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Technical Information for R 60 / R 60.1 keyless chuck

Observe the safety instructions of the drive unit!


Safety instructions


- Disconnect the mains plug before securing R 60 / R 60.1 keyless chuck and stirring element to the stirrer.
- Check R 60 / R 60.1 keyless chuck and stirring element is held in position securely prior to each use and also at regular intervals. The chuck and stirring element must only be tightened and changed when the equipment is stationary and the power supply is disconnected.
- Please note and observe the operating instructions of the used overhead stirrer.
- R 60 / R 60.1 keyless chuck can only be used for overhead stirrers with max. torque of the stirring shaft less than 200 Ncm.

Assembly

Securing R 60 / R 60.1 keyless chuck to overhead stirrer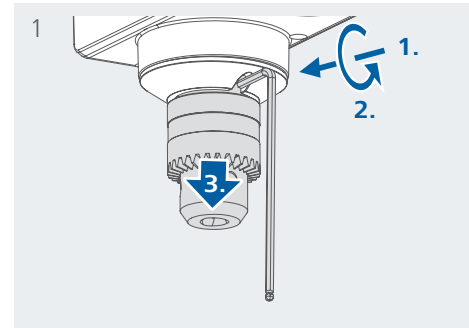


Disassemble the included chuck from overhead stirrer.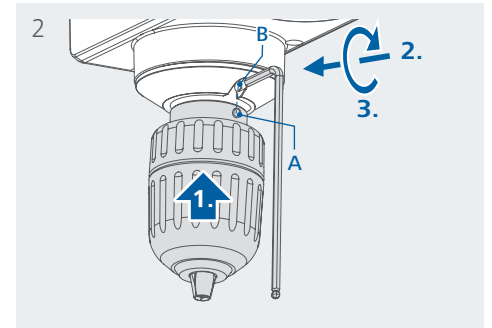

Slide R 60 / R 60.1 keyless chuck along the output shaft and align the hexagonal socket screw (A) with the hole (B). Tighten the hexagonal socket screw using the cranked wrench. Until at last hexagonal socket screw is screwed into the screw hole completely.

Note: before installing the chuck please loose the hexagonal socket screw (A) firstly so that the output shaft could be inserted into.


Securing the stirring element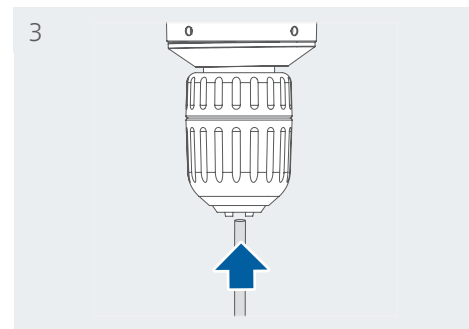

Slide the stirring tool into the chuck.

Note: before attaching the stirring tool please loose the R 60 / R 60.1 keyless chuck firstly so that the stirring tool can be inserted into.

Technical specifications may be changed without prior notice.

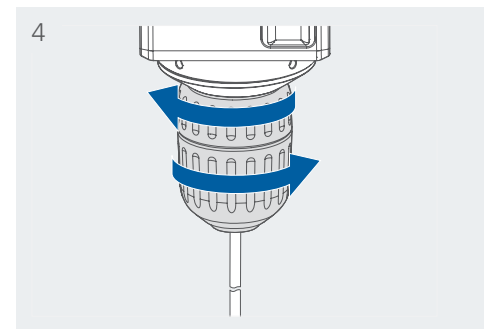Grip the upper and lower part of R 60 / R 60.1 keyless chuck with hands and tighten the chuck according to above shown direction until stop position. For loosening the R 60 / R 60.1 keyless chuck, turn the upper and lower part in reverse direction.

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