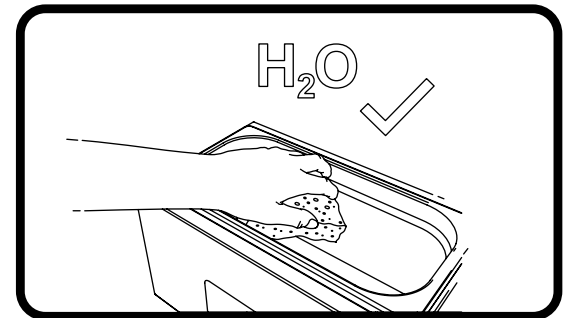
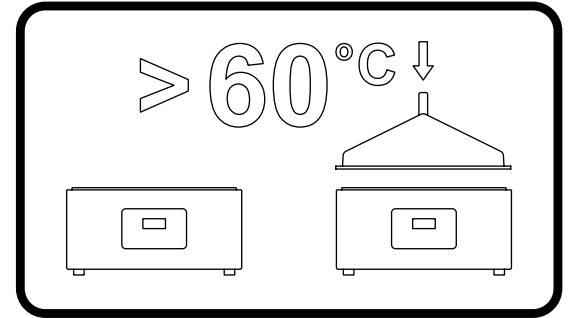
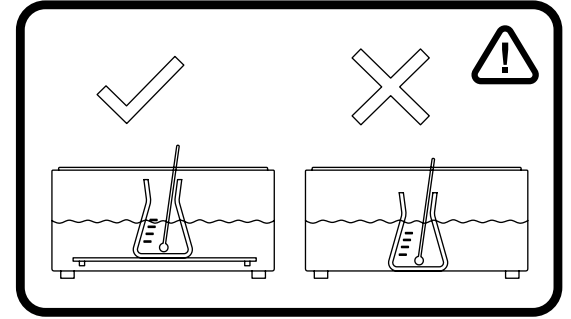
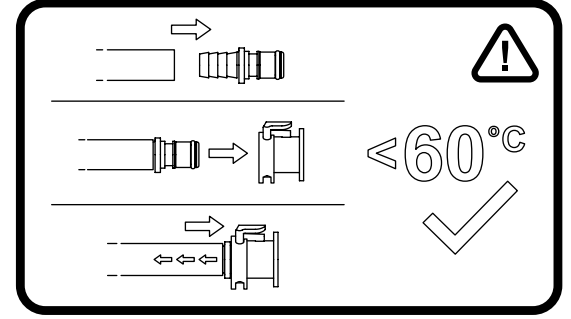
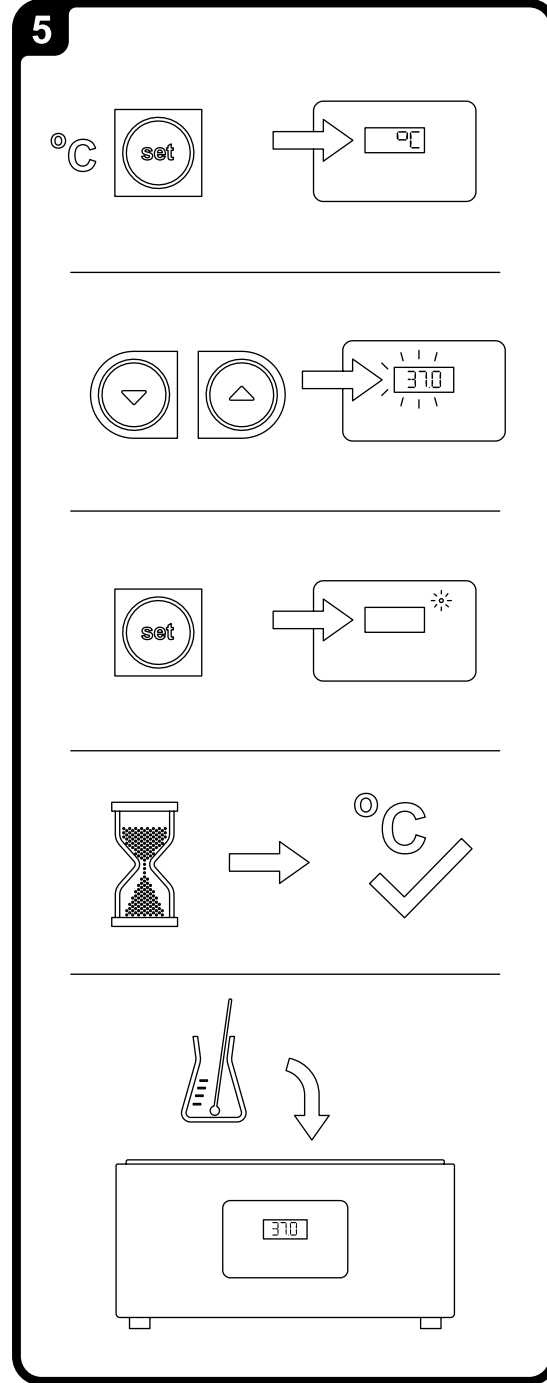
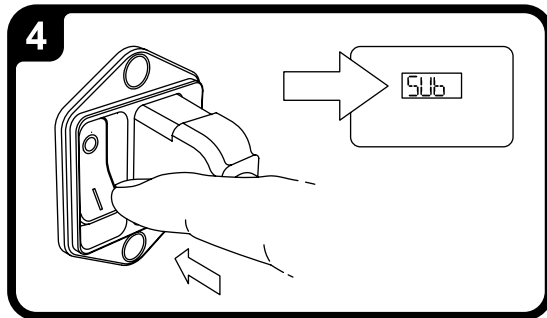
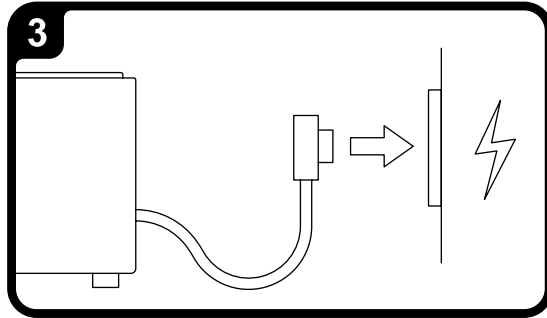
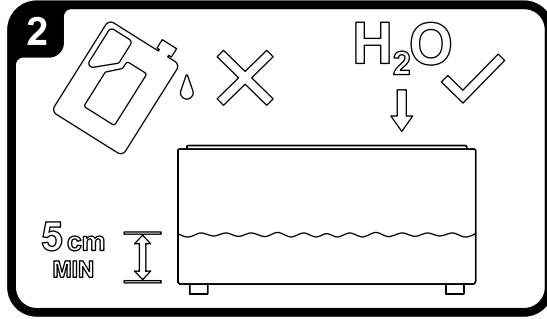
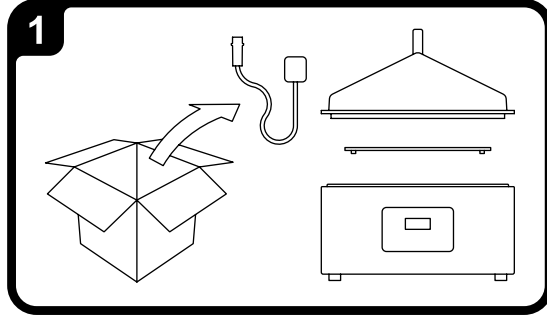
Safety information

GB

FR

DE

ES

IT

This product is a general purpose thermostatically controlled water bath designed for indoor laboratory use by or supervised by a professional user.

Read the whole of the instructions. Safety may be impaired if they are not followed.

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Surfaces and water can be hot during and after use.

This bath is only intended for use with water. Use of other fluids may invalidate warranty and present a risk of fire or explosion.

Place on a stable flat surface to reduce the risk of accidental spillage.

No user serviceable parts. Risk of electric shock after disassembly or operation with covers removed.

Not for use in environments with a risk of flammable or explosive gases. To be operated within the limits listed below.

Only use the mains cord provided. Ensure that the mains plug and the switch are easily accessible.

A clearance of >10cm around the bath is required to ensure adequate air flow.

For optimum performance prevent tray from touching bath sides.

If a potentially hazardous liquid is spilt onto the equipment, disconnect it from the power supply and have it checked by a competent person. It is the user's responsibility to carry out appropriate decontamination if hazardous material is spilt on the equipment.

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To preserve your water bath in peak condition consult the extra guidance listed below. Failure to do so may affect your warranty. Consult online resources for additional important information.

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Online resources

How to videos

- Care and maintenance of your bath
- Calibrating your bath
- Configuring advanced features

Full user manuals available for download at:
www.grantinstruments.com

Warranty

When used in accordance with the user instructions this product is guaranteed for 3 years against faulty materials and workman ship
Register your warranty online
<http://www.grantinstruments.com/support/warranty-registration/>

Specification

Suitable for use with water and heat transfer beads (excluding SAP2 & SAP2S)

Mains supply:
220-230V@50/60Hz
110-120V@50/60Hz
Earth connection required

Ambient operating temp:
+5°C to 40°C
Ambient storage temp:
+20°C to +60°C

Current consumption

	230v	120v
SAP2	1.0A	2.0A
SAP2S	1.5A	3.0A
SAP5	1.5A	3.0A
SAP12	3.5A	6.5A
SAP18	6.0A	8.5A
SAP26	6.0A	8.5A
SAP34	8.0A	11.0A
SAPD	5.0A	9.0A