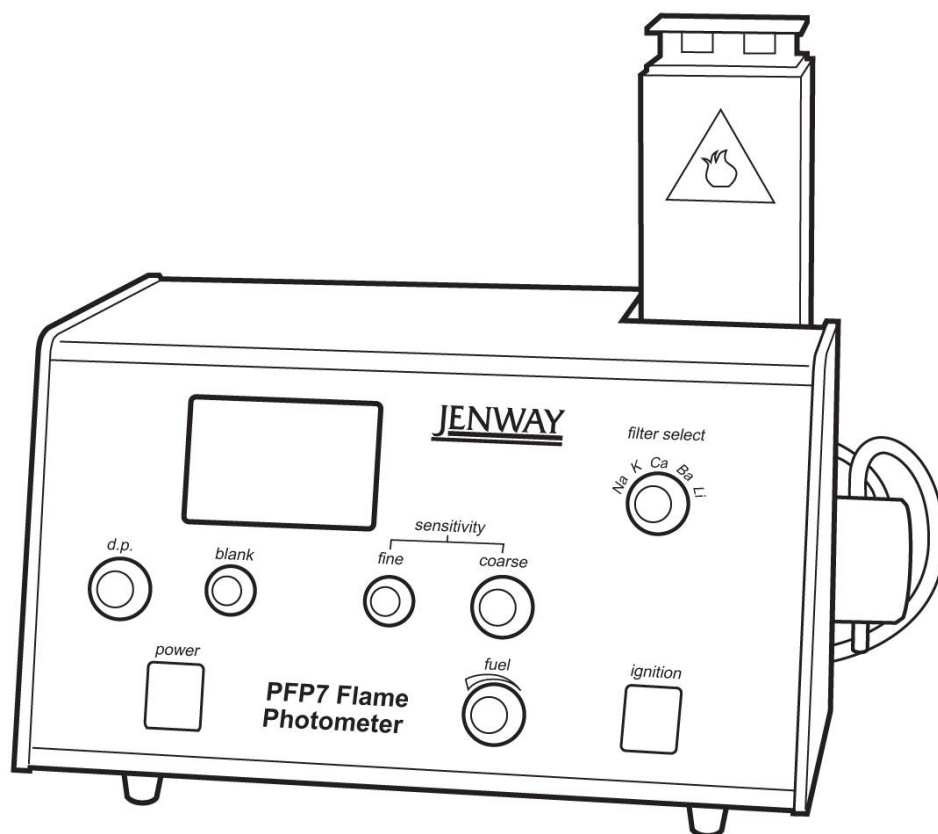
# Flame Photometer

Models PFP7 and PFP7/C

## Operating Manual

# English

## Safety Information

Please read this information carefully prior to installing or using this equipment.

1. The unit described in this manual is designed to be operated only by trained personnel. Any adjustments, maintenance and repair must be carried out as defined in this manual, by a person qualified to be aware of the hazards involved.
2. It is essential that both operating and service personnel employ a safe system of work, in addition to the detailed instructions specified in this manual.
3. Other than for those items defined in the maintenance procedures herein there are no user serviceable items in this instrument. Removal of covers and attempted adjustment or service by unqualified personnel will invalidate the warranty and may incur additional charges for repair.
4. References should always be made to the Health and Safety data supplied with any chemicals used. Generally accepted laboratory procedures for safe handling of chemicals should be employed.
5. If it is suspected that safety of the unit has been impaired in any way, the unit must be made inoperative and secured against any intended operation. The fault condition should immediately be reported to the appropriate servicing authority.

**NOTE:** Failure to use this equipment in accordance with the manufacturer's manual may compromise your basic safety protection afforded by the equipment and may invalidate the warranty. The warranty does not cover damage caused by faulty installation or misuse of the equipment.

### Prevention of Fire and Electric Shock.

To prevent a risk of fire or electric shock, **DO NOT** open your product case without authorisation. Only qualified service personnel should attempt to repair this product.

Replace fuses only with the type as listed in section 4.1.2.

Ensure the mains power supply conforms to the rating found on the serial plate located on the back of this product.

**Never** operate this equipment without connection to earth / ground. Ensure the mains supply voltage is correctly earthed / grounded in accordance with current area legislation.

### General Safe Operating Practice

Always follow good laboratory practice when using this equipment. Give due recognition to your company's safety and legislative health & safety procedures and all associated legislation applicable to your areas of operation. Check laboratory procedures for substances being used and ensure all hazards (e.g. explosion, implosion or the release of toxic or flammable gases) that might arise have been suitably addressed before proceeding. When using certain

substances the liberation of hazardous gases may require the use of a fume cupboard or other means of extraction.

Be aware of biological contamination. Ensure sterilization of the capillary tube, waste, nebuliser and mixing chamber is performed after contamination.

Ensure equipment is used on a clean, dry, non-combustible, solid work surface with at least 300mm suitable clearance all around and at least 1m clearance above the chimney.

Ensure that the outer chimney is properly installed before operating the equipment.

**Attention:** Ensure that the gas and air pipes have been connected correctly and that there are no gas leaks present before igniting.

Ensure that the unit is properly connected to the gas cylinder making sure there are no gas leaks. Ensure that the correct gas regulator is used and connected properly. The responsible body must ensure that the type of connector used on the outlet of the gas-pressure regulator conforms to the applicable national requirements.

Take care when fitting and removing the capillary tube to the nebulizer needle.

It is not recommended to leave the flame photometer unattended during operation.

**Do not** operate or handle any part of the product with wet hands.

**Do not** look down the flame chimney when igniting or operating the equipment.

**Do not** handle the flame chimney whilst in operation.

**Do not** lean or stretch over the equipment whilst in operation.

**Do not** cover the chimney whilst in use. **Do not** block or obstruct ventilation slots / airways.

**Do not** spill substances onto the unit. If spillage does occur, disconnect unit from mains supply. If using hazardous substances please adhere to standard laboratory operating procedures.

**Do not** position the product so that it is difficult to disconnect from the mains supply.

**Do not** move or carry the unit when in use or connected to the mains electricity supply.

**Important:** Allow sufficient time for the chimney to cool before handling.

**Do not** use acetylene with the flame photometer.

**Warning:** Over adjustment of the fuel valve will cause excess flame.

Avoid installing the flame photometer in areas that are susceptible to drafts, but ensure that there is adequate ventilation to prevent any build-up of gas.

Before using any cleaning or decontamination methods except those specified by the manufacturer, **responsible bodies** should check with the manufacturer that the proposed method will not damage the equipment.

In order to completely disconnect power to the flame photometer the mains power cord must be detached from the back of the unit.

Ensure a suitably rated detachable mains cord is used.

This hazard symbol is present on the flame chimney outer casing and indicates that the surface may be **hot to touch**.

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### Introduction

#### 1.1 Instrument description

The PFP7 and PFP7/C are low temperature, single channel emission flame photometers designed for the routine determination of sodium (Na), potassium (K) and lithium (Li). Additional filters for the determination of calcium (Ca) and barium (Ba) are supplied with the PFP7. Both versions are fitted with an automatic flame failure detection for user safety, making them ideal for use in clinical, industrial or educational applications. Model PFP7/C is specifically designed for use in clinical applications. The in-built lineariser circuitry enables readings of both Na and K, at normal clinical serum concentrations to be displayed directly in mmol/l. Serum samples must be diluted 200:1 or 100:1 prior to presentation to the flame photometer.

#### 1.2 Specification

|                       | PFP7                                                                                                                                         |         | PFP7/C                                 |         |
|-----------------------|----------------------------------------------------------------------------------------------------------------------------------------------|---------|----------------------------------------|---------|
| Ranges                | 0 to 199.9ppm                                                                                                                                |         | 120-160 mmol/L Na (linearized)         |         |
|                       |                                                                                                                                              |         | 0-10.0 mmol/L K                        |         |
| Limits of Detection   | Na                                                                                                                                           | 0.2ppm  | Na                                     | 0.2ppm  |
|                       | K                                                                                                                                            | 0.2ppm  | K                                      | 0.2ppm  |
|                       | Li                                                                                                                                           | 0.25ppm | Li                                     | 0.25ppm |
|                       | Ca                                                                                                                                           | 15ppm   | -                                      |         |
|                       | Ba                                                                                                                                           | 30ppm   | -                                      |         |
| Reproducibility       | 1% Coefficient of variation (C.V.) for 20 consecutive samples using 10ppm Na set to read at 50.0 (PFP7) or 140 (PFP7/C).                     |         |                                        |         |
| Linearity             | Better than 2% when concentration of 3ppm Na and K and 5ppm Li are set to read 100.                                                          |         |                                        |         |
| Specificity           | Interference from Na, K and Li when equal in concentration to the test element will be less than 0.5%                                        |         |                                        |         |
| Stability             | Better than 2% over 5 minutes when continuously aspirating 10ppm, sample set to read 50.0. Zero drift better than 2% per hour after warm-up. |         |                                        |         |
| Sample requirements   | Between 2 and 6 ml/minute                                                                                                                    |         |                                        |         |
| Recorder Output       | Nominal 1.00 volt for readout of 100.0.                                                                                                      |         |                                        |         |
| Warm Up               | The flame must be alight for at least 15 minutes to ensure achievement of the above stated specification.                                    |         |                                        |         |
| Services              | Electrical:                                                                                                                                  |         | 90-125V or 190-250V @ 50/60Hz          |         |
|                       | Air:                                                                                                                                         |         | Moisture and oil-free.                 |         |
|                       | Fuel:                                                                                                                                        |         | Propane, butane, natural gas or L.P.G. |         |
| Operating Environment | 15°C to 35°C                                                                                                                                 |         |                                        |         |
| Size                  | 420 x 360 x 300mm                                                                                                                            |         |                                        |         |
| Weight                | 8kg                                                                                                                                          |         |                                        |         |

### 1.3 Principles of operation

Flame photometry relies upon the fact that the compounds of the alkali and alkaline earth metals can be thermally dissociated in a flame and that some of the atoms produced will be further excited to a higher energy level. When these atoms return to the ground state they emit radiation which lies mainly in the visible region of the spectrum. Each element will emit radiation at a wavelength specific for that element. The table below gives details of the measurable atomic flame emissions of the alkali and alkaline earth metals in terms of the emission wavelength and the colour produced.

| Element       | Emission Wavelength (nm) | Flame Colour |
|---------------|--------------------------|--------------|
| Sodium (Na)   | 589                      | Yellow       |
| Potassium (K) | 766                      | Violet       |
| Barium (Ba)   | 554                      | Lime Green   |
| Calcium (Ca)  | 622*                     | Orange       |
| Lithium (Li)  | 670                      | Red          |


\*Note: Calcium is measured by using the calcium hydroxide band emission at 622nm as the Calcium main atomic emission occurs at 423nm.

Over certain ranges of concentration the intensity of the emission is directly proportional to the number of atoms returning to the ground state. This is in turn proportional to the absolute quantity of the species volatilized in the flame, i.e. light emitted is proportional to sample concentration.

It can be seen that if the light emitted by the element at the characteristic wavelength is isolated by an optical filter and the intensity of that light measured by a photo-detector, then an electrical signal can be obtained proportional to sample concentration. Such an electrical signal can be processed and the readout obtained in a digital form.

A simple flame photometer consists of the following basic components:

- The burner: a flame that can be maintained in a constant form and at a constant temperature.
- Nebuliser and mixing chamber: a means of transporting a homogeneous solution into the flame at a steady rate.
- Simple colour filters (interference type): a means of isolating light of the wavelength to be measured from that of extraneous emissions.
- Photo-detector: a means of measuring the intensity of radiation emitted by the flame.

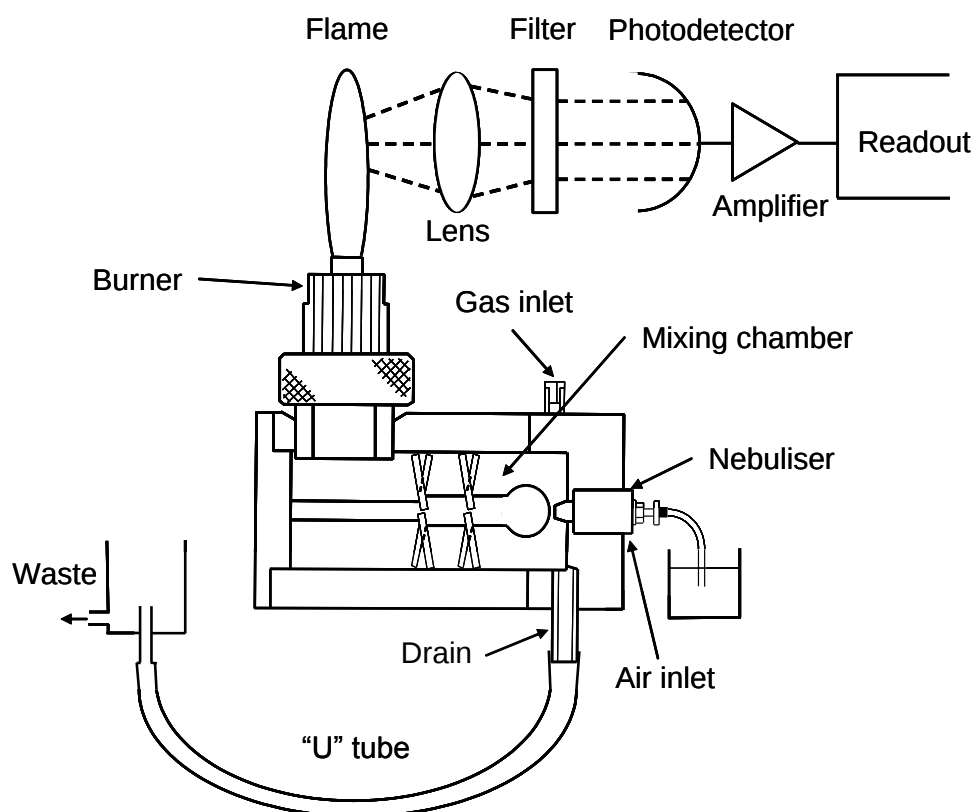**Figure 1.3.1:** Schematic diagram showing the component parts of a flame photometer.

The analysis of alkali and alkaline earth metals by flame photometry has two major advantages:

- i. Their atoms reach the excited state at a temperature lower than that at which most other elements are excited.
- ii. Their characteristic wavelengths are easily isolated from those of most other elements due to wide spectral separation.

The analysis of Na, K, Li, Ba and Ca are typically determined at low temperatures, i.e. 1500-2000°C, therefore suitable fuel mixtures are propane/air, butane/air and natural gas/air.

## 1.4 Good practice guidelines

1. It is most important that the nebuliser, mixing chamber and burner are kept clean by carrying out the correct shutdown procedure and by periodic maintenance. If high salt solutions are aspirated, correspondingly longer periods should be spent aspirating deionised water prior to shut down.
2. It is recommended that blank and standard solutions should have a wetting agent (e.g. Triton X-100<sup>1</sup> or Decon 90) added to promote good stability and self-cleaning. Any such wetting agent should be non-ionic and used at a concentration of less than 3ppm. It should be added to the blank, standards and samples at the same concentration.
3. Take care when preparing standards. The performance of the instrument depends upon the accuracy and purity of the calibration standards.
4. If standard solutions are required to be stored for any length of time or at an elevated temperature, a suitable mould inhibitor e.g. azide should be added. However if this contains the element to be measured (e.g. sodium) it is important that the samples also contain an equivalent amount.
5. Always sample from the top half of the sample container. The bottom half may contain sediment or particulate matter which could easily block the fine tubing used in the nebuliser.
6. Always use recommended spares. Even where an alternative part may be obviously suitable there may be good reasons for not using it.
7. Never use glass containers to store calibration standards.

---

<sup>1</sup> Triton X-100 is a registered trademark of Union Carbide Chemicals and Plastics Co. Inc.

## Section 2

### Installation

#### 2.1 Services required

The fuel and air supplies to the instrument must be clean and dry and supply pressures regulated within the limits specified. Any contamination, moisture or variation in supply pressure will directly affect the performance of the instrument.

**NOTE: The instrument will only operate with the correct type of air compressor and gas regulator.**

##### 2.1.1 Voltage

This unit is designed to operate at 90 - 250V and at 50 or 60Hz.

##### 2.1.2 Fuel

- Propane or butane regulated to 10-15 inches water gauge (0.36-0.54psi or 0.025-0.038kg/cm<sup>2</sup>).
- Natural gas at mains pressure between 3 and 10 inches water gauge (0.11-0.36psi or 0.0076-0.025kg/cm<sup>2</sup>).
- L.P.G. regulated to 10-15 inches water gauge (0.36-0.54psi or 0.025-0.038kg/cm<sup>2</sup>).
- Suitable regulators are available from Jenway; refer to Section 2.5.

##### 2.1.3 Air

A supply of dry, clean and pulse-free air at a pressure at 15 - 30psi (approx. 1 - 2 kg/cm<sup>2</sup>) at 6 litres/minute is required. A suitable compressor and water separator are available from Jenway; refer to Section 2.5.

##### 2.1.4 Drain

The instrument will need to be sited near a drain or sink to enable disposal of waste liquid. A suitable receptacle can be utilised if provision is made for easy disposal of its contents. **WARNING:** The waste liquid will still contain any hazardous materials that were in the original samples and should be handled and disposed of with the same care. Waste liquid should always be considered to be of a pathogenic nature where the instrument is used in a clinical environment.

## 2.2 Unpacking

Remove the instrument from the packaging and ensure the following items are present:


| Part                                                  | Part Code          |
|-------------------------------------------------------|--------------------|
| Model PFP7 or PFP7/C Flame photometer                 | 500 701 or 500 801 |
| Auxiliary power plug                                  | 009 035            |
| Mains cable with connector (plug optional as ordered) | 013 046            |
| Nebuliser inlet tube (500mm)                          | 500 193            |
| Gas tube, low pressure (2 metres)                     | 500 191            |
| Silicon rubber tube (drain)                           | 023 003            |
| Air tube (2 metres)                                   | 500 192            |
| Nozzle 1: gas                                         | 026 012            |
| Nozzle 2: air                                         | 026 013            |
| Allen key, 2.5mm                                      | 060 037            |
| Adjustable hose clip                                  | 060 083            |
| Drain trap                                            | 500 018            |
| Drain trap clip                                       | 500 114            |
| Nebuliser cleaning wire                               | 500 194            |

**Any shortages or damage should be reported immediately to the manufacturer or local distributor.**

## 2.3 Mixing Chamber Assembly


Place the instrument on a flat, stable bench and proceed as follows (see **Figures 2.3.1** and **2.3.2** below):

1. Take the mixing chamber assembly and ensure that the fluted burner (1) is in place (remove the retaining tape). Place the mixing chamber and burner assembly inside of the flame photometer and line up the burner with the bottom of the chimney. Look down the chimney and ensure that the burner locates centrally in the inner chimney. Screw the locking ring (2) onto the threaded boss provided and tighten until finger tight, ensuring that the mixing chamber is at right angles to the side of the instrument.
2. Fit the nebuliser (3) into the end cap (8) and connect the air tubing to the nebuliser inlet (4).
3. If natural gas is to be used, remove the restrictor screw (5) fitted into the end of the fuel inlet (6) on the mixing chamber. The restrictor screw should remain in place if butane, propane or L.P.G. is to be used as fuel. Keep the restrictor screw in case of any future change in gas supply. Push the fuel tubing onto the fuel inlet connector (6) – this tube can be secured with a metal hose clip.

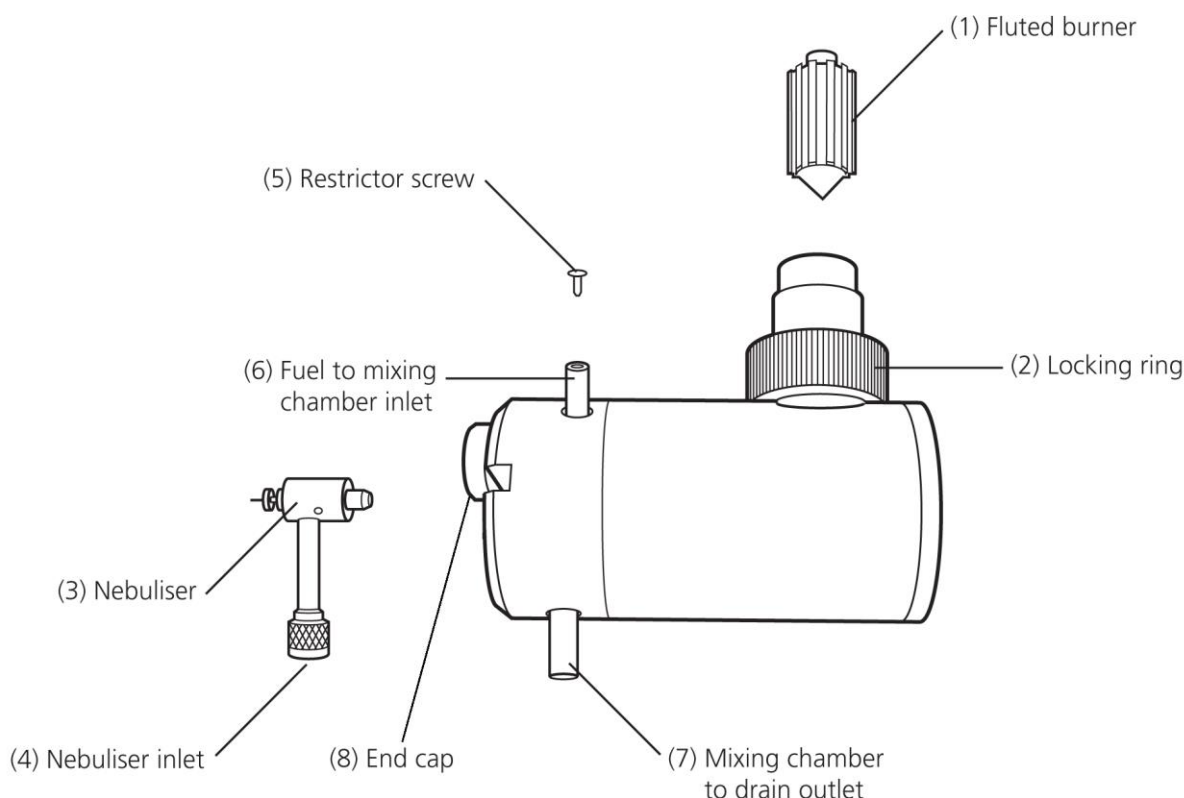**Figure 2.3.1:** Assembly of the mixing chamber, burner and nebuliser.

4. Take the drain trap clip (9) and, with the screw provided, fix to the hole situated at the chimney end of the rear panel. Fix a length of silicon rubber tubing (10) onto the side port on the drain trap such that it is long enough to reach either the sink or the waste receptacle to be used. Fit the drain trap into the clip on the rear panel and run the short piece of tubing (11) from the bottom of the drain trap to the mixing chamber waste outlet (7) on the bottom of the mixing chamber. This tube can be pushed on and does not need securing.

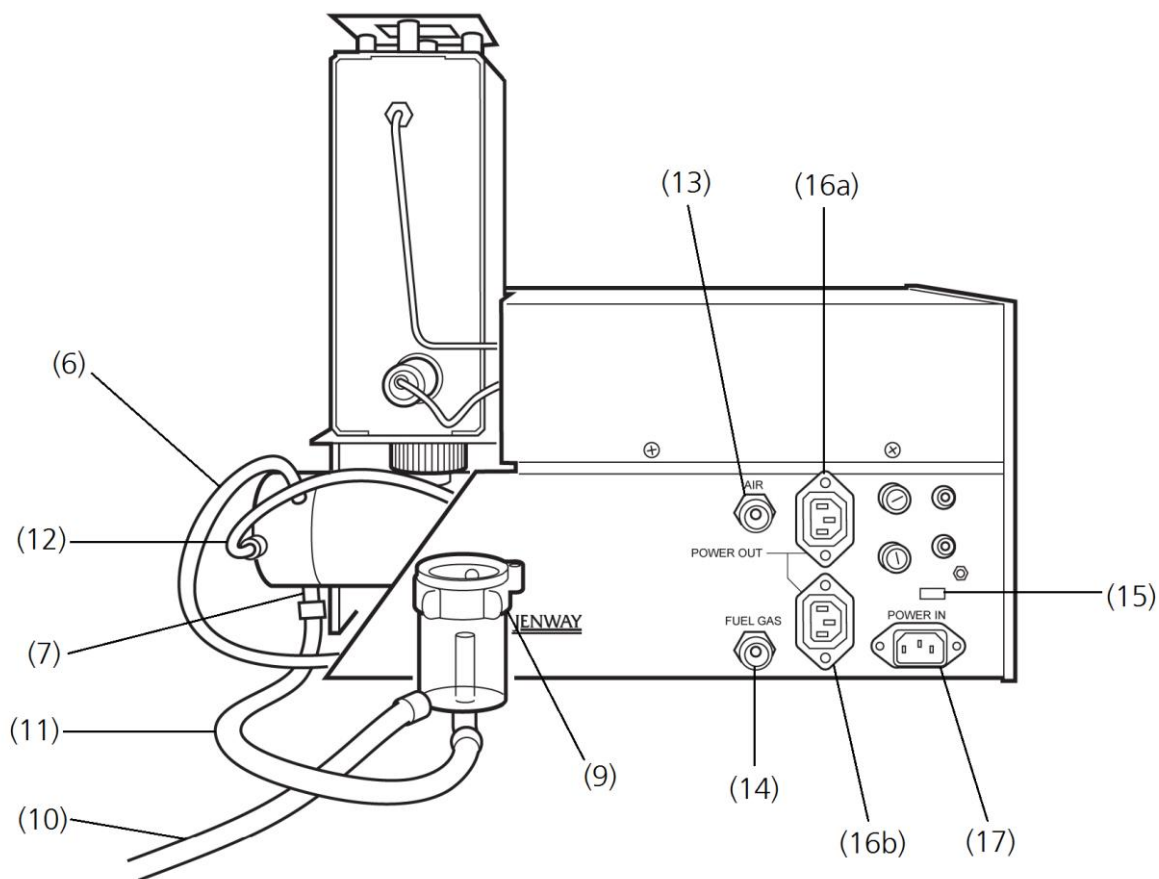**Figure 2.3.2:** Rear panel showing the drain trap, air and fuel inlets.

- 1) Fluted burner
- 2) Locking ring
- 3) Nebuliser
- 4) Nebuliser inlet
- 5) Restrictor screw
- 6) Fuel to mixing chamber inlet
- 7) Mixing chamber to drain outlet
- 8) End cap
- 9) Drain trap clip
- 10) Drain trap to waste container tubing
- 11) Mixing chamber to drain trap tubing
- 12) Compressed air to nebuliser tubing
- 13) Compressed air input connector
- 14) Fuel gas input connector
- 15) Adjustable voltage fuse
- 16) POWER OUT sockets
- 17) POWER IN socket

## 2.4 Installation

### WARNING

The exhaust gases from the chimney are very hot. No obstruction should be placed above the instrument and the instrument should be located in a position that makes accidental contact with the chimney or its exhaust unlikely.

The instrument needs to be operated in a well-ventilated room (although very strong draughts should be avoided).

Ensure the air tubing is routed away from sources of heat.

Fit the air and fuel inlet connectors to the ports provided on the rear of the instrument.

1. Fit the appropriate lengths of 8mm and 6mm tubing to the connectors 13 and 14 (figure 2.3.2). Run the tubing to the air and fuel supplies to be used.

**NOTE:** The maximum inlet pressure and regulation requirements are defined in Section 2.1.

Turn on the fuel supply at the cylinder or source and check all fuel joints for leaks by using soap solution or suitable gas detector.

2. Plug the power cable into the POWER IN socket 17 (figure 2.3.2). If necessary connect a suitable plug to the 3-pin AC supply lead. The colours of the wires conform to the internationally recognised standard such that:

|                     |                |
|---------------------|----------------|
| <b>BROWN</b>        | <b>LIVE</b>    |
| <b>BLUE</b>         | <b>NEUTRAL</b> |
| <b>GREEN/YELLOW</b> | <b>EARTH</b>   |

**IMPORTANT – THE INSTRUMENT MUST BE EARTHED.**


The green/yellow wire in the AC supply must be connected to a properly grounded terminal.

If a compressor is to be run directly from the instrument then it should be connected to the plugs provided and connected to the POWER OUT sockets 16a and 16b (figure 2.3.2).

Press the power switch on the front panel. The digital display should be illuminated.

3. Fill the centre tube of the drain trap and the connecting silicon rubber tubing to the mixing chamber with deionised water. Check that no air bubbles are trapped in this tube and that it flows and runs to waste freely. Ensure that the drain trap is pushed completely down on its clip.

When the mixing chamber and drain trap are fully connected your flame photometer should look similar to the illustration below:

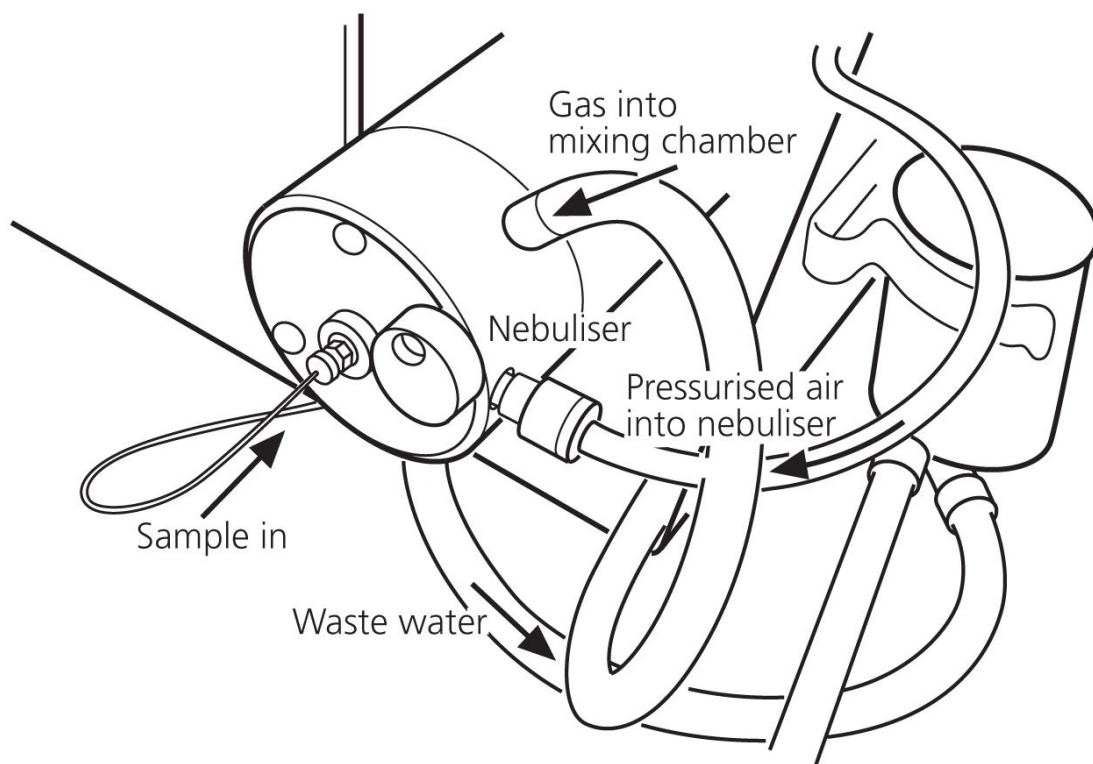**Figure 2.4.1:** An alternative view of the mixing chamber showing inlet and outlet tubing

## 2.5 Accessories

Always use recommended spares and accessories. Even if an alternative part may appear obviously suitable there may be some minor variations in specification that could degrade the performance of the instrument.

| Accessories                           | Part Code   |
|---------------------------------------|-------------|
| Model 8515 air compressor (220V 50Hz) | 535 001     |
| Model 8516 air compressor (110V 60Hz) | 535 002     |
| Butane regulator                      | 500 178     |
| Propane regulator                     | 500 179     |
| Natural gas regulator                 | 500 180     |
| Water separator (small)               | 500 176     |
| Water separator (large)               | 500 177     |
| Dust cover                            | 500 134     |
| Cleaning solution (1 litre)           | 025 171     |
| Minor spares kit                      | PFP7/MINKIT |

## Section 3

### Analysis Preparation

#### 3.1 Calibration standards

A comprehensive range of aqueous calibration standards is available from Jenway. These must be diluted to a suitable concentration for aspiration into the flame (see Sections 3.3 and 4.3)

| Industrial Standards (500ml) | Part Code |
|------------------------------|-----------|
| 1000ppm K                    | 025 023   |
| 1000ppm Li                   | 025 024   |
| 3000ppm Ba                   | 025 025   |
| 1000ppm Na                   | 025 021   |
| 1000ppm Ca                   | 025 009   |

When preparing standards always observe the following:

1. Standards must always contain the constituents that are present in the samples in the same concentration ratios; i.e. if samples are prepared in 0.05M HCl then the standards should also contain 0.05M HCl.
2. Always ensure that the standards encompass the expected range of the sample concentrations.
3. Standards should be prepared so as to ensure that the region in which measurements are made coincide with the concentrations that produce the optimum performance from the flame photometer, i.e...
  - ...when measuring sodium, the top standard is ideally 10ppm,
  - ...when measuring potassium, the top standard is ideally 10ppm,
  - ...when measuring calcium, the top standard is ideally 100ppm,
  - ...when measuring barium, the top standard is ideally 1000ppm,
  - ...when measuring lithium, the top standard is ideally 10ppm.

NOTE: The blank used should contain all the constituents of the standard solutions except the element being measured.

Since a flame photometer measures the concentration of the element itself in solution, standard solutions prepared from the salts of sodium, potassium, lithium, calcium and barium must be made up to contain the concentrations required in terms of the quantity of the elements.

Below are two examples of how to prepare standards of 1mg Na/100ml (10ppm Na) and 1mg K/100ml (10ppm K).

### 3.1.1 Sodium

Accurately weigh 0.634g of dry “Analar” quality NaCl, dissolve in pure deionised water and wash into a 500ml volumetric flask. Fill to the mark with pure deionised water. To prepare the standard solution for use with the flame photometer, this stock solution should be diluted 1 in 50.

Calculation:

Atomic weight Na = 23.0

Molecular weight NaCl = 58.46

Therefore, 0.634g NaCl contains

$$\frac{0.634 \times 23}{58.46} = 0.25 \text{g Na}$$

Thus in 500ml of solution there is 250mg Na or 50mg Na/100ml.

Diluting 1 in 50 gives a standard of 1mg Na/100ml = 10ppm Na.

### 3.1.2 Potassium

Accurately weigh 0.477g of dry “Analar” quality KCl, dissolve in pure deionised water and wash into a 500ml volumetric flask. Fill to the mark with pure deionised water. To prepare the standard solution for use with the flame photometer, this stock solution should be diluted 1 in 50.

Calculation:

Atomic weight K = 39.1

Molecular weight KCl = 74.56

Therefore, 0.477g KCl contains

$$\frac{0.477 \times 39.1}{74.56} = 0.25 \text{g K}$$

Thus in 500ml of solution there is 250mg K or 50mg K/100ml.

Diluting 1 in 50 gives a standard of 1mg K/100ml = 10ppm K.

### 3.1.3 Storage

Store solutions away from direct sunlight in a cool place, ideally at temperatures below 25°C. Glass containers should not be used for storage as they can affect the sodium concentration levels. Standards should be stored in sealed, plastic vessels and in high concentrations, (e.g. as a stock 1000ppm solution) and dilutions prepared as required. The long-term storage of low concentration standards is not recommended due to degradation of ionic species.

## 3.2 Sample preparation

There are several practical points regarding sample preparation, which should be adhered to in order to achieve the required accuracy in your analysis:

1. Avoid handling samples with fingers. This leads to serious contamination, e.g. if a finger is immersed in 20ml of deionised water the resulting Na concentration will exceed that of a 10ppm standard.
2. All analyses involve the use of a diluent, which is almost always deionised water. This should be of the highest quality for accurate flame analysis. Sodium, potassium and calcium are present in high concentrations in tap water and thus efficient deionisation is essential.
3. Species that cause interference should be removed from samples or the equivalent concentration of the interfering ion should be present in the standards so as to avoid erroneous results, e.g. if a sample of approximately 10ppm Na contains approximately 1000ppm Ca, then Na analysis can only be achieved by removing the Ca with oxalate/oxalic acid or ensuring all standards contain 1000ppm Ca.
4. Always try to follow a well-documented analytical procedure, which should contain information pertaining to interference removal when applicable.
5. Standards and samples should not be exposed to the atmosphere for long periods due to contamination from airborne particles and the evaporation of the solvent that could lead to elevated concentrations.
6. When in doubt about the equipment or application, the operator should contact Jenway at [jenwayhelp@bibby-scientific.com](mailto:jenwayhelp@bibby-scientific.com) for advice.

### 3.2.1 Sample extraction

A number of methods for extracting sodium, potassium, lithium, calcium and barium from a wide variety of raw materials may be obtained by contacting the technical support helpline at [jenwayhelp@bibby-scientific.com](mailto:jenwayhelp@bibby-scientific.com) or by visiting [http://jenway.com/notes\\_Flame\\_Photometers.asp](http://jenway.com/notes_Flame_Photometers.asp)

The sample must be in the form of an aqueous solution, with no solid matter present, to be suitable for direct introduction into the flame photometer. This is achieved by:

- Extracting the salts from solid samples using deionised water or suitable extractants e.g. saturated  $\text{CaSO}_4$  for sodium in soil. Extraction is more successful using a blender, macerator or shaking machine.
- If the sample is organic then the organic material should be removed by ashing. The remaining oxides are then dissolved using strong acids.
- Filtration/centrifugation is used to remove solid debris.

When aqueous, the sample can then be diluted to a known, accurately measured volume using deionised water. If it is a concentrated sample then the dilution ratio should be increased. If the sample concentration is low then a small volume of diluent and initial extractant should be used. Whichever method of extraction is used, the resultant solution must always be free of any particulate matter that may cause blockages in the nebuliser capillary tube.

### **3.3 Dilution**

In order to obtain samples and standards of the right concentration for aspiration into the flame, various levels of dilution will often be necessary. Good quality deionised water should normally be used for carrying out these dilutions and it is recommended that the same batch of water should be used for diluting the samples and standards. More information on typical dilution ratios is given in Section 4.3.

## Section 4

### Operation

#### 4.1 Front and rear panel controls

##### 4.1.1 Front panel controls


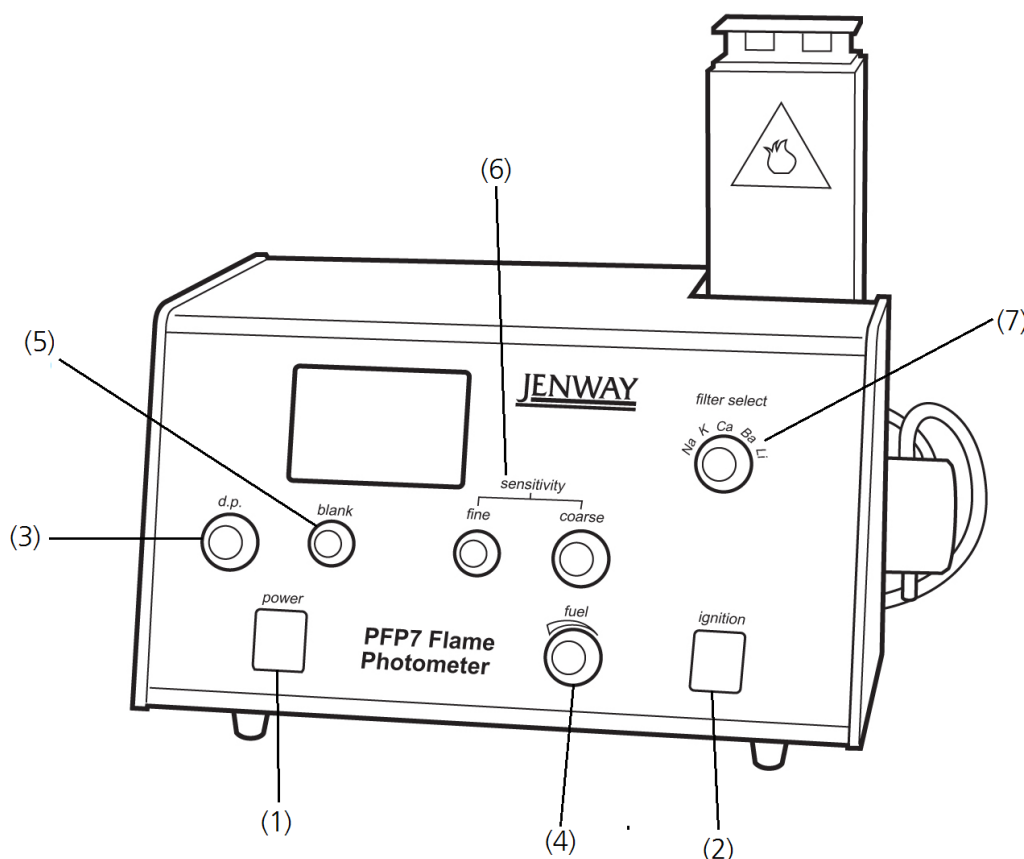
Figure 4.1.1.1: Front panel

- |                    |                                                                                                                                                                                                                                |
|--------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>1. Power</b>    | A two-position rocker switch which controls the AC supply of the instrument. Any accessories connected to the auxiliary <b>POWER OUT</b> sockets on the rear panel are also controlled by the front panel <b>power</b> switch. |
| <b>2. Ignition</b> | A spring loaded switch which, when depressed, will cause an electrical discharge between the ignition electrode and the burner unit, thereby causing fuel ignition.                                                            |
| <b>3. D.P.</b>     | This dial controls the position of the decimal point on the digital readout.                                                                                                                                                   |
| <b>4. Fuel</b>     | A fine needle valve that controls the flow of fuel and enables optimum flame conditions to be set.                                                                                                                             |
| <b>5. Blank</b>    | This dial control sets the zero point when a blank standard is aspirated.                                                                                                                                                      |

6. **Sensitivity:**  
**Fine and Course** Two controls which are used in conjunction to set the digital readout to an appropriate number when a calibration standard is aspirated.
7. **Filter Select** A five position control which will select the appropriate optical filter for the element being determined.

**NOTE:** For the PFP7/C positions 3 and 4 will normally be blank

#### 4.1.2 Rear panel controls

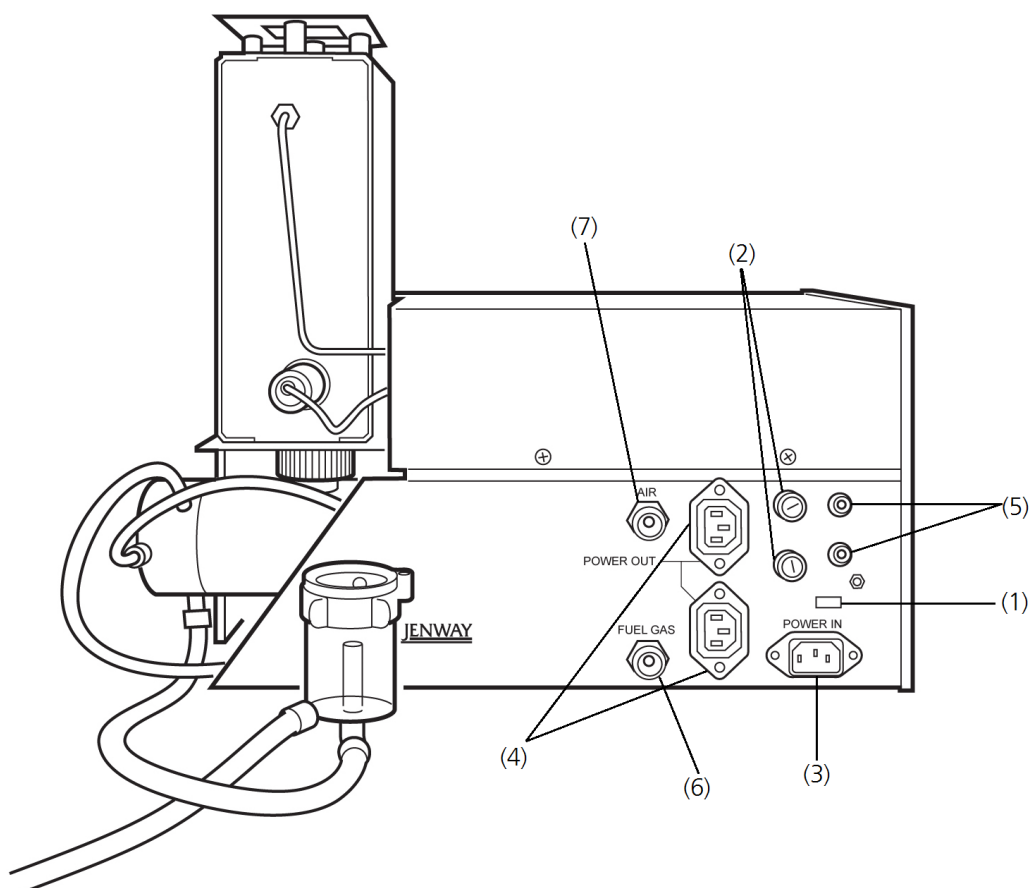**Figure 4.1.2.1:** Rear panel

1. **VOLTAGE SELECTION** A two-position switch marked 230 and 115. These positions allow operation from the voltage supplies 190-250 volts and 90-125 volts respectively at either 50 or 60Hz.  
**WARNING:** when adjusting the operating voltage it may be necessary to change the top fuse as indicated on the rear panel. Please note that the unit is set to the required voltage in the factory depending on the part code ordered.

2. **FUSES** Two fuses are fitted to the instrument. **FS1** is the primary fuse provided to protect the instrument and any accessories connected to the **power out** sockets. **FS2** is to protect the electronic circuitry in the PFP7. Both FS1 and FS2 should be of an anti-surge type.
3. **POWER IN** A three pin receptacle for the AC mains supply.
4. **POWER OUT** Two three pin sockets for flame photometer accessories such as a compressor. The power from these sockets is controlled by the instrument **power** switch on the front panel.
5. **RECORDER** Two 4mm sockets which will provide an analogue signal of approximately 1 volt when the readout is 1000 digits or 100.0, 10.00, 1.000.
6. **FUEL GAS** 1/4" connector for fuel tubing.  
  
**IMPORTANT: Fuel supply must be regulated to 10-15 inches water gauge (0.36-0.54psi or 0.025-0.38kg/cm<sup>2</sup>). See section 2.1.2 for information on different fuel types.**
7. **AIR** 5/16" connector for air tubing.  
  
**IMPORTANT: Air pressure on the inlet to the instrument must not exceed 30psi (2kg/cm<sup>2</sup>). See section 2.1.3.**

## 4.2 Operation – Flame Ignition and Optimisation

1. Ensure that the drain trap is pushed fully down on its clip. Ensure that the drain trap has solution in it and that no air locks are present. If necessary, purge by adding deionised water and allowing the surplus to run away.
2. Close the **fuel** valve by turning fully clockwise. To avoid damaging the valve, it should not be forced.
3. Turn the **fuel** valve the required number of turns anti-clockwise depending on the fuel being used.

| Fuel        | Number of turns |
|-------------|-----------------|
| Propane     | 3               |
| Butane      | 4               |
| Natural gas | Fully open      |
| L.P.G.      | 3.5             |

4. Turn on the fuel supply at source i.e. cylinder.
5. Switch on electrical power by depressing the **power** switch. If the air compressor is powered separately, switch on the air compressor. Ensure that air is present by listening for the hissing created as it passes through the nebuliser.

6. Depress the ignition switch and hold down. Watch the **FLM** indicator in the display window. When this indicator is illuminated the flame is a light and the ignition switch can be released. If the **FLM** indicator does not light within approximately 20 seconds, release the switch and open the **fuel** valve one additional turn. Allow the gas to disperse before continuing. Depress the ignition switch for a further 20 seconds. This process may be repeated until successful ignition occurs. If the **fuel** valve has to be opened more than 5 turns more than recommended above, refer to Section 7.5.
7. Set the **filter select** control to the desired position.
8. Aspirate deionised water and set the readout to zero by adjusting the **blank** control.
9. Aspirate a standard solution of slightly higher concentration than expected in the samples to be tested. Adjust the **fine and coarse** control until a positive reading is obtained. This is a nominal value to be used in optimising flame conditions.
10. Adjust the **fuel** valve in a clockwise direction until a peak reading is obtained.

**NOTE:** There is a time delay between adjusting the flow of fuel and seeing the effect of the adjustment. A pause of a few seconds is therefore necessary between making every fine adjustment.


**The optimum flame conditions are different for Na and K and retuning is necessary when switching between elements.**

11. For optimum performance the instrument should be allowed 15 minutes to warm-up. During this warm-up period a blank deionised water sample should be aspirated. It should be noted that results, adequate for most purposes, can be obtained from switch on, although standardisation settings will need frequent checking during the warm-up period.

### 4.3 Calibration: PFP7

It is important to understand that the principles of flame photometry are such that, over **certain concentration ranges**, light emitted from the flame is **directly** proportional to the concentration of the species being aspirated.

The graph below shows that the direct relationship between the flame emission and concentration is only true at relatively low concentrations. Above these low levels the flame begins to saturate and the flame emission ceases to increase in a linear relationship to concentration.

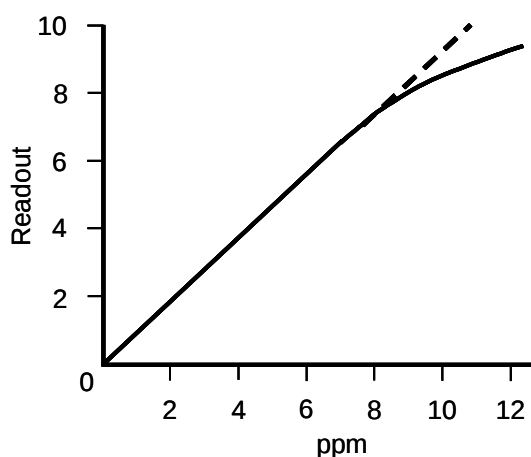**Figure 4.3.1:** Relationship between sample concentration and flame emission.


If the samples being analysed lie on the linear part of the curve then the user can take direct concentration readings from the digital display. If, however, the concentration of samples are above the levels shown on the graph then the user has to dilute the samples so that they lie on the linear part of the curve.

The user can manually construct a calibration curve using standard solutions containing known concentrations of the elements to be determined and if necessary, other materials to ensure that the standard and sample backgrounds match. The concentration range covered by the calibration curve will depend upon the expected concentration of the samples so that the sample readings fall somewhere in the middle of the calibration curve.

Once the calibration curve has been plotted, the readings for the sample solutions are compared with the curve to allow the sample concentrations to be established.

It is important to realise that each element has its own characteristic curve and separate calibration curves must be constructed.

Instrument re-calibration is easily achieved by setting the blank solution to read zero and remeasuring the top standard. The graph in **Figure 4.3.2** shows a typical curve obtained when measuring in parts per million (ppm).

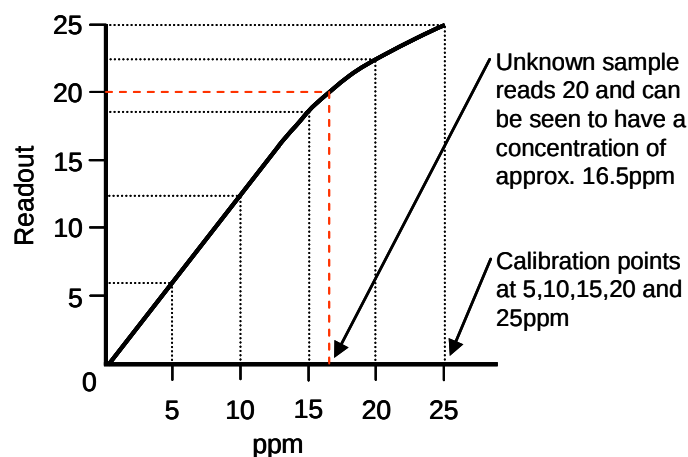**Figure 4.3.2:** Typical calibration curve measuring ppm.

Customers working in medical environments are quite likely to be using the S.I. unit of mmol/l to report their results. The relationship between mmol/l and ppm is defined below:

|           |    |                     |                 |
|-----------|----|---------------------|-----------------|
| Sodium    | Na | 1ppm = 0.0435mmol/l | 1mmol/l = 23ppm |
| Potassium | K  | 1ppm = 0.0256mmol/l | 1mmol/l = 39ppm |
| Lithium   | Li | 1ppm = 0.1441mmol/l | 1mmol/l = 7ppm  |
| Calcium   | Ca | 1ppm = 0.0250mmol/l | 1mmol/l = 40ppm |

This relationship means that Na and K samples in the normal clinical range of 136-145mmol/l Na and 3.5-5.0mmol/l K should be pre-diluted 1 in 100 or 1 in 200 to get optimum results from the flame photometer.

Aspirate a blank solution and set the readout to 000 using the **blank** control.

Aspirate the highest standard solution and set the readout to an appropriate reading using the **fine** and **coarse** sensitivity controls. Re-check the blank setting and adjust if necessary.

Aspirate the remaining standard solutions (if used) to construct the calibration curve and note the results.

When the blank and standards are set, unknown samples can be aspirated and the results noted directly from the instrument readout.

Calibration needs to be checked periodically by aspirating the blank and standard solutions. Initially this check should be carried out after every 10 samples. Experience and increased confidence in the PFP7 will enable you to best judge the frequency of this check.

The decimal point (**d.p.**) switch can be set to illuminate the decimal point in any significant position. This should be chosen to give sufficient resolution for the test required.

#### 4.4 Calibration: PFP7/C

When using the PFP7/C, a calibration curve is not required as the display is calibrated in direct concentration units. Therefore only one top standard is required to enable this calibration to be performed.

Installation and set-up procedures should be carried out as for the standard Model PFP7.

**NOTE: The Model PFP7/C allows direct readout of Na and K concentration in serum samples. Na levels in urine samples may also be read directly from the display.**

**Samples of both serum and urine should be diluted 1 in 100 or 1 in 200 to obtain optimum results from the flame photometer.**

To calibrate the unit for clinical use the following steps should be carried out:

1. Aspirate deionised water and set the display to read zero.
2. Select the appropriate element (Na or K) on the **filter select** control and the position of the decimal point using the **d.p.** switch.
3. Aspirate the required standard, e.g. 140mmol/l Na, 5.0mmol/l K which has been pre-diluted at least 1 in 100 and set the display reading accordingly using the **fine** and **coarse** sensitivity controls.
4. Adjust the **fuel** valve in a clockwise direction until a peak reading is obtained. Re-set the display reading accordingly for the standard being aspirated.

**NOTE: There is a time delay between adjusting the flow of fuel and seeing the effect of the adjustment. A pause of a few seconds is therefore necessary between making every fine adjustment.**

**The optimum flame conditions are different for Na and K and retuning is necessary when switching between elements.**

5. Re-check the zero setting and adjust if necessary.
6. Aspirate the pre-diluted sample and note the reading.
7. Calibration needs to be checked periodically by aspirating the blank and standard solutions. Initially this check should be carried out after every 10 samples. Experience and increased confidence in the PFP7/C will enable you to best judge the frequency of this check.

#### 4.5 Shutdown

1. Aspirate deionised water for at least ten minutes.
2. If the shutdown is short term i.e. the instrument is to be used again the same day/shift, it is adequate to remove electrical power from the unit by depressing the **power** switch. This will safely extinguish the flame.

3. If the shutdown is longer term or if the laboratory is likely to be left unattended during the shutdown, then the fuel supply should be turned off at source; wait for the **FLM** indicator to extinguish and then turn off the power. This ensures that there is no gas left in the tubing to the unit.

#### **4.6 Operating precautions**

1. The fuel gases used in the flame photometers are inflammable and therefore potentially hazardous. Cylinders of fuel gas should always be stored and used in line with the supplier's recommendation.
2. It is possible that a small quantity of fuel will escape from the instrument during the ignition sequence. The amount of fuel is harmless although may smell slightly. If the smell of fuel gas persists the instrument should be immediately shut down and the source of the leakage determined by using a soap solution on the hose joints.
3. Do not leave the instrument running unattended while the flame is alight.
4. The top of the instrument chimney unit becomes very hot when running and can cause severe burns if touched.
5. The exhaust gases from the flame are very hot and the area approximately 1 metre above the chimney must be avoided. Never attempt to look down the chimney whilst the flame is running. Always use the inspection window.
6. The instrument uses potentially hazardous electrical supplies. Never remove covers from the instrument without first ensuring that it has been isolated completely from the AC mains supply.
7. If the instrument is used in a pathology laboratory, all samples should be handled with the caution normally accorded to those known to contain pathogenic organisms. Care should also be taken when undertaking maintenance on instruments that have been used in these environments. A bactericidal agent should be used when cleaning parts during routine maintenance.

**For detailed safety precautions please refer to the beginning of the manual, page 1 onwards.**

## Section 5

### Accessories

#### 5.1 Water separator

To avoid unstable readings it is essential that the air supply to the mixing chamber is clean and dry. Jenway recommends that a water separator be used with all flame photometers as when air is compressed, condensation occurs. The amount of condensation will depend on the humidity in the air, so in high humidity areas, water separators are essential. The water separator fits on the air-line between the compressor and the flame photometer. This is where condensation forms. Any water collecting in the tube will cause unstable readings and if it enters the instrument, will lead to the eventual corrosion of internal components in the air-line.

##### 5.1.1 Features

1. Available in two sizes: large (normally recommended) (part code **500 177**) or small (for areas of low humidity) (part code **500 176**).
2. Complete with all parts to connect to the flame photometer.
3. Provides optimum performance in minimum space.
4. Provides effective liquid and solid separation.
5. No maintenance required.
6. Finger operated plungers for easy draining.

##### 5.1.2 Specifications

| Model                       | Large                         | Small                         |
|-----------------------------|-------------------------------|-------------------------------|
| Part code                   | 500 177                       | 500 176                       |
| Port size:                  | ¼", ⅜"                        | ⅛", ¼"                        |
| Maximum operating pressure: | 100psi                        | 100psi                        |
| Working temperature range:  | -5 to 80°C (with no freezing) | -5 to 80°C (with no freezing) |
| Nominal filtration rating:  | 5µm                           | 5µm                           |
| Bowl material:              | Polycarbonate                 | Polycarbonate                 |
| Bowl guard:                 | Yes                           | No                            |
| Drain capacity (ml):        | 25                            | 8                             |
| Weight (kg):                | 0.22                          | 0.18                          |

#### 5.2 Compressor

The compressor is a reliable, quiet air compressor of oil-less design requiring minimal maintenance. The compressor is constructed from a lightweight but rigid aluminium extrusion and is housed in a protective case to ensure maximum user safety and convenience. Load bearings are permanently lubricated and sealed for the life of the compressor. A PTFE composition cup offers excellent sealing and wear properties. The integral motor has its own cooling fan and is thermally protected against overload.

##### 5.2.1 Specification

The compressor gives a supply of dry, clean and pulse free air at a pressure between 15 and 30 psi (approximately 1-2 kg/cm<sup>2</sup>) at 6 litres/minute. There are two models available depending on the required voltage.

| Model     | 8515          | 8516       |
|-----------|---------------|------------|
| Part Code | 535 001       | 535 002    |
| Voltage:  | 230V, 50/60Hz | 110V, 60Hz |

### 5.2.2 Installation

The compressor should be positioned on a stable flat surface within 2 meters of the PFP7 or PFP7/C. Connection to the PFP7 or PFP7/C should be carried out as follows:

1. Ensure the power supply to the flame photometer is switched on.
2. Connect the air tubing (supplied with the flame photometer) to the compressor outlet nozzle and secure with a hose clip.
3. Connect the other end of the air tubing to the air inlet located.
4. Plug the compressor mains lead into either auxiliary power socket located on the rear of the panel of the flame photometer or directly into the mains power supply.

## Section 6

### Maintenance

#### 6.1 General

The design of the Model PFP7 & PFP7/C is such that maintenance requirements are minimal. To maintain good performance and prolong the life of the instrument it is important that the procedures defined in the manual are carried out regularly. The performance of the instrument depends upon an adequate supply of compressed air. Recommended compressor maintenance procedures should be carried out to ensure that compressor performance does not deteriorate.

#### 6.2 Weekly maintenance

The nebuliser should be cleaned weekly to ensure efficient nebuliser operation.

**Equipment required:** 10ml graduated measuring cylinder  
Stop watch  
Cleaning wire (500 194)

1. Switch on the air supply to the flame photometer but do not light the flame.
2. Fill the measuring cylinder with deionised water and present to the nebuliser inlet tube for one minute.
3. The consumption rate should be between 2 and 6ml/min. If correct, no further action is required.
4. If the consumption rate is too low, this is likely to be caused by a blockage in the fine capillary tubes and can usually be cleared by passing the cleaning wire through the nebuliser. If, after taking this action and re-checking the consumption rate it is still too low, then the inlet tube should be discarded and the test repeated using fresh tubing.
5. If the nebuliser operation is still unsatisfactory, remove the nebuliser from the mixing chamber, disconnect from the air line, remove the inlet tubing and soak the nebuliser in hot deionised water.
6. Dry the nebuliser before refitting and repeating the test.
7. If operation is still unsatisfactory, a new nebuliser should be fitted.

**NOTE: Under no circumstances should the nebuliser be adjusted.**

#### 6.3 Monthly maintenance

**Equipment required:** 10ml graduated measuring cylinder  
Stop watch  
Cleaning wire (500 194)  
Cleaning solution (025 171)

1. Carry out the **Weekly maintenance** schedule.

2. With the flame alight aspirate a 1 in 100 diluted sample of cleaning solution for 30 minutes, followed by deionised water for a further 30 minutes.

## 6.4 Six-monthly maintenance

Burner cleaning.

|                            |                                                                                                                                 |
|----------------------------|---------------------------------------------------------------------------------------------------------------------------------|
| <b>Equipment required:</b> | 10ml graduated measuring cylinder<br>Stop watch<br>Cleaning wire (500 194)<br>Cleaning solution (025 171)<br>Small, stiff brush |
|----------------------------|---------------------------------------------------------------------------------------------------------------------------------|

1. Carry out the **Weekly maintenance** schedule.
2. If the instrument has been running with the flame alight, allow 30 minutes for the burner system to cool down.
3. Disconnect the mixing chamber from the chimney unit by unscrewing the knurled locking ring and dropping the chamber clear of the chimney.
4. Remove the fluted burner from the top of the burner tube.
5. Rinse the burner in deionised water and remove any deposits or encrustations by brushing with a small stiff brush. If deposits are persistent, heating or boiling the deionised water will help.
6. Dry and replace the burner, pointed end downwards and re-assemble the mixing chamber onto the chimney.
7. Carry out the flushing procedure as described in **Monthly maintenance**.

## Section 7

### Trouble shooting

#### 7.1 General

This section is a step-by step guide that should allow an operator to take appropriate actions to clear simple faults. Any action not defined in this trouble shooting guide should only be undertaken by qualified personnel.

#### 7.2 Unstable results

| Possible cause                       | Solution                                                                                                                                                        |
|--------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Condensation in the air supply.      | Drain the tubing. If the problem persists a moisture-free air supply should be arranged by using a water separator.                                             |
| Nebuliser blocked.                   | Check the nebuliser performance as defined in Section 6.2, Weekly maintenance.                                                                                  |
| Flame temperature not correctly set. | Ensure that the <b>fuel</b> valve is correctly set for the element being determined; refer to Section 4.2, Operation.                                           |
| Fuel cylinder nearly empty.          | Replace cylinder with a fully charged one.                                                                                                                      |
| Instrument in strong draught.        | Relocate the instrument.                                                                                                                                        |
| Instrument in very bright sunlight.  | Relocate the instrument.                                                                                                                                        |
| Instrument not draining correctly.   | Ensure the "U" tube is clear of blockages and that the drain trap is at the correct level i.e. pushed down fully so the rim at the top is resting on the clip.  |
| Atmosphere contaminated.             | Remove source of contamination or improve ventilation.                                                                                                          |
| Mixing chamber/burner contaminated.  | Carry out the six-monthly maintenance procedures as defined in section 6.4<br><br>Ensure the instrument is always flushed with deionised water before shutdown. |

#### 7.3 Unable to set standard reading

| Possible cause                                                          | Solution                                               |
|-------------------------------------------------------------------------|--------------------------------------------------------|
| Calibration standard at incorrect concentration or incorrectly diluted. | Make new standards to check the calibration procedure. |
| Blank calibration standard contaminated or incorrectly set.             | Re-make standard and re-set blank.                     |

|                                      |                                                                                                       |
|--------------------------------------|-------------------------------------------------------------------------------------------------------|
| Nebuliser blocked.                   | Check the nebuliser performance as defined in Section 6.2, Weekly maintenance.                        |
| Flame temperature not correctly set. | Ensure the <b>fuel</b> valve is correctly set for the element being determined, refer to Section 4.2. |

#### 7.4 Non-linear results

| Possible cause                                             | Solution                                                                                              |
|------------------------------------------------------------|-------------------------------------------------------------------------------------------------------|
| Nebuliser blocked.                                         | Check the nebuliser performance as defined in Section 6.2, Weekly maintenance.                        |
| Contaminated deionised water.                              | Make up new standards in fresh water.                                                                 |
| Incorrect dilution or incorrect calibration standard used. | Prepare fresh dilutions of the calibration standard.                                                  |
| Flame temperature not correctly set.                       | Ensure the <b>fuel</b> valve is correctly set for the element being determined, refer to Section 4.2. |
| Instrument not draining adequately.                        | Ensure the “U” tube is clear of blockages and that the drain trap is at the correct level.            |

#### 7.5 Flame will not light

| Possible cause                                              | Solution                                                                                                                    |
|-------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------------------|
| Drain “U” tube empty.                                       | Fill with deionised water.                                                                                                  |
| Fuel cylinder empty or regulator pressures incorrectly set. | Replace cylinder or reset regulator.                                                                                        |
| Inadequate air supply.                                      | Check air connection to the rear panel and the operation of the compressor.                                                 |
| Propane/butane fuel restrictor blocked.                     | Carefully “blow out” the restrictor. Do not use wire to clear the blockage as this will result in damage to the restrictor. |

#### 7.6 No electrical power

| Possible cause               | Solution                                    |
|------------------------------|---------------------------------------------|
| AC supply not fully engaged. | Check the power lead is correctly inserted. |
| AC supply not available.     | Check mains supply.                         |

## 7.7 Unable to set blank

| Possible cause              | Solution                                                                                                                                   |
|-----------------------------|--------------------------------------------------------------------------------------------------------------------------------------------|
| Blank contaminated.         | Re-make the blank and re-check.                                                                                                            |
| Burner encrusted with salt. | Clean the burner as defined in Section 6.4, Six monthly maintenance.<br>Ensure that the cleaning procedure is carried out more frequently. |

## 7.8 No reading on display

| Possible cause                         | Solution                                                                       |
|----------------------------------------|--------------------------------------------------------------------------------|
| Incorrect filter selected.             | Select correct filter.                                                         |
| Calibration standard incorrectly made. | Re-make standard.                                                              |
| Nebuliser completely blocked.          | Check the nebuliser performance as defined in Section 6.2, Weekly maintenance. |

## Section 8

### Spares

#### 8.1 Minor Spares Kit (PFP7/MINKIT)

| Description                                                 | Part code | Quantity |
|-------------------------------------------------------------|-----------|----------|
| Nebuliser                                                   | I/500019  | 1        |
| Cleaning Wire (3x150mm)                                     | I/500194  | 1        |
| Cleaning Solution                                           | 025016    | 2        |
| Ignition Probe Assembly                                     | PFP7/IPA  | 1        |
| Mk2 Small Knob (Blank & Fine) Assembly                      | PFP7/SKB  | 1        |
| Mk2 Large Knob (d.p., Coarse & Filter Select) Assembly      | PFP7/LKB  | 1        |
| Mk2 Fuel Knob Assembly                                      | PFP7/FKB  | 1        |
| Fuse Kit                                                    | PFP7/FKIT | 1        |
| Drain Trap Kit                                              | I/500551  | 1        |
| Ignition Switch                                             | 017025    | 1        |
| Power Switch                                                | 017050    | 1        |
| Nebuliser to Solution Tube                                  | I/500193  | 500mm    |
| Air Regulator to Nebuliser Tube                             | 023009    | 500mm    |
| Air Regulator to Nebuliser Tube (Plastic Olive)             | 026023    | 1        |
| Air Regulator to Nebuliser Tube (Brass Nut)                 | 026007    | 1        |
| Air Regulator to Nebuliser Tube (Coupling to Air Regulator) | 026024    | 1        |
| Allen Key                                                   | 060196    | 1        |
| Foot Assembly                                               | PFP7/FA   | 1        |

# Declaration of Conformity

## Flame Photometer, Model PFP7

This product complies with the requirements of the EU Directives listed below:

**2004/108/EC**      **EMC Directive.**  
**2006/95/EC**      **Low voltage Directive (LVD)**

Compliance with the requirements of these Directives is claimed by meeting the following standards:

**EN 61326-1:2006 (Electrical Equipment for Measurement, Control and Laboratory use).**  
**EN 61010-1: 2001**  
**(Safety Requirements Electrical Equipment for Measurement, Control and Laboratory use)**  
**EN 61010-2-010:2003 (Particular requirements for laboratory equipment for heating of materials)**

**CE mark affixed 1996**

Signed:  (Mr C. Warren)

Date: 06/09

Authority: Technical Manager  
Bibby Scientific Ltd

Bibby Scientific Ltd - Stone - Staffs - ST15 0SA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

# Declaration of Conformity

## Flame Photometer, Model PFP7/C

This product complies with the requirements of the EU Directives listed below:

2004/108/EC      EMC Directive.  
2006/95/EC      Low voltage Directive (LVD)

Compliance with the requirements of these Directives is claimed by meeting the following standards:

EN 61326-1:2006 (Electrical Equipment for Measurement, Control and Laboratory use).  
EN 61010-1: 2001  
(Safety Requirements Electrical Equipment for Measurement, Control and Laboratory use)  
EN 61010-2-010:2003 (Particular requirements for laboratory equipment for heating of materials)

CE mark affixed 1996

Signed:                      (Mr C. Warren)

Date:           027 09          

Authority: Technical Manager  
Bibby Scientific Ltd

**Bibby Scientific**

Bibby Scientific Ltd - Stone - Staffs - ST15 0SA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

**Bibby Scientific Ltd**

Beacon Road Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)1785 812121  
Fax: +44 (0)1785 813748  
e-mail: [info@bibby-scientific.com](mailto:info@bibby-scientific.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**Bibby Scientific France SAS**

ZI du Rocher Vert - BP 79  
77793 Nemours Cedex  
France  
Tel: +33 1 64 45 13 13  
Fax: +33 1 64 45 13 00  
e-mail: [bsf@bibby-scientific.fr](mailto:bsf@bibby-scientific.fr)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**Bibby Scientific Ltd. – Singapore**

Prudential Tower, Level 26,  
30 Cecil Street 049712  
Singapore  
Tel: +65 6631 2976  
Fax: +44 (0) 1785 810405  
e-mail: [bibby@bibby-scientificasia.com](mailto:bibby@bibby-scientificasia.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**Bibby Scientific Middle East Ltd**

BPO Box 27842  
Engomi 2433  
Nicosia Cyprus  
Tel: +357 22 660 423  
Fax: +357 22 660 424  
e-mail: [sales@bibby-scientificme.com](mailto:sales@bibby-scientificme.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**Bibby Scientific US Ltd**

3 Terri Lane Suite 10  
Burlington NJ 08016  
USA  
Tel: 800-225-9243  
Fax: 609-589-2571  
[www.bibby-scientific.com](http://www.bibby-scientific.com)