


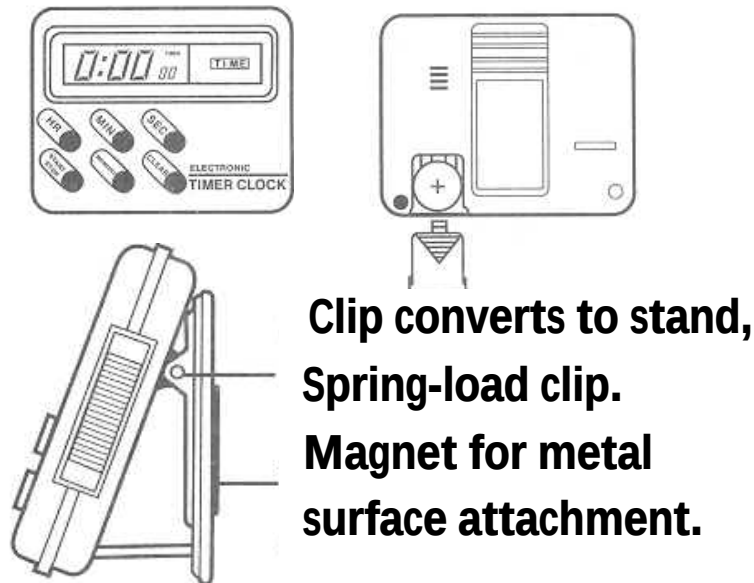

OPERATING INSTRUCTIONS (Be sure battery is properly inserted with ``+`` and ``-`` terminals) Model E 920 630

1. TO SET CLOCK

Slide switch to ``Clock``. Press and hold ``HR``, ``MIN`` or ``SEC`` button to actual time.

2. TO SET TIMER

Slide switch to ``Timer``-- press and hold ``HR``, ``MIN`` or ``SEC`` for timing. Press ``START``. To restart, press ``START``. A 60 second alarm when the timer has completed its operation.

3. TIME`S UP FEATURE

The timer will count up from 0:00:00 to tell you how long since the alarm sounded. To stop count up and/or alarm, press ``STOP`` and then clear.

4. MEMORY

Press and hold ``HR``, ``MIN`` or ``SEC`` button for desired time then press ``MEMORY`` to enter preset time (``MEMORY`` flag on display on). ``PRESET TIME`` can be recalled Instantly in ``TIMER`` mode by pressing ``MEMORY`` and operated by START/STOP button as normal to save time in setting for repeated process with same timing requested. Press ``MEMORY`` to recall preset time and then press ``CLEAR`` to cancel preset time.

5. COUNT UP

Slide switch to ``COUNT UP`` position, press ``START`` to start counting upward. Press ``STOP`` to temporarily stop timing/to restart, press ``START``. Press ``CLEAR`` to reset digits to 00HR:00MIN:00SEC.

6. 12HOUER AM7PM FORMAT OR 24 HOUR INTERNATIONAL TIME USER-SELECTABLE:

Slide switch to ``CLOCK`` position and press and hold ``CLEAR`` button for 2 sec to change 12 Hour mode to 24 Hour mode or vice versa.