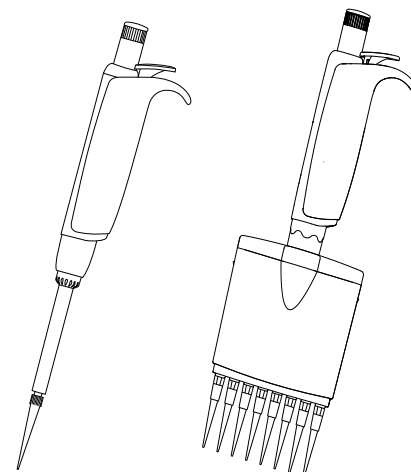
# ACURA<sup>®</sup> manual 815 / 825 / 835 / 855

## DIGITAL READING PIPETTES PIPETTEN MIT DIGITAL-DISPLAY PIPETTES À LECTURE DIGITALE

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference.
- Observe manufacturer's instructions and guidelines for operation and maintenance.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals (such as acetone) to clean instrument.
- This product should be used only for its intended purpose.
- Mind possible hand-fatigue during serial pipetting and its medical consequences (such as repetitive strain injuries RSI).

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#### DESCRIPTION (Fig. 1)

The Acura<sup>®</sup> manual is an air displacement micropipette line intended for the accurate, precise and easy pipetting. The single channel models cover volumes ranging from 0.1 µL to 10 mL. Available in eight models with 8 or 12 channels, the multichannels are perfectly adapted for a continuous microplate pipetting in the 0.5 to 350 µL range.

Materials have been carefully selected to ensure durability and optimal resistance to aggressive media. The unit is autoclavable at 121°C fully assembled.

Major features include digital display window (E) for permanent volume visibility and integrated, adjustable tip ejector Justip<sup>™</sup> (H) activated by separate push button (C). Ergonomics and pipette activation provide for an ultra-soft operation. Calibration easily performed by laboratory personal with *swift-set* system (F).

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#### OPERATION (Fig. 1)

##### Holding the pipette

Place finger rest (D) on the phalanx of forefinger. Ease of access of both plunger button (B) and ejector push button (C) allow natural and comfortable thumb movement. The multichannel casing (K) revolves to allow the selection of the best working position.

##### Volume setting

Set desired volume of variable models by rotating push button (B) clockwise to decrease the volume and vice versa, until it appears in display window (E). Set volume is exactly indicated by the top of the pointer. Precise click-stops of the micrometric mechanism and a free rotating smartie colour cap (A) on top of button prevent any volume alteration during pipetting.

**Note:** If the letter E appears beside digits in the display window, the set volume is no longer within working range of the pipette. Overturning the micrometric screw may impair the volume setting mechanism.

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### Tip selection

Disposable Qualitips® – with or without filter – were carefully designed to provide the best pipette and tip combination for all types of applications. Regular tips are autoclavable at 121°C. Filter tips are either autoclavable or supplied in a sterile version. Use tips supplied by instrument manufacturer, or similar (see tip selection on p II). Detailed flyer available on request.

Apply new tip as often as deemed necessary in regard with change of sample nature, droplet accumulation after blow out, or foam formation. When using tip racks, place the instrument on tips, press firmly to ensure a good tightness. This is particularly important for multichannel models.

**Note:** Enhance precision by filling new tip or Pasteur pipette once before sampling.

### Pasteur pipette (835 model only)

Glass Pasteur pipettes are of advantage in handling PP affecting solvents. Optional adapter-nozzles (cat. No. 1.835.631, 2 mL or 1.835.633, 5 mL), accommodate standard 2 mL Pasteur pipettes (ext. Ø 6.5-7.2 mm) in addition to Polypropylene tips. Refer to instructions for disassembling volumetric module for changing the adapter nozzle. Performance shown on QC certificate are obtained with PP tips, recalibrate if necessary with Pasteur pipette.

**Note:** Do not set volume higher than 2 mL.

### Tip ejection (Fig. 2 and 3)

The patented Justip™ ejection system allows instant height adjustment ( $\pm 2$  mm) of ejector for best fitting of tip used (optimal distance). Set correct position ( $\sim 0.5$  mm space between tip and ejector) by rotating ejector screw (< LO – HI >). Click stops prevent any unwanted change while pipetting.

A special design ensure sequential ejection on multichannel models.

### Protection filter (835 models)

Macropipettes (2, 5 and 10 mL models) accommodate a nozzle protection filter, which drastically reduces risks of barrel and plunger contamination. Its use is particularly recommended with large volumes and/or when handling biohazardous, radioactive or highly aggressive liquids. Change filter if wet or contaminated. Filter must not be autoclaved.

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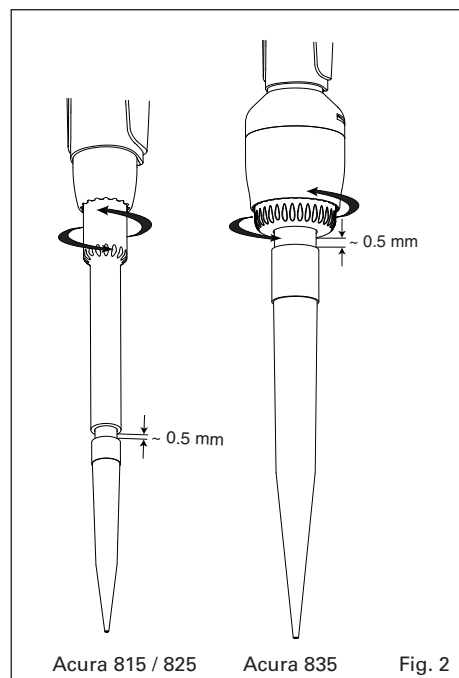
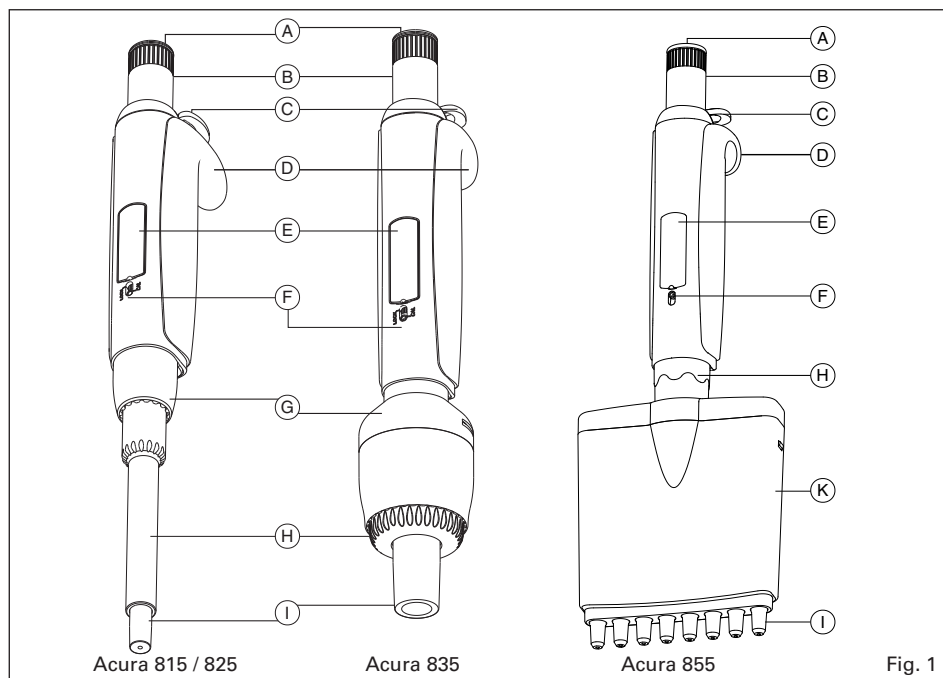
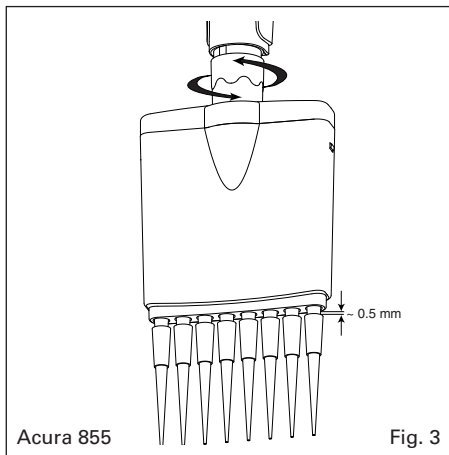
## PIPETTING OPERATION

Before pipetting, tightly fit new, clean tips on the nozzle.

### Forward pipetting (Fig. 4)

Exact volume desired is aspirated, then dispensed.

- Phase 1** Depress plunger button to intermediate stop ending calibrated stroke.
- Phase 2** Immerse tips vertically (2 - 3 mm) into liquid, wait 2 sec. and smoothly release plunger button. Withdraw tips from liquid, without touching container wall.
- Phase 3** Place tips against wall of receiving vessel, depress plunger button to first stop and let liquid out.
- Phase 4** Depress button to second stop: blow out will expel remaining droplets. Withdraw tips from receiving vessel, sliding it over 10-15 mm along container wall.
- Tip ejection** Remove used tips by pressing ejector button (C). Change tips as required.

### Reverse pipetting (Fig. 5)

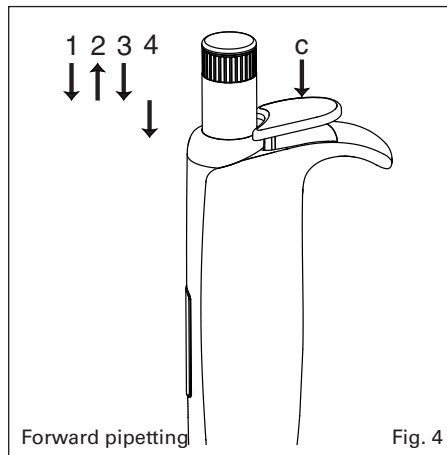
Liquid is aspirated in excess of the selected volume but only the set volume is dispensed.

- Phase 1** Depress plunger button in full.
- Phase 2** Immerse tips vertically (2 - 3 mm) into liquid, wait 2 sec. and smoothly release plunger button. Withdraw tips from liquid, without touching container wall.
- Phase 3** Place tips against wall of receiving vessel, depress plunger button to first stop only and let liquid out.
- Phase 4** Withdraw tips from receiving vessel, sliding it over 10 -15 mm along container wall and start operation to phase 2 again, or discard residual liquid by depressing button to second stop.
- Tip ejection** Remove used tips by pressing ejector button. Change tips as required.

### Measuring unknown volumes (825/835 models only)

The fine volume resolution of the micrometric mechanism allows for precise liquid measurement by aspiration. Remember set volume before starting, dip pipette tip into liquid and turn adjustment button counter-clockwise until unknown quantity is completely aspirated. Measured volume consists in read volume minus initial set volume.

**Example:** Volume before aspiration: 10.0 µL. Volume read-out after aspiration: 64.7 µL. Liquid aspirated: 64.7-10.0 = 54.7 µL.

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## MAINTENANCE

The Acura® manual pipettes are in conformity with Council directive IVD 98/79 EEC (in vitro diagnostic medical devices). Consequently, any change on the pipettes (i.e. parts, calibration) is done under responsibility of the user. The models were designed to sustain long-term operation with minimal maintenance and service requirements. However it is recommended to control instrument performance in accordance with laboratory internal procedures (SOP, GLP, etc) or at least once a year. Defect parts must be replaced, or instrument returned to dealer for repair. Order original spare parts from authorised dealers only.

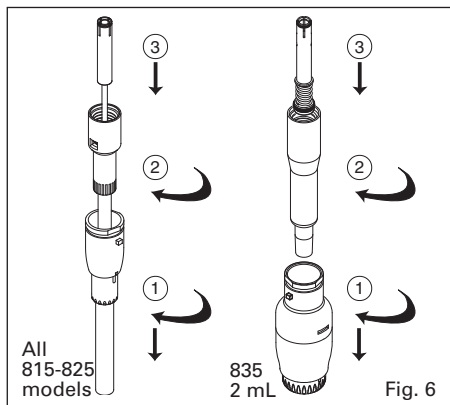
### Tightness control

A tight air chamber is paramount for the instrument to provide accurate and precise results. Leaks can be detected by observing a droplet forming and falling from the extremity of the tips or unequal liquid level in the tips of a multichannel model. Special attention must be paid to the following points in case of leakage:

- Matching of tips and nozzle?
- Damaged nozzle?
- Worn-out PTFE sleeve and/or O-ring?

### Cleaning

Elements of volumetric module (lower assembly) can be cleaned with alcohol, or soaked in appropriate decontamination or disinfecting solutions. This operation is particularly important if liquid accidentally entered the pipette. Rinse and dry instrument before reassembly. If ultra-sonic bath is used to remove sticking residues, slightly grease O-ring and / or plunger before reassembly.

### Disassembling volumetric module and O-ring change, 815/825/835 - 2 mL models (Fig. 6)

- Removing volumetric module (lower assembly):
1. Press ejector button (2), turn part (20) or (24) for 2 mL to the left and pull.
  2. Unscrew barrel (18).
  3. Depress plunger button in full then unclip (pull) plunger.

#### O-ring and PTFE sleeve replacement (815/825):

- After removing volumetric module, take barrel (18), press both clips of ring (14) with a pipette tip or a pointed tool.
- Remove ring from barrel, then remove spring (15) to access O-ring/PTFE sleeve assembly.
- Change parts if needed. Clean plunger (13) to eliminate any grease residue.

**Note:** To prevent damages when reassembling, first place PTFE sleeve (16) on plunger, then O-ring (17) on sleeve.

- Slightly grease plunger, sleeve and O-ring and reassemble volumetric module.

**Note:** Tightness parts are not available separately on 2, 10, 10Y and 20 µL models. In case of tightness problem, the whole barrel must be changed.

#### O-ring replacement (835 - 2 mL):

- Unscrew plunger rod (21) and remove washers (22, 23) and spring (15).
- Change parts if needed. Uniformly grease O-ring (17) and barrel (18).

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### Disassembling volumetric module and O-ring change, 835 5 and 10 mL models (Fig. 7)

Removing volumetric module (lower assembly):

1. Slightly turn ejector nut (20) to unclip from ejector cap (24).
2. Unscrew barrel (18), depress push button in full then pull barrel to unclip.

O-ring replacement:

3. Slightly and simultaneously press both barrel clips (18) with fingers to separate from bonnet (25).
- Pull out plunger assembly.
- Unscrew plunger rod (21) and remove washers (22, 23) and spring (15).
- Change parts if needed. Uniformly grease O-ring (17) and barrel (18).
- Reassemble plunger assembly, barrel and bonnet, screw barrel on pipette, then clip plunger rod by pressing in full plunger button (B).

To use both Pasteur pipettes and PP tips, replace regular barrel by Pasteur pipette adapter nozzle (1.835.633).

### Disassembling volumetric module and barrel change, 855 model (Fig. 8)

Removing volumetric module (lower assembly):

1. Turn ejector nut clockwise till its lowest position.
2. Press ejector button (2), turn ejector nut to the left, unclip ejector ring and release ejector button.
3. Hold volumetric module firmly with one hand and, while pulling down casing, rotate slowly to unscrew.
4. Press plunger button then unclip plunger rod (28).

Barrel change (Fig 9 to 11):

**Note:** Tightness O-ring cannot be removed from barrel. Change barrel if tightness is deficient.

1. Press with pointed tool on both clips of cover (27) and remove casing (31).
2. Press bottom clips a) of barrel holder and separate.
3. Pull out barrel assembly without bending the plunger head.

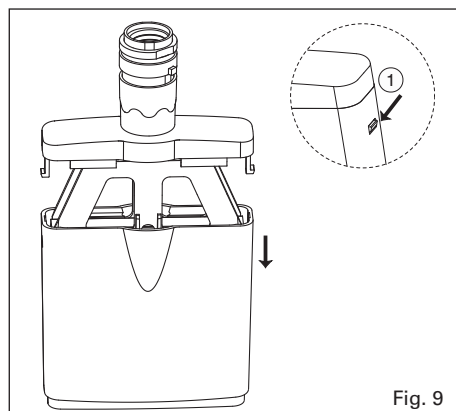
Fig. 9

### Reassembling volumetric module, 855 model:

- Apply thin, regular grease layer on all surface length of plunger.
- Separate both plate of barrel holder.
- Introduce barrel (30) on plunger (29) without bending it.
- Clip barrel plate, all barrels must be sitting properly and aligned.
- Introduce casing (31) and clip on cover (27).

**Note:** Markings on casing (volume) and on cover (Justip) should not appear simultaneously.

- Hold casing with fingers pulling against barrel extremities.
- Slowly screw volumetric module.
- Press ejector button (2) and introduce tooth of ejector ring in recess.
- Turn counter clockwise to click in ejector rod (11) and release ejector button.
- Press firmly plunger button until plunger rod (28) clips in.
- Control proper functioning of plunger and ejector buttons.
- Set ejector height to desired position.

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## TROUBLE SHOOTING

Observation	Possible cause	Action
<b>Tips not fitting the nozzle</b>	Inadequate position of tip ejector Inadequate tips	Adjust position of tip ejector Use original or other compatible tips
<b>Pasteur pipette does not hold or too «tight»</b>	Pasteur not compatible	Check/change O-rings Grease slightly O-rings
<b>Plunger sticking or dragging</b>	Dirty plunger	Disassemble and clean plunger (835, 855 grease as well)
<b>Liquid not aspirated</b>	Dry residues block pipette nozzle Volumetric module not properly assembled Dirty or wet nozzle filter on 835 models	Disassemble and clean Reassemble according to instructions Change nozzle filter
<b>Uneven liquid level in tips of multichannel model</b>	Lack of tightness	Check tip fitting Change for damaged nozzle or tightness parts
<b>Poor instrument performance</b>	Lack of tightness Inadequate tips  Dirty or wet nozzle filter on 835 models Instrument out of calibration Instrument used with viscous or volatile liquids or at particular temperatures	Check for damaged nozzle or tightness parts Check tip compatibility and fitting Pasteur not compatible, check tightness with adapter Change nozzle filter Perform new calibration Perform new calibration with liquid or at temperature used

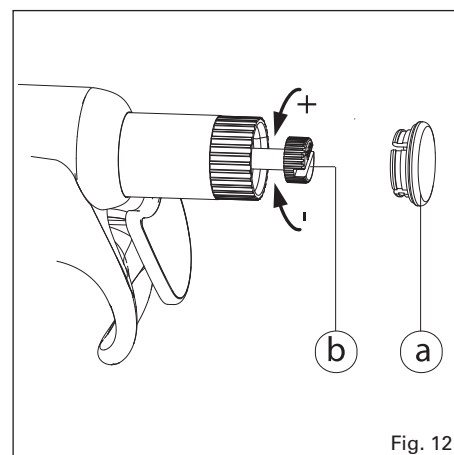
Fig. 12

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## Sterilisation

Instrument has been designed for repeated sterilisation in the autoclave at 121°C (20 minutes), fully assembled. Before autoclaving, remove nozzle filter on 835 models. Place instrument horizontally in the autoclave, avoiding any direct contact with metal.

Allow instrument to dry and cool down before use. Insert new nozzle filter on 835 models. Check tightness and accuracy regularly but at least after 50 autoclaving cycles. Check volumetric module, tighten barrel if loose. Repeated autoclaving may affect material colours and free rotation of smartie colour caps. Correct autoclaving and resulting sterility are the responsibility of the user.

## CALIBRATION

Each Acura® *manual* has been factory calibrated and individually controlled. The QC certificate accompanying the instrument includes control data, serial number and operator identification. Calibration is carried out gravimetrically with aqua-bidest. at constant temperature ( $\pm 0.5^\circ\text{C}$ ) comprised between 20 and 25 °C according to ISO 8655.

### In-lab calibration

The swift-set calibration system with integrated key and locking mechanism makes tedious procedures obsolete. If performance results are no longer within recommended values, for instance after QC check or replacing parts, or if changes occur in physical parameters (liquid density, temperature), recalibration is easily performed using an analytical balance. Ten measurements should be first performed on small volume after checking seal tightness and perfect working condition of the instrument.

### Adjustable volume models (Fig. 12 and 13)

- Remove smartie cap (a) on plunger button and pull out calibration screw (b).
- Remove "Calibration seal" (c) and move calibration cursor (d) on cal position (use pointed item such as tip or clip).
- Turn calibration screw until average weighed volume appears in the display window, without holding push button.
- Place calibration cursor on *lock* position. Cover plunger button with Smartie cap again.
- Perform few plunger movements and check resulting volume.
- Place a new "Calibration seal" over the calibration cursor.

### Fixed volume models (Fig. 12)

Same as above, except

- Turn calibration screw (b) clockwise to decrease and counter-clockwise to increase resulting volume.
- Check resulting volume after each fine tuning, until pipette performance are within acceptable values.

## WARRANTY

Your Acura® *manual* pipette is guaranteed against any material or manufacturing defects for the period of time specified in its QC certificate. Damages due to non-respect of manufacturer's instructions, safety precautions or autoclaving conditions, as well as material colour alteration are excluded from the warranty. Repair and replacement of parts do not extend warranty time.

Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased after obtaining return authorization.

**Decontaminate instrument prior to returning it.**

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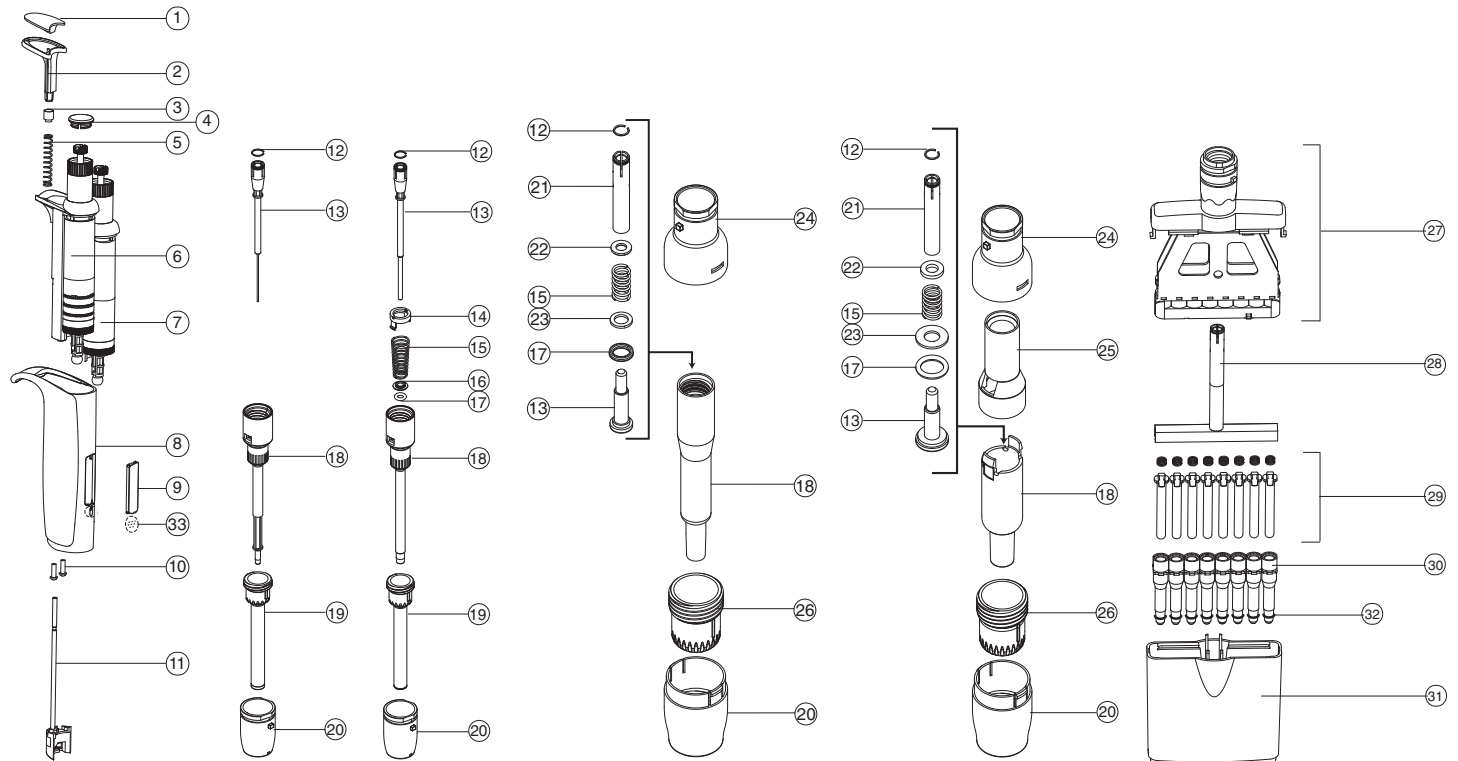
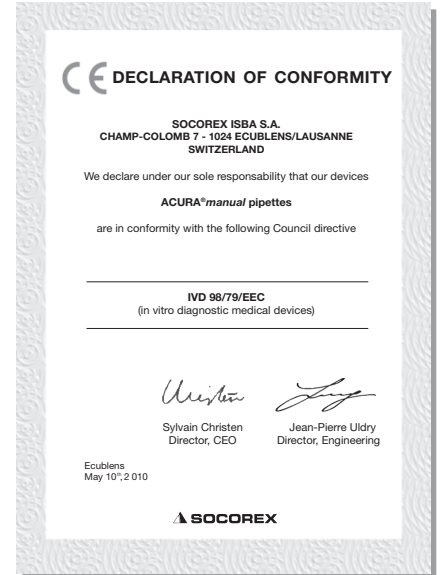
SAS Accredited  
Service Laboratory  
ISO/IEC 17025

**G.M. SYSTEM**  
ISO 9001/13485  
CERTIFIED

**PERFORMANCE AND ORDERING INFORMATION**

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<sup>1)</sup> Measured at 0.5  $\mu$ L. Indicative data at 0.2  $\mu$ L: E<= 12%, CV<8%. <sup>2)</sup> Measured at 1  $\mu$ L. Indicative data at 0.5  $\mu$ L: E <+5%, CV<3%.

**Acura 835, adjustable volume macropipettes**

Volume mL	Division mL	Inaccuracy (E%)			Imprecision (CV%)			Tip style	Cat. No.
		Min. vol.	Mid. vol.	Max. vol.	Min. vol.	Mid. vol.	Max. vol.		
0.2 – 2	0.002	< $\pm$ 1.5 %	< $\pm$ 1.0 %	< $\pm$ 0.5 %	< 0.5 %	< 0.3 %	< 0.2 %	2 mL	835.02
0.2 – 2 <sup>1)</sup>	0.002	< $\pm$ 1.5 %	< $\pm$ 1.0 %	< $\pm$ 0.5 %	< 0.5 %	< 0.3 %	< 0.2 %	2 mL/Pasteur	835.02PP
0.5 – 5	0.01	< $\pm$ 1.5 %	< $\pm$ 1.1 %	< $\pm$ 0.6 %	< 0.6 %	< 0.5 %	< 0.3 %	5 mL	835.05
0.5 – 5 <sup>1)</sup>	0.01	< $\pm$ 1.5 %	< $\pm$ 1.1 %	< $\pm$ 0.6 %	< 0.6 %	< 0.5 %	< 0.3 %	5 mL/Pasteur	835.05PP
1 – 10	0.01	< $\pm$ 1.5 %	< $\pm$ 0.7 %	< $\pm$ 0.5 %	< 0.5 %	< 0.3 %	< 0.2 %	10 mL	835.10

Measurements done with nozzle protection filter. Above values obtained using PP tips.

<sup>1)</sup> Model includes Pasteur adapter nozzle.

**Acura 855, adjustable volume multichannel pipettes**

Volume $\mu$ L	Division $\mu$ L	Inaccuracy (E%)			Imprecision (CV%)			Tip style	Cat. No.
		Min. vol.	Mid. vol.	Max. vol.	Min. vol.	Mid. vol.	Max. vol.		
<b>8-channel</b>									
0.5 – 10	0.01	< $\pm$ 3.5 % <sup>1)</sup>	< $\pm$ 2.5 %	< $\pm$ 1.5 %	< 3.0 % <sup>1)</sup>	< 2.0 %	< 1.0 %	Ultra 10 $\mu$ L	855.08.010
5 – 50	0.1	< $\pm$ 1.0 %	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< 1.0 %	< 0.7 %	< 0.4 %	200 $\mu$ L	855.08.050
20-200	0.2	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< $\pm$ 0.7 %	< 0.6 %	< 0.5 %	< 0.3 %	200 $\mu$ L	855.08.200
40 – 350	0.4	< $\pm$ 1.0 %	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< 0.6 %	< 0.5 %	< 0.3 %	350 $\mu$ L	855.08.350
<b>12-channel</b>									
0.5 – 10	0.01	< $\pm$ 3.5 % <sup>1)</sup>	< $\pm$ 2.5 %	< $\pm$ 1.5 %	< 3.0 % <sup>1)</sup>	< 2.0 %	< 1.0 %	Ultra 10 $\mu$ L	855.12.010
5 – 50	0.1	< $\pm$ 1.0 %	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< 1.0 %	< 0.7 %	< 0.4 %	200 $\mu$ L	855.12.050
20-200	0.2	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< $\pm$ 0.7 %	< 0.6 %	< 0.5 %	< 0.3 %	200 $\mu$ L	855.12.200
40 – 350	0.4	< $\pm$ 1.0 %	< $\pm$ 0.9 %	< $\pm$ 0.8 %	< 0.6 %	< 0.5 %	< 0.3 %	350 $\mu$ L	855.12.350

<sup>1)</sup> Measured at 1  $\mu$ L. Performance values obtained with bi-dist. water at constant temperature ( $\pm$  0.5°C) comprised between 20 and 25°C in accordance with ISO 8655.

**Acura 835, fixed volume macropipettes**

Volume $\mu$ L	Inaccuracy (E%)	Imprecision (CV%)	Socorex Cat. No.	Tip style
2	<+/- 0.8 %	< 0.5 %	835.F02	2 mL <sup>1)</sup>
2.5	<+/- 0.8 %	< 0.5 %	835.F02.5	5 mL
5	<+/- 0.7 %	< 0.3 %	835.F05	5 mL
10	<+/- 0.5 %	< 0.2 %	835.F10	10 mL

<sup>1)</sup> 5 mL tip up to serial number 17091050

**Smarties plunger caps**

Colour	Socorex Cat. No.	Colour	Socorex Cat. No.
Smartie plunger caps, alternative colour selection (6 / pack)			
White	825.710	Rose	825.719
Grey	825.715	Pink	825.720
Vanilla	825.716	Purple	825.721
Lemon	825.712	Ice Blue	825.722
Yellow	825.717	Blue	825.713
Orange	825.718	Mint	825.723
Red	825.714	Green	825.711
Smartie mix assorted colours (14 / pack)			1.825.700

**ACCESSORIES**

**ZUBEHÖR**

**ACCESSOIRES**

Description	Packaging	7 positions	6 positions
		Model 337	Model 336
<b>Universal work station – fits single channel pipettes of most brands</b>			
Grey	1 / pack	320.337G	320.336Q
Blue	1 / pack	320.337B	320.336B
Yellow	1 / pack	320.337Y	320.336Y
Green	1 / pack	320.337M	320.336G
Pink / Red	1 / pack	320.337R	320.336R
Orange	1 / pack	–	320.336O
White	1 / pack	–	320.336W

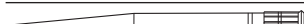
Description	Packaging	Socorex Cat. No.
Work station 340 for 3 multi- or single-channel pipettes	1 / pack	320.340
Shelf pipette holder for 2 instruments (transparent blue)	1 / pack	320.332
	4 / pack	320.332.4

V-shape, 70 ml, autoclavable PP reservoir	20 / pack	330.01
Large capacity with lid, 125 mL, autoclavable PP reservoir	20 / pack	330.01L
8-channel, 48 mL, sterile, single wrapped PS reservoir	10 / pack	330.08.9
12-channel, 48 ml, sterile, single wrapped PS reservoir	10 / pack	330.12.9

Nozzle protection filters, PP fibres, for 2 and 5mL models	250 / pack	322.05
Nozzle protection filters, PP fibres, for 10mL model	100 / pack	322.10
Pasteur pipette adapter, 2 mL model	1 / pack	1.835.631
Pasteur pipette adapter, 5 mL model	1 / pack	1.835.633
Protective calibration seal stickers	25 / bag	1.835.025
	100 / bag	1.835.100

**QUALITIPS® PIPETTE TIPS**
**QUALITIPS® PIPETTENSPITZEN**
**EMBOUTS QUALITIPS®**

Superior pipette tips cover a range between 0.2 µl and 10 ml. All models autoclavable at 121 °C.

Tips without filter	Description	Vol. (µL)	Racks	Socorex Cat. No.	Bags	Socorex Cat. No.
	Natural ultra-microtip	10	10 x 96/rack	309.0010R	1 x 1000/bag	309.0010B
	Natural gel load tip	10	4 x 204/rack	302.0010GR		
	Natural microtip	10	10 x 96/rack	302.0020R	1 x 1000/bag	302.0020B
	Natural universal tip	200			2 x 500/bag	307.0200B
	Yellow universal tip	200	10 x 96/rack	327.0200R	2 x 500/bag	327.0200B
	Natural tip	200	10 x 96/rack	308.0200R	1 x 1000/bag	308.0200B
	Natural gel load tip	200	4 x 204/rack	308.0200GR		
	Natural tip	200	10 x 96/rack	309.0200R	2 x 500/bag	309.0200B
	Yellow tip	200	10 x 96/rack	329.0200R	2 x 500/bag	329.0200B
	Natural tip	350	10 x 96/rack	308.0350R	1 x 1000/bag	308.0350B
	Natural universal tip	1000	10 x 60/rack	307.1000R	4 x 250/bag	307.1000B
	Blue tip	1000	8 x 60/rack	318.1000R	4 x 250/bag	318.1000B
	Natural tip	1000	8 x 60/rack	309.1000R	4 x 250/bag	309.1000B
	Blue tip	1000	8 x 60/rack	319.1000R	4 x 250/bag	319.1000B
	Natural macrotip	2 mL			2 x 250/bag	312.02
	Natural macrotip	5 mL	2 x 50/rack	312.05R	2 x 250/bag	312.05B
	Natural macrotip	10 mL	4 x Rack alone	312.10ER	3 x 100/bag	312.10

Efficient protection against aerosols. Include high density PE filter for complete, uncontaminated sample retrieval. Certified DNase and RNase free. Sterile tips are guaranteed pyrogenfree

Filters tips	Description	Vol. (µL)	Racks	Socorex Cat. No.	Bags	Socorex Cat. No.
	Sterile ultra-microtip	10	10 x 96/rack	309.0010FR		
	Sterile microtip	10	4 x 100/rack	302.0010FR		
	Autoclavable microtip	10			1 x 1000/bag	302.0010AFB
	Sterile tip	20	10 x 96/rack	308.0020FR	5 x 100/bag	308.0020FS
	Sterile gel load tip	20	4 x 204/rack	308.0020GFR		
	Sterile tip	100	10 x 96/rack	308.0100FR		
	Sterile gel load tip	100	4 x 204/rack	308.0100GFR		
	Sterile extended tip	200	8 x 204/rack	308.0200LFR	5 x 100/bag	308.0200LFS
	Sterile tip	300	10 x 96/rack	308.0300FR		
	Autoclavable tip	300			1 x 1000/bag	308.0300AFB
	Sterile tip	1000	10 x 100/rack	309.1000FR	5 x 100/bag	309.1000FS
	Autoclavable tip	1000			1 x 1000/bag	309.1000AFB

B = bag, R = rack, F = filter, G = gel, L = long, S = sterile, A = autoclavable

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SOCOREX ISBA S.A. - CH. CHAMP-COLOMB 7 - CH-1024 ECUBLENS/LAUSANNE - SWITZERLAND