# Thermostatschränke TC-Serie

## TC 140 G / TC 256 G

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Seite 2 - 9

**GB Instruction Manual**

Page 10 - 17

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Page 18 - 25

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Página 26 - 33

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Página 34 - 41

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43 82 35

## Disposal notes

The appliance contains reusable materials and should be disposed of properly - not simply with unsorted household refuse. Appliances which are no longer needed must be disposed of in a professional and appropriate way, in accordance with the current local regulations and laws.

When disposing of the appliance, ensure that the refrigerant circuit is not damaged to prevent uncontrolled escape of the refrigerant it contains (data on type plate) and oil.

### Do not vent R134a into atmosphere.

**R134a is a fluorinated greenhouse gas, covered by Kyoto Protocol, with a Global Warming Potential (GWP) = 1300.**

- Disable the appliance.
- Pull out the mains plug.
- Cut through the connection cable.
- Disable safety lock.

### ⚠ WARNING

Danger of suffocation due to packing material and plastic film! Do not allow children to play with packaging material. Take the packaging material to an official collection point.

## Range of appliance use

The thermostat-controlled incubators are designed to allow continuous regulation of temperature for various fields of application, e.g.

- 20 °C BOD<sub>5</sub> measurement and research purposes at:
- 4 °C storage of sewage samples
- 25 °C enzyme activity (TTC-test)
- 37 °C colony count
- Microbiology
- Food industry
- Laboratories

All temperature control problems in the usual range of 2 °C to 40 °C are solved by the factory set universal plug-in controller which is maintenance-free and reliable. This appliance is not suitable for storing flammable liquids, acids and lyes.

## Function

The inner temperature of the fully insulated incubator is controlled exactly by an integrated corrosion-proof temperature probe which connects or switches-off either the cooling unit or the heating element.

The measured inner temperature is directly indicated by a display. The temperature range of 2 °C to 40 °C can be set in increments of 0,1 °C with the aid of two tactile response keys. The keyboard is protected by a rugged foil. Aircirculation is provided by 2 axial flow fans with a discharge width of 160 cm<sup>2</sup> and an output of 300 m<sup>3</sup>/h.

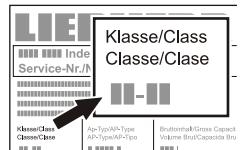
## Noise emissions from the appliance

The noise level while the appliance is operating is below 70 dB(A) (relative noise level 1 pW).

## Climate rating

The climate rating indicates the room temperature at which the appliance may be operated in order to achieve full refrigeration performance.

The climate rating is indicated on the type plate.

### Climate rating      Room temperature

4 (SN)	+10°C to +32°C
4 (N)	+16°C to +32°C
4+ (ST)	+16°C to +38°C
4+ (SN-ST)	+10°C to +38°C
5 (T)	+16°C to +43°C
5 (SN-T)	+10°C to +43°C

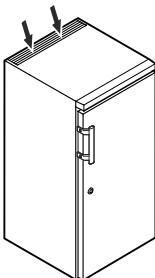
**Do not operate the appliance outside the specified room temperature range.**

## Safety instructions and warnings

- To prevent injury or damage to the unit, the appliance should be unpacked and set up by two people.
- In the event that the appliance is damaged on delivery, contact the supplier immediately before connecting to the mains.
- To guarantee safe operation, ensure that the appliance is set up and connected as described in these operating instructions.
- Disconnect the appliance from the mains if any fault occurs. Pull out the plug, switch off or remove the fuse.
- When disconnecting the appliance, pull on the plug, not on the cable.
- Any repairs and work on the appliance should only be carried out by the customer service department, as unauthorised work could prove highly dangerous for the user. The same applies to changing the mains power cable.
- Do not allow naked flames or ignition sources to enter the appliance. When transporting and cleaning the appliance, ensure that the refrigerant circuit is not damaged. In the event of damage, make sure that there are no ignition sources nearby and keep the room well ventilated.
- Do not stand on the plinth, drawers or doors or use them to support anything else.
- This appliance can be used by children of 8 years old and over, and also by persons with restricted physical, sensory or mental capacity or lack of experience and knowledge, if they are supervised or have been instructed on safe use of the appliance and understand the resulting risks. Children must not be allowed to play with the appliance. Cleaning and user maintenance must not be carried out by children without supervision.
- Avoid prolonged skin contact with cold surfaces or chilled/frozen food. This could cause pain, numbness and frostbite. In the case of prolonged skin contact, protective measures should be taken, e.g. gloves should be worn.
- Do not store explosives or sprays using combustible propellants such as butane, propane, pentane, etc. in the appliance. Electrical components might cause leaking gas to ignite. You may identify such sprays by the printed contents or a flame symbol.
- If you have a lockable appliance, do not keep the key near the appliance or within reach of children.
- The appliance is designed for use in enclosed areas. Do not operate the appliance outdoors or in areas where it is exposed to splash water or damp conditions.
- Do not place items with too high temperature in the interior.

## Setting up


- Do not place the appliance in direct sunlight or near cookers, radiators and similar sources of heat.
- The floor on which the appliance stands should be horizontal and level. Compensate for uneven floors with the adjustable feet.
- Do not cover ventilation openings or grille. 
- Always install the appliance directly against the wall.
- Well ventilated rooms

### Important

When unpacking and setting up use protective gloves.
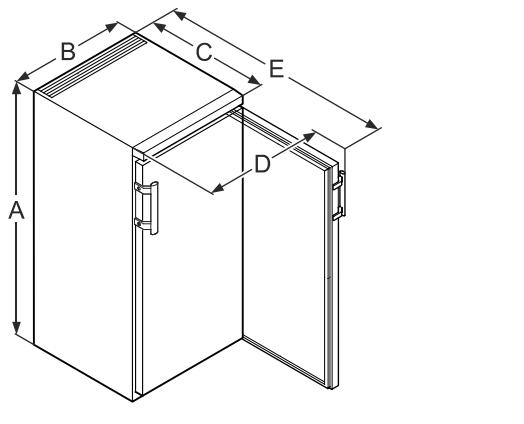
## Appliance dimensions


<b>TC 140 G</b>	850 mm	600 mm	600 mm	648 mm	1153 mm
<b>TC 256 G</b>	1640 mm	600 mm	610 mm	650 mm	1180 mm

## Installing beneath a worktop TC 140 G

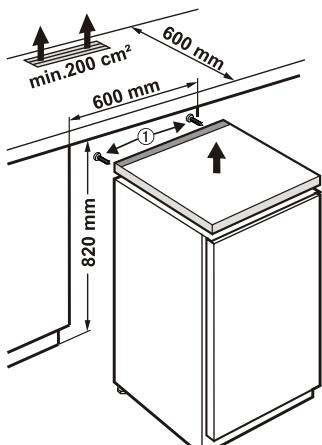For continuous worktops the refrigerator's tabletop can be detached. The refrigerator can then be pushed under the existing worktop.

- To take off tabletop remove screws ①.
- A cross-section of min. 200 cm<sup>2</sup> is required in the worktop for ventilation of the rear of the appliance.


## Electrical connection

Only operate the appliance with **alternating current (AC)**. The permissible voltage and frequency are indicated on the type plate. The socket must be properly earthed and protected by a fuse. The tripping current of the fuse must be between 10 A and 16 A. The socket must not be situated behind the appliance and must be easily accessible.

Do not connect the appliance using an extension cable or extension socket.

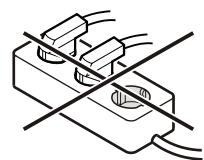Do not use stand-alone inverters (conversion of direct current to alternating current/three-phase current) or energy-saving plugs. Risk of damage to the electronic control system!


## Safety lock TC 140 G

The lock in the appliance door is equipped with a safety mechanism.

### Locking the appliance

- Insert the key as shown by arrow 1.
- Turn the key 90°.

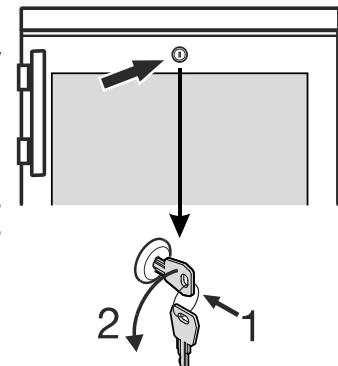To unlock the appliance, the same procedure must be repeated in the same order.


## Safety lock TC 256 G

The lock in the appliance door is equipped with a safety mechanism.

### Locking the appliance

- Insert the key as shown by arrow 1.
- Turn the key 90°.

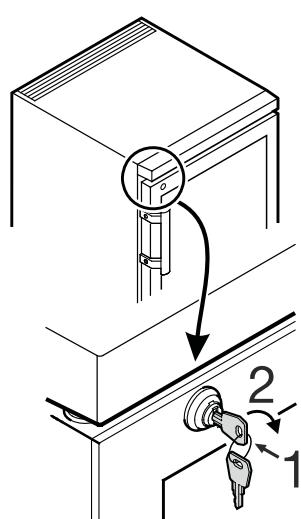To unlock the appliance, the same procedure must be repeated in the same order.

### Important

#### TC 140 G, TC 256 G

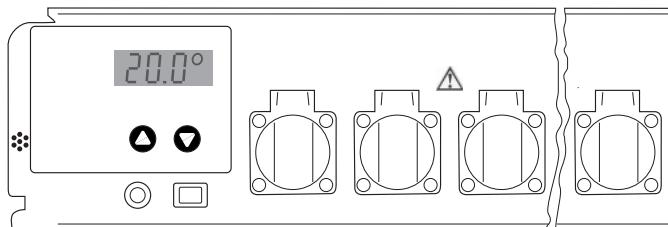
The maximum load per grid shelf is 45 kg.

## Adjustment and operation

The unit is switched on and off via the On/Off switch.

The „Up“ key is used to increase the temperature

The „Down“ key is used to decrease the temperature

The illustration shows the actual temperature of 20.0 °C.

## Setting the temperature values

When the unit is switched on, the display will initially show briefly the type of oven and a version number. It will then display the preset temperature, together with the symbol for the specified figure (3 horizontal lines). After 5 seconds, this will be replaced by the current, actual temperature.

The unit is preset in the factory to a working temperature of 20 °C. If this is required to be 25 °C, for example:

Press the „ $\uparrow$ “ key once. The display will show the specified figure which has been set, together with the symbol for the specified figure (3 horizontal lines). Now press the „ $\uparrow$ “ key again, or hold it down until the display shows the new figure required (25 °C in this example). Then release the key. The display will show the new setting for 5 seconds and will then display the actual, current temperature.

If a lower working temperature is required, the same procedure applies, using the „ $\downarrow$ “ key. The temperature can be adjusted in steps of 0.1 °C between 2 °C and 40 °C.

If you wish to read the specified temperature but without changing it, simply press the „ $\uparrow$ “ or „ $\downarrow$ “ key once and release it, so that the display shows the specified figure instead of the current temperature. The specified figure cannot be changed unless one of the keys is pressed while the display is already showing the specified figure. The preset temperature will not be lost if there is a power failure or if the unit is switched off. The data are stored in an internal electronic memory and are called up automatically when power is restored or the unit is switched on.

## Detecting a power failure

When the unit is first switched on, the display will flash. Pressing the „ $\uparrow$ “ or „ $\downarrow$ “ key will cause it to stop flashing and remain permanently lit.

If there is a power failure, the display will start to flash again. This may indicate that the temperature in the oven can no longer be held at the working temperature, if the power is off for an extended period. The controls will continue to operate normally, even if the display is flashing. The flashing merely indicates that the unit has been switched off and then back on again, or that the power supply failed for a time and no key has been pressed since then.

It is important to note that the display will flash, even if the power supply has failed for a very short period. It does not necessarily mean that the temperature in the oven has deviated significantly from the specified level.

## Temperature control and accuracy

The thermostatically controlled incubators are not specified according ASTM E 145 (9-point-measurement) or DIN 12880, part 2 (27-point-measurement).

Specifications are based on a one-point-measurement.

### Test procedure

1. Place a sample bottle filled with 500 ml water with a stirring rod on a stirring place and place it in the middle of a grid (5th insert from the bottom).
2. Completely immerse the probe of the meter in the water sample (according to the specifications of the meter). The probe should not touch the bottle wall.  
Specifications of the meter:  
Accuracy at least 0.5°C  
Resolution at least 0.1°C
3. The temperature is read after 4 hours.

## Trouble shooting

If the verification of the temperature accuracy is not successful the following should be taken under consideration:

- The ambient temperature of the thermostatically controlled incubator must match the climate rate (type plate).
- A certain circulation of air is necessary for the compressor to work properly.
- The incubator should not be opened to often and / or for longer period of time.
- Check the function of the two fans by visual inspection.
- Check the cooling device by setting 10 °C and wait for a significant change of the temperature in the sample towards 10 °C.
- Check the heating device by setting 40 °C and wait for a significant change of the temperature in the sample towards 40 °C.
- Depending on the sample temperature it can take up to 7 hours for temperature stability at the set-value.

### Important

The air temperature in the cabinet may fluctuate by several degrees centigrade. Nevertheless the actual temperature of the sample fluctuates only by a maximum of  $\pm 0.5$  °C.

## Other operation controls

On the front side of the temperature control unit there are covered shock-proof sockets for the connection of BOD's appliances or stirring apparatuses. The total load on the sockets must not exceed 4,5 amps.

The thermostatically controlled oven and laboratory sockets are protected by a fuse (6.3 A T) which is accessible from the front panel of the unit.

The control unit is connected by means of a 4-pole plug on the bottom of the incubator.

## Defrosting

When operating at a very cold thermostat setting, the interior rear wall may ice up.

In this case, the appliance must be defrosted manually.

### • Pull out the mains plug!

- Leave the door of the appliance open while defrosting. Defrosted water collects in an evaporator tray and is evaporated automatically. Attention must be paid that the defrosted water can drain off freely through the drainage opening in the back panel of the incubator. The drainage opening should, when necessary be cleaned with the aid of a suitable device (e.g. bottle brush). After defrosting mop up the remaining water and clean the appliance.

**Do not use any mechanical devices or other artificial aids for defrosting other than those recommended by the manufacturer.**

## Cleaning

Clean the appliance at least twice per year.

**Before cleaning always switch off the appliance. Pull out the mains plug or switch off or unscrew the fuse.**

- Clean the inside, equipment and outer walls with lukewarm water and a little detergent. Do not use abrasive or acid cleaners or chemical solvents.
- Ensure that no cleaning water penetrates into the electrical components or ventilation grille.
- The dust should be removed from the refrigeration unit and heat exchanger - metal grid at the back of the appliance - once a year.
- Do not damage or remove the type plate on the inside of the appliance. It is very important for servicing purposes.
- **Important!**  The compressor and other parts may still be hot from operation.

**Do not use steam cleaners because of the risk of injury and damage.**

## Malfunctions

If a malfunction occurs during operation, check whether it is due to an operating error.

**You may be able to rectify the following faults yourself:**

### • Appliance does not function:

- Is the appliance switched on?
- Is the plug correctly fitted in the mains socket?
- Is the fuse intact? If necessary, replace with a new 6.3 A (slow).

### • Loud running noise:

- Is the appliance set up firmly on the floor?
- Does the appliance cause nearby items of furniture or objects to vibrate? Please note that noises caused by the refrigerant circuit cannot be avoided.

### • Correct display of inside temperature but no regulation to the temperature setpoint:

- Is the temperature setting correct (see "Setting the temperature")?
- Is the ventilation system working properly?
- Is the appliance set up too close to a heat source?

If none of the above causes apply and you cannot rectify the fault yourself. Contact your dealer or supplier, if necessary send in the control unit or thermostat-controlled incubator for repair (original packing if possible!).

## Important

**Before undertaking any repair work disconnect plug from mains! Only trained specialists should work on electrical devices.**

## Shutting your appliance down

If your appliance is to be shut down for any length of time, switch it off and disconnect the plug or switch off or unscrew the fuse. Clean the appliance and leave the door open in order to prevent unpleasant smells.

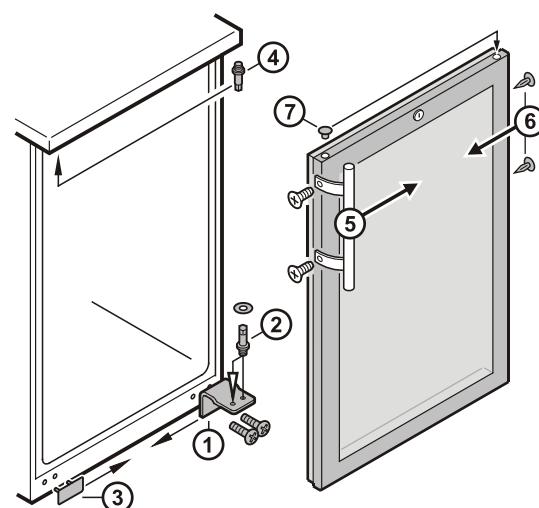
**The permissible room temperature for storing the appliance after decommissioning is between -10°C and +50°C.**

**Storing the appliance outside this temperature range can damage the appliance and cause it to malfunction.**

The appliance complies with the relevant safety regulations and EC Directives 2004/108/EC and 2006/95/EC.

## Changing over door hinges TC 140 G

- Unscrew hinge bracket ①.
- Pull the door out at the bottom and lift off.
- Transfer cover ③ to the opposite side.
- Transfer hinge pin ② on hinge bracket ① to the opposite side.
- Transfer hinge pin ④ to the opposite side.
- Transfer door handle ⑤ and plugs ⑥ to the other side.
- Mount door on hinge pin ④ and close.
- Insert hinge bracket ① in lower door mounting and screw into place.


## Changing over door hinges TC 256 G

Door hinges should only be changed by a trained expert.

Changing the door hinges must be done by two people.

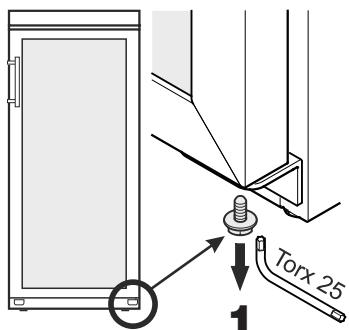1. Remove screw from lower hinge bracket.


2. Open the door approx. 45° and lift upwards.


**Important**

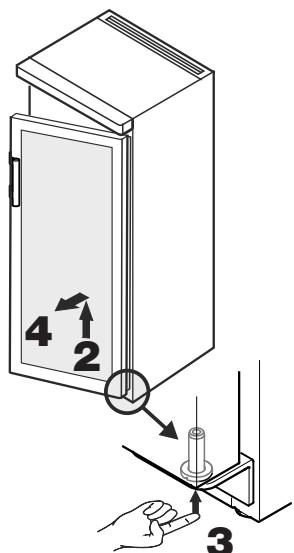As a result of its heavy weight, the door must be held secure by one person.

3. Push the pin upwards using your fingertip.

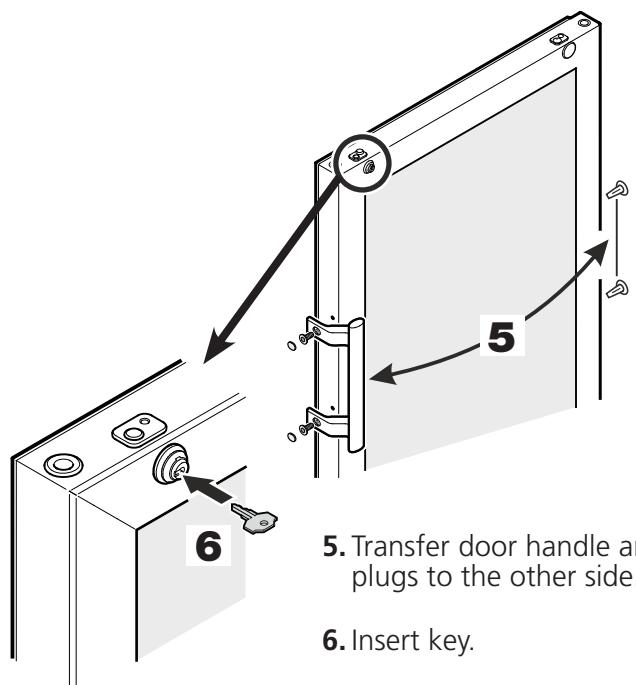
4. Pull the door to the front and remove it downwards.

5. Transfer door handle and plugs to the other side.


6. Insert key.

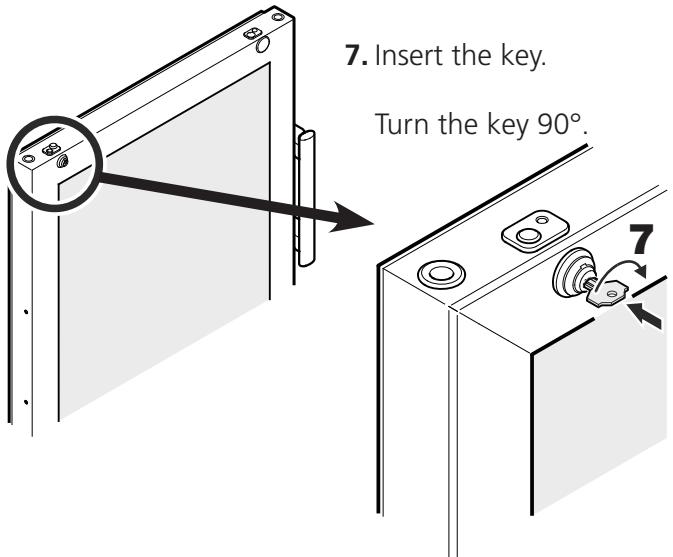7. Insert the key.


8. Remove cover.

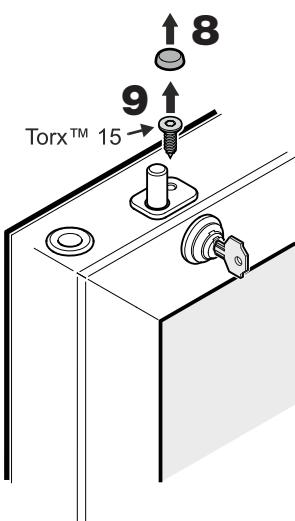9. Remove screw.


10. Turn the lock approx. 3° clockwise.

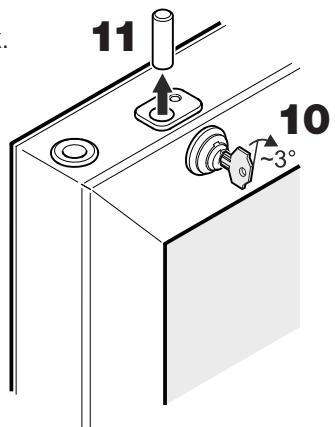11. Remove pin.


12. Remove bushing.

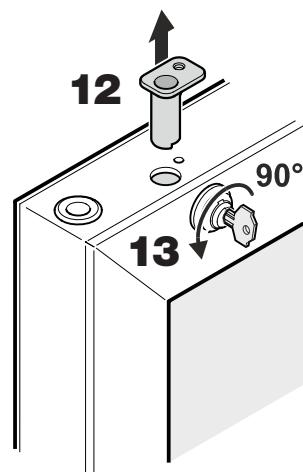13. Turn back the lock 90° anticlockwise.


**14.** Remove lock.


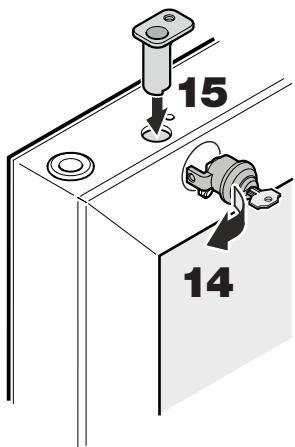
**15.** Reinsert hinge bushing.

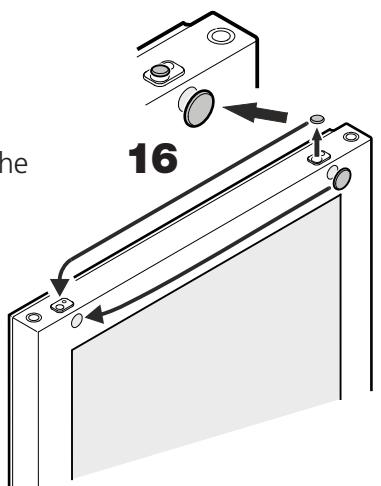**16.** Transfer covers to the opposite side.


**17.** Transfer plug and bushing.

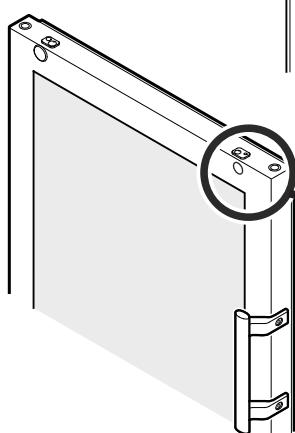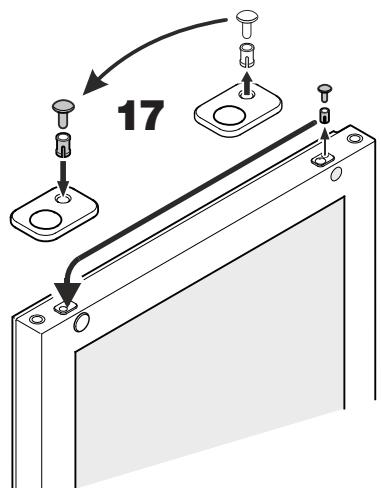
**18.** Remove bushing.

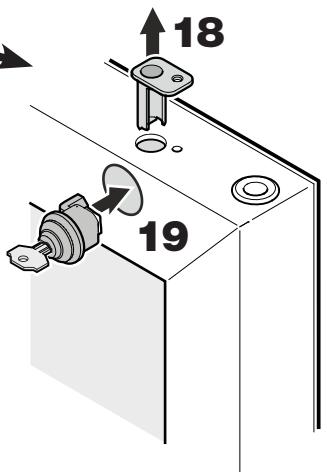**19.** Insert the lock as shown in the illustration.


**20.** Turn the lock 90° clockwise.

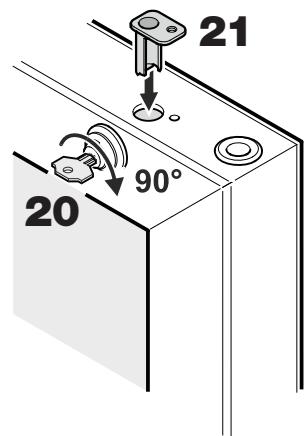**21.** Reinsert hinge bushing.


**22.** Turn the lock further - approx. 3° clockwise.

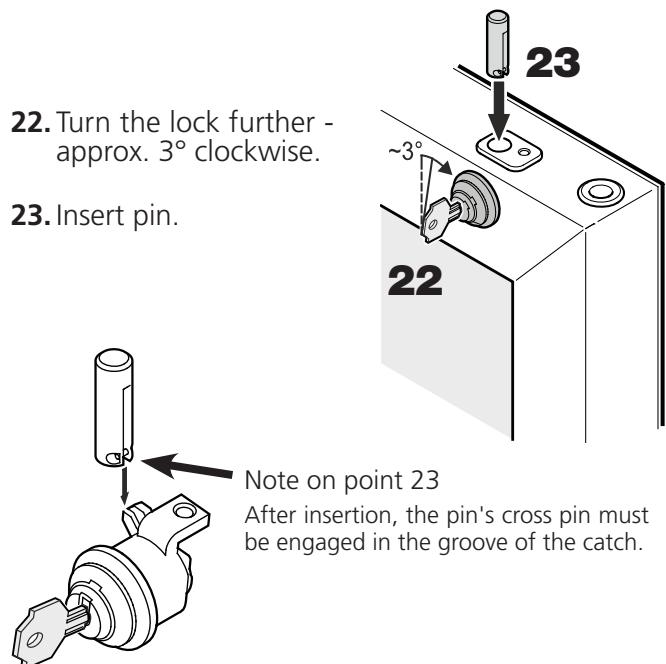**23.** Insert pin.


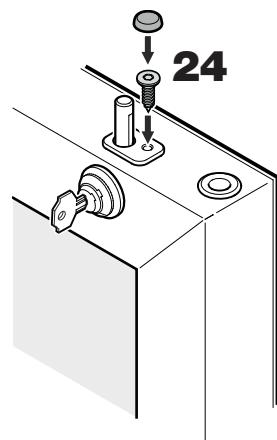
**24.** Screw in the screw and attach the cover cap.

**25.** Insert the key.

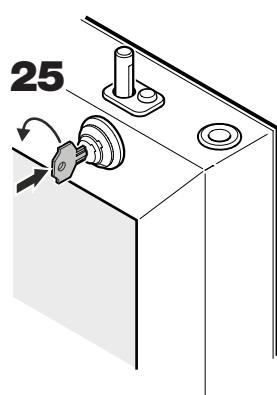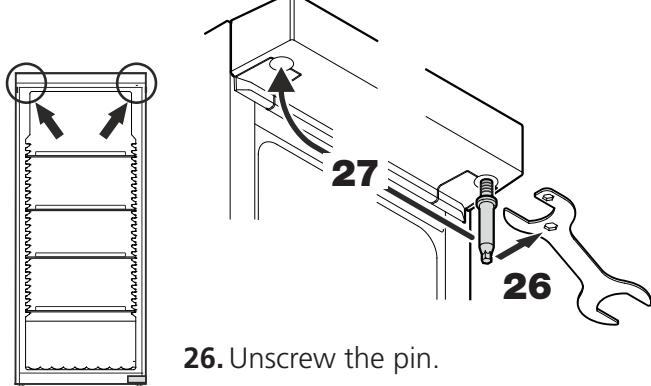Turn the key 90°.



**26.** Unscrew the pin.

Use the spanner provided.

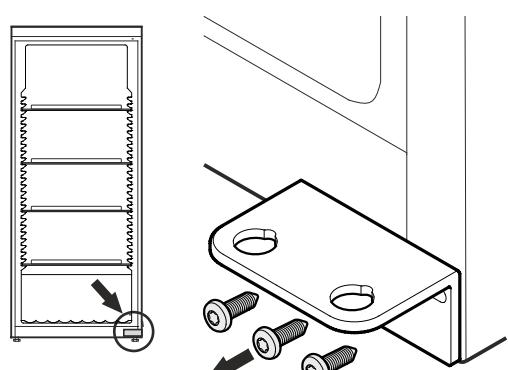
**27.** Screw in the pin on the opposite side.


**28.** Unscrew hinge bracket.

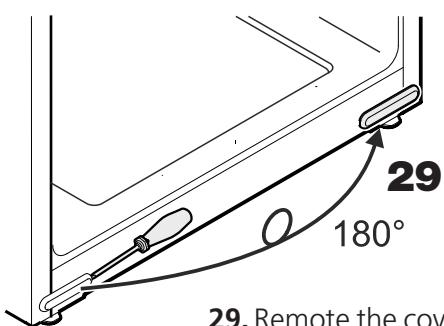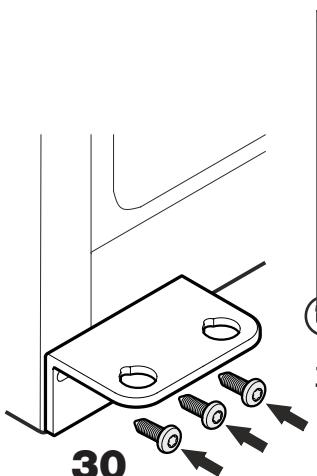**29.** Remove the cover, turn it through 180° and re-fit it on the opposite side.

**30.** Screw hinge bracket into place.

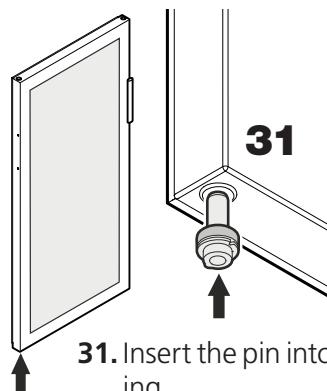**31.** Insert the pin into the left-hand door mounting.

Use your index finger to hold the pin while you are assembling the door.

**32.** Locate the door on the upper pin as shown in the illustration and guide it to the lower hinge bracket.


**Important**

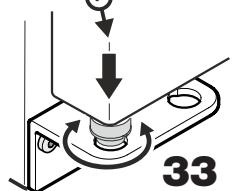As a result of its heavy weight, the door must be held secure by one person.


**33.** Lift the door upwards and turn the pin until it fits into the bore hole of the hinge bracket.

Slowly lower the door downwards. The lower part of the pin must be fully engaged in the bore hole of the hinge bracket.

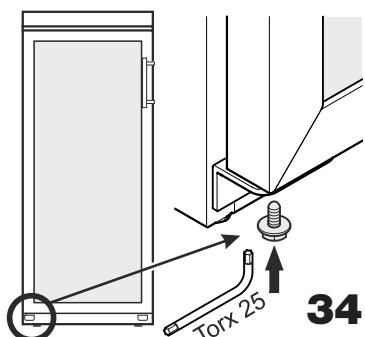
**34.** Screw the screw into the pin again and tighten.


## Adjusting the lateral tilt of the door

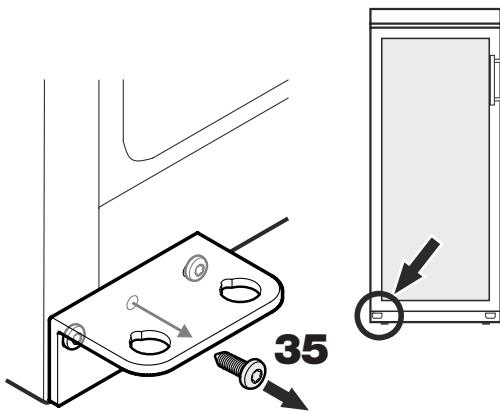If the door is at an angle, adjust the angle.

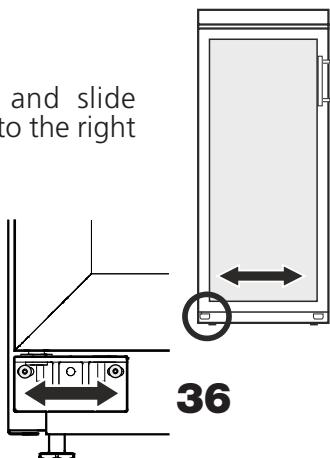- 35.** Remove middle screw from hinge bracket.  
This screw is no longer needed.

- 36.** Undo screws and slide hinge bracket to the right or left.

Tighten screws.

## Technical Data

### Construction

fully insulated incubator with universal plug-in temperature. Door, double glazed insulated in ABS frame

### Control unit operation

robust, water protected foil front panel, 2 keys with tactile acknowledgement  
2 °C to 40 °C, adjustable in 0.1 °C-increments

### Constancy of temp.

20 °C, ± 1.0 °C  
the climate rating is indicated on the type plate

### Climate class

3 digit for actual and target temperature, resolution 0.1 °C  
compressor-system

### Display

heating element

### Fan

axial, discharge capacity 320 m³/h

### Power requirement

220 - 240 V / 50Hz

### Sockets

CEE 7/5, type E, with hinged lid, 230 V / 16 A, 2p + E, IP 44

### Coolant

R134a

### TC 140 G (4 sockets)

140 Ltr.

### Net capacity

275 VA

1.77 kWh / 24 h  
(ambient temperatur 25 °C,  
target temperature 20 °C with interior lighting switched on)

### Overall dimensions

600 W x 600 D x 850 H mm with worktop

600 W x 600 D x 819 H mm without worktop

513 W x 441 D x 702 H mm (with 3 retractable grids and 1 bottom grid)

### Weight

48.0 kg

### Order Code

43 82 10

### TC 256 G (7 sockets)

255 Ltr.

### Net capacity

340 VA

1.56 kWh / 24 h  
(ambient temperatur 25 °C,  
target temperature 20 °C with interior lighting switched on)

### Overall dimensions

600 W x 610 D x 1640 H x mm

470 W x 440 D x 1452 H mm (with 4 retractable grids and 1 bottom grid)

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