



Short Instructions HandyStep® touch and HandyStep® touch S


Connect battery

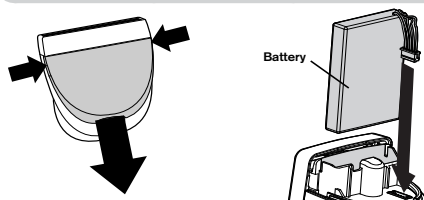


Select work mode

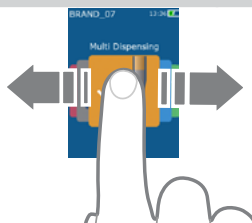


Set values

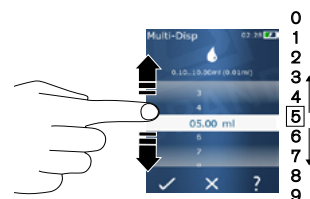


Charge battery

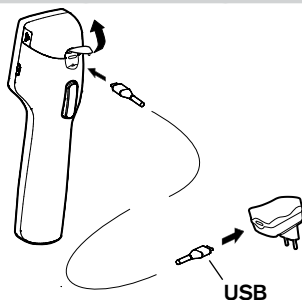


Tap on work mode to open it

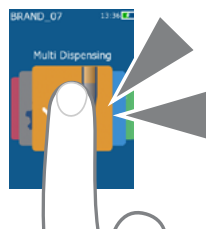


Switch digits

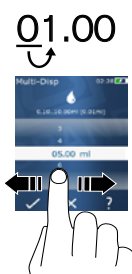

Turn on/off

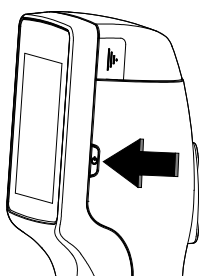

Insert tip

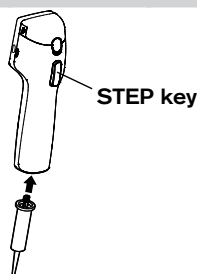

Confirm/abort settings



Safety

For safety information refer to operating manual HandyStep® touch and HandyStep® touch S. For current operating manuals refer to www.brand.de.

Menu reference

MULTI-DISP



Dispensing stepwise.

AUTO-DISP



Automatic dispensing. Interval settings between steps manually or via learn function.

PIP



Pipetting predefined volumes.

SEQ-DISP (+)



Dispense volumes based on predefined steps.

MULTI-ASP (+)



Aspirate volumes manually, stepwise or predefined with different step volumes.

TITRATION (+)



Dispense liquid manually or stepwise for titration.

SETTINGS



Device
Time/Date
Display
Energy
Sound
Firmware
Factory settings


GENERAL




MULTI-DISP




AUTO-DISP

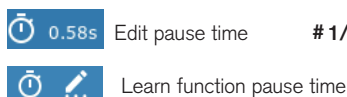


SEQ-DISP(+)

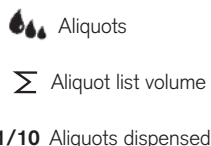

MULTI-ASP (+)

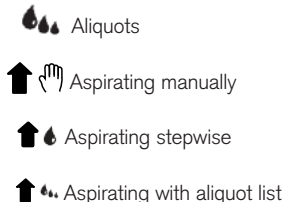

TITRATION (+)

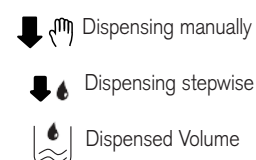

+ = HandyStep touch S only



888-522-2726 | www.brandtech.com

