Manual

CertoClav Classic 125°C/140°C

CertoClav Classic 115°C/121°C

Laboratory autoclave

Autoclave using steam for the treatment of material and laboratory processes e.g. sterilization, as an aid in research and development

English

German version page 13

Contents

CE Declaration of Conformity	3
Introduction.....	4
Warranty	4
Safety devices.....	5
What is the autoclave CertoClav Classic	6
<i>Names and functions of each part.....</i>	<i>6</i>
<i>Product uses.....</i>	<i>7</i>
<i>Product Features</i>	<i>7</i>
List of items	7
Installation instructions	7
Transportation	8
Operation.....	8
<i>Loading the autoclave</i>	<i>8</i>
<i>Closing and locking</i>	<i>8</i>
<i>Heating up</i>	<i>9</i>
<i>Sterilization / Holding time</i>	<i>9</i>
<i>Cooling down.....</i>	<i>9</i>
<i>Opening</i>	<i>9</i>
Cleaning and maintenance	10
Spare parts:.....	10
Technical data.....	11
<i>Overview.....</i>	<i>11</i>
<i>Stickers</i>	<i>11</i>
Electrical cable plan	12

CE Declaration of Conformity

PRODUCT: CERTOCLAV AUTOCLAVE CLASSIC

This is to certify, that the **CERTOCLAV** Product Range CertoClav Classic complies with the following regulations of the European Community:

2014/35/EU Low Voltage Directive

Associated Standards:

EN 61010-2-040:2005 Part 2: Particular requirements for sterilizers and washer-disinfectors used to treat medical materials

EN 61010-1:2001 Safety requirements for electrical equipment for measurements, control and laboratory use – Part 1: General requirements

The Certoclav pressure equipment is constructed according Low Voltage Directive.

2014/30/EU Electromagnetic compatibility

Associated Standards:

EN 61326-1:2013 Electrical Equipment for measurements, control and laboratory use - EMC Requirements- Part 1: General requirements

EN 61000-3-2:2006 + A1:2009 + A2:2009 Electromagnetic Compatibility (EMC) — Part 3- 2: Limits

Limits for harmonic current emissions (equipment input current equal to or less than 16 A per phase)

EN 61000-3-3:2013 Electromagnetic compatibility — Part 3- 3: Limits — Limits for voltage fluctuation and flicker on the public low-voltage system (equipment having an input current equal to or less 16 A per phase) and not subject to conditional connection.

2011/65/EU Restriction of the use of certain hazardous substances (RoHS) Directive

Associated Standards:

EN 50581:2012 Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances

Quality of Production:

CertoClav produces according modern production methods and with high quality materials. Experience for many years and a competent team ensure high quality of products and professional service. All units are 100 % tested on function and quality. Certoclav manufactures in Austria only.

Manufacturer:

CERTOCLAV Sterilizer GmbH

Georg Grinninger-Str. 37

A-4050 Traun / Austria

Traun, 21.11.2016

Mag. Moritz Gruber
CEO

Introduction

Thank you for choosing CertoClav! Since 50 years we stand for quality made in Austria. Should the product not fulfil your expectations please contact your retailer or CertoClav directly via email to support@certoclav.com.

The CertoClav Classic laboratory autoclave regulates heat through a combination of a thermostat and a precision valve. The regulation is calibrated to pure steam, required proper evacuation of air. Time can be conveniently regulated via the built in timer.

Please obey the safety remarks

Please register your unit via email to support@certoclav.com.

Please send name, contact data and serial number of the unit.

Your

CertoClav team

Warranty

The unit has been checked from our quality control by 100 %. We give guarantee for 2 years on all parts which are not consumables from the date of purchase.

An extension of the warranty period to 5 years is available for purchase.

No warranty will be given in case of inappropriate use e.g. corrosive loads.

Safety devices

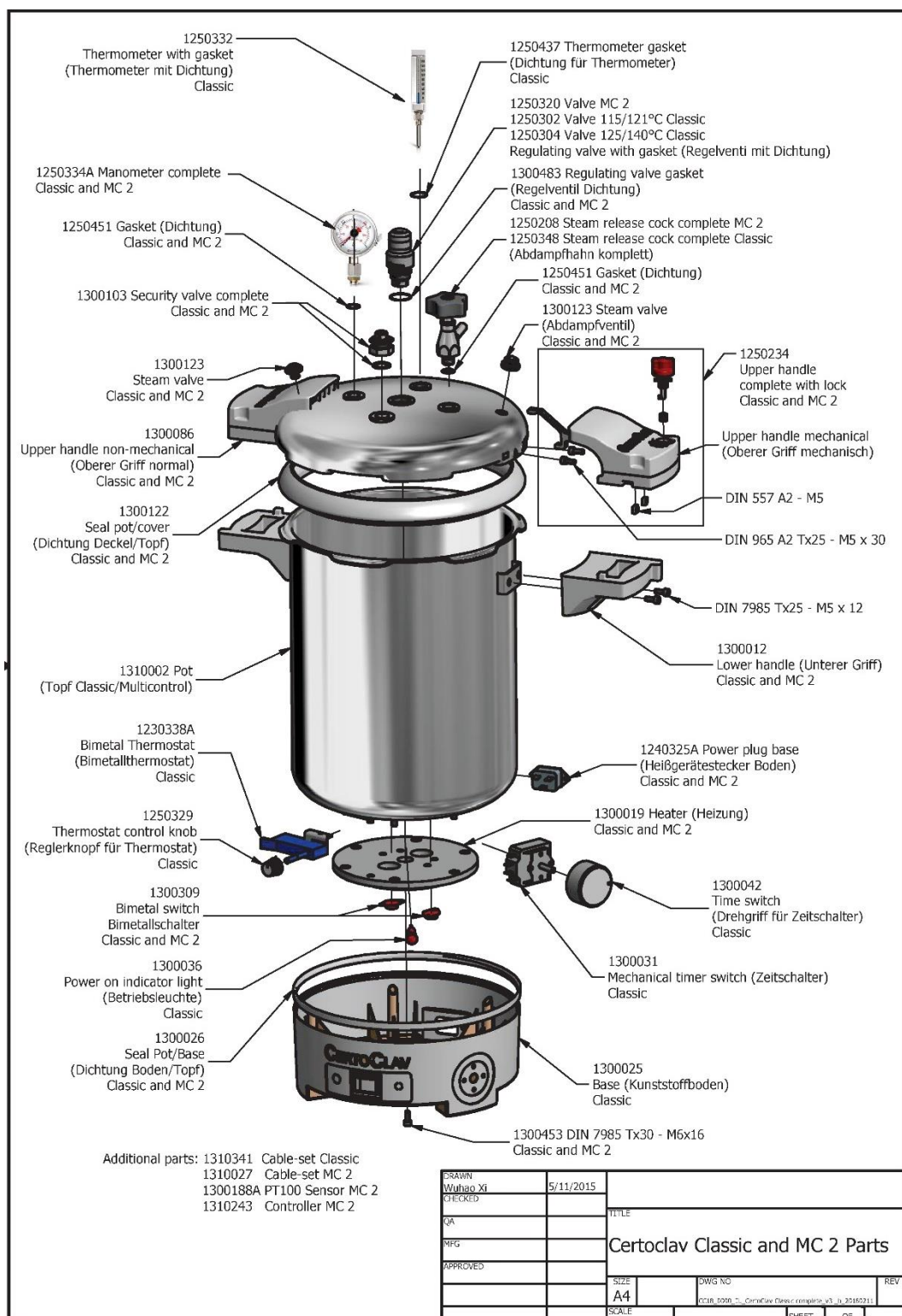
In this manual, the following headings are applied to items to which attention should be given:

	Warning	Precaution indicating an imminent dangerous situation which if not avoided may lead to death or serious injury.
	Caution	Precaution indicating a dangerous situation which if not avoided may lead to moderate or slight injury.
	Important	Indicates items you are strongly advised to obey.
	Note	Items that will aid in proper operation of the equipment.

-  Never manipulate on safety devices. Operate only controller and chamber of the same series.
-  Do not mix lid and chamber of different types of autoclaves.
-  Surfaces are hot while in operation (metal and plastic parts). Wear protective gloves.
-  Use only in accordance with this manual. Store the manual in a space accessible to all users.
-  Must not be installed in a space that is subject to the danger of explosions
-  Use only with rated voltage.
-  Only for use in laboratories. Operate only by trained personnel. Do not use in domestic settings.
-  Never operate without water. Target water amount: 1,5 liters (min.1 liter)
-  Do not autoclave tightly sealed containers (especially glass bottles) since there is no pressure compensation.
-  Use only water for steam creation.
-  Do not release steam in the direction of a person – danger of burning. We recommend to connect a temperature resistant hose (inner diameter 10mm) to release hot steam into a suitable container or sink. We offer this hose and container as accessory: Art. no. 8582016 – hose 1.0m
-  If the pressure is released quickly when autoclaving liquids, there is a possibility that the liquid will boil. I this should be avoided we recommend cooling the autoclave to 70°C before opening the valve.
-  If the running time too long it is danger of cooking dry and overheating. The autoclave could be damaged. Check regularly water level.
-  Do not transport when filled. Do not open with force.
-  This autoclave must only be maintained by CertoClav or authorized retailers. Always use original spare parts.

What is the autoclave CertoClav Classic

Names and functions of each part

Product uses

The product is used to sterilize substances which can withstand high temperature and high pressure steam such as tools of glass, ceramic, metal or rubber, water, media and reagents.

Product Features

- Integrated timer
- Precision thermometer
- Precision pressure gauge
- Process control via thermostat and precision valve
- Steam release cock
- Volume: 18 litre
- Max. Temp.: 140°C
- Max. pressure: (rel.): 2.7 Bar
- 220-240VAC, 50Hz, 1300W
- Certificates: CE, TÜV
- Operating time: S2 max. 30 min

List of items

The following parts are delivered with the autoclave, please check:

1. Autoclave CertoClav Classic with lid
2. Manometer and steam release cock with gasket and nut
3. Thermometer with gasket
4. Mains cable
5. In a poly bag: manual and warranty certificate
6. Instrument tray with water level indicator

Should one of these parts be missing, please contact your dealer or call us directly.

Additionally, we recommend the following accessories:

Wire basket diam. 25, stainless steel ø 25 cm. Art. No. 1300149

For further information about our product range please visit: www.certoclav.com/shop

Installation instructions

1. Either put the autoclave on a table or the floor, so that you get a convenient work height.
2. Unpack the manometer and the steam release cock and mount them with gaskets by putting in the holes and fix it with nut from inside of lid. These appliances have been packed separately to avoid damage in transportation.
3. Mount the thermometer with gasket according to the instructions packed with.
4. Mount a tube/hose on the steam release valve, inner diameter 10 mm.
5. Plug in the mains cable to the socket on the back side.

Transportation

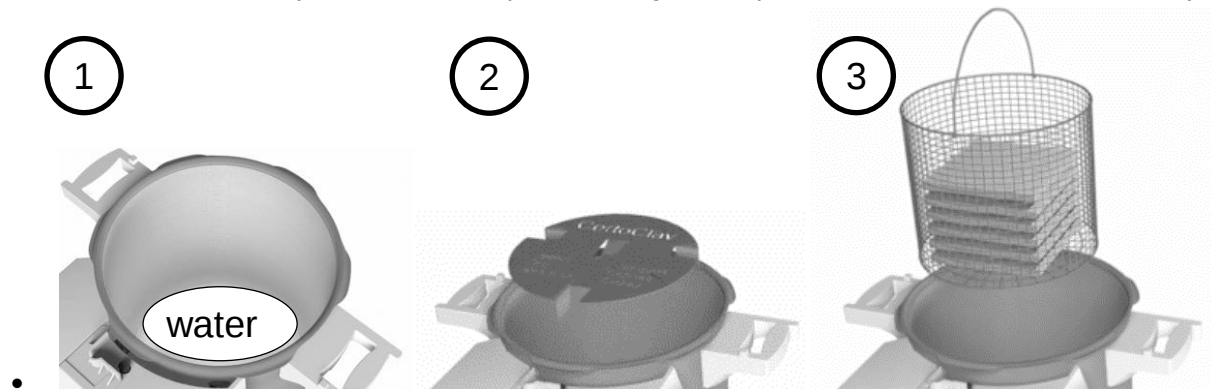
1. Check if the autoclave is cooled down and the chamber is completely empty.
2. When transporting the equipment, unmount the manometer and steam release cock and put them into a separate shock resistant box.
3. Close the lid and slide the open/close lever into locked position to prevent the lid from opening

Operation

- ⚠ Read instructions before first use
- ⚠ The first run should be without load; only with water.
- ⚠ Do not autoclave tightly sealed containers (especially glass bottles) since there is no pressure compensation.

Loading the autoclave

- Fill in 1.0 – 1.5 litres of water.
- Put in instrument tray and basket and place loading. It is important that steam can circulate freely.

Closing and locking

- Take the lid and make sure that the red knob is turned to open.
- Lid is put on 30° turned counter clockwise. Then turn the lid clockwise until the handles are in line with each other.
- Turn the red knob to closed = locked.
- Open the steam release cock (turn black knob anti-clockwise).
- Optional: Connect tube to steam release cock.
- Turn the red maximum indicator of the manometer (pressure gauge) to "0".
- Pressure control valves with two temperatures have to be set to the desired temperature by rotating cup until it comes to a stop. Anticlockwise to 125°C / 115°C – and clockwise to 140°C / 121°C. Do NOT open the small screw on the regulating valve.
- Set the black thermostat knob to the same temperature as sterilization valve at the lid.

Heating up

Switch on the autoclave by turning the white timer knob clockwise to its always-on position.

If timer knob is set a position for continuous operation in this case the autoclave has to be switched off manual. The time can be set after temperature is reached with timer. In this case the autoclave switched off automatically after time run off.

15 – 20 minutes after switch on steam escapes through the steam release cock.

To insure sufficient air removal let the steam release cock open for 10 minutes than close it by turning the black knob clockwise. Saturated steam will be built.

Now pressure and temperature rise according adjustment.

Sterilization / Holding time

- The sterilization time starts when temperature is reached. Steam could escape through the pressure control valve. The indicator on the manometer should stay within the specified limits (red or blue marking).
- You can now use the timer to set the holding time you need (e.g. 15 minutes to sterilize 500ml culture media at 121°C)

Load of fluid in ml	Initial timer switch position
3000 ml	30 minutes
2000 ml	25 minutes
1000 ml	20 minutes
500 ml	15 minutes

- Pressure regulation is done by the pressure control valve, steam escapes in regular intervals.
- After the end of sterilization time switch off by turning the timer knob to OFF. If you have used the timer function, the unit switches off automatically.
- If timer function is not active, you have to switch off manual. Turn the knob to OFF.

⚠ If timer function is not active don't forget to switch off, the autoclave could cook dry and overheat. Make sure that enough water is in the chamber.

⚠ Energy is controlled by the thermostat. Make sure that the same temperature on valve and knob for thermostat on the base is set, they work together.

⚠ On CertoClav Classic autoclaves energy is controlled by the thermostat. Make sure that the same temperature on valve and knob for thermostat on the base is set, they work together.

Cooling down

Check that the manometer whether the black indicator shows to Zero. Check temperature at thermometer. For solid loads, please wait until temperature is lower than 80°C. For fluid loads, please wait until temperature is lower than 70°C.

⚠ SAFETY HINT: If loading contains glass bottles wait until chamber temperature is below 80°C, for liquids wait until temperature has reached 70°C = approx. 40 min. Danger of boiling over of fluids.

Opening

- Open the steam release cock (turn anticlockwise).
- Push & turn red knob to open position. Steam release valve will be opened. Steam escapes.
- Turn lid anticlockwise and remove.
- To get better drying of the loading remove the basket from the chamber and let it cool down to room temperature before taking goods out.

⚠ SAFETY HINT: INCOMPLETE CYCLES

Should the cycle interrupted caused by switch off or electrical failure or other reason – the sterilization is not completed correct and has to be repeated. For opening please follow the safety hint for cooling down.

Cleaning and maintenance

For CLEANING the appliance wipe inside and outside with a damp cloth. Never immerse the lid and chamber of the unit in water.

Before start of cleaning or disinfection plug off mains cable and make sure that appliance has cooled down.

For DISINFECTION of surface disinfection liquids with following substances are suitable:

Ethanol, Propyl alcohol, Isopropyl alcohol, Glutaraldehyde, quaternary ammonium compounds.

Please follow the application notes of the manufacturer for the cleaning and disinfection detergents.

Discolouration and stains occur when not using distilled water. These stains may block the pressure control valve. Therefore, always use distilled water or demineralised water with tap water added.

To remove discolouration, add some citric acid to the aqua dest. and boil for about 10 minutes, do not close the CertoClav. Empty the chamber immediately.

Never manipulate on safety devices. Operate only lid and chamber of the same series (Classic or MultiControl). Do not mix.

Repairs on valves, safety features and electric must only be carried out by an appointed CertoClav partner or by the manufacturer. Use original CertoClav spare parts only.

Every two years the unit should be checked by an appointed retailer or CertoClav.

Spare parts:

CERTOCLAV SERVICE SET 2 – Art.Nr. 1310186.

It contains all spares which are under wear and tear and have to be replaced.

Technical data


Overview


Technical data			
Material	Stainless steel	Certificates	CE, TÜV
Dimensions			
Volume	18 litre	Height	60 cm
Chamber diameter	26cm	Width	44 cm
Chamber height	35cm	Depth	40 cm
Useable height chamber	30 cm	Weight	10 kg
Working data			
Max. working temp	140°C	Current	230V, 50/60 Hz
Max. working pressure	2,7 Bar	Power	1300 Watt
Rated pressure	4,1 Bar	Operating time	S2 max. 30min

Stickers

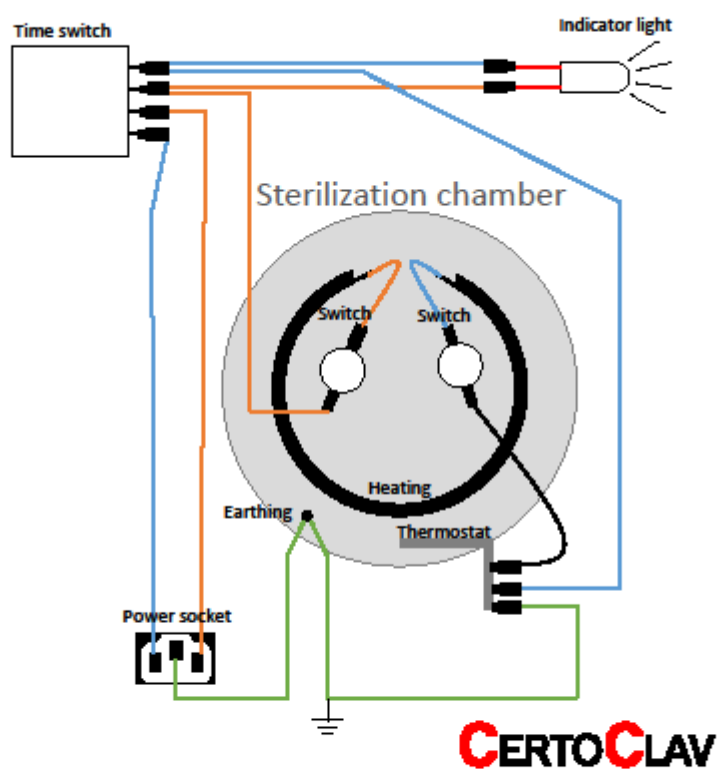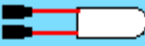	Directive 2002/96/EC (WEEE): The supplier will take the supplied commodity back on the expense of customer after use completion and ensure best available treatment, recovery and recycling according to the laws duly.
	ATTENTION: Consult manual before use
	ATTENTION: Hot surface
	CE certified according to: Low voltage directive 2014/35/EU + EMC 2014/30/EU.

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