Temperature

Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Test Tube Shakers Hei-MIX Reax top/Hei-MIX Reax control
Operating manual must be read before initial start-up.
Please follow the safety instructions provided.
Please keep for future reference.

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About this Document	4
Version and Variants	4
About this manual	4
Terms and signs used.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / Electrical safety	7
Personnel qualification.....	7
Operating company's obligations.....	7
Installation site	7
Changes to the unit	8
Safety for the personnel	8
Safety during use	8
Disposal	8
Device Description	9
Device overview	9
Hei-MIX Reax top	9
Hei-MIX Reax control	9
Start-Up	10
Setting up device	10
Connect / disconnect power cord	10
Switch unit on and off	11
Operation	12
Shaking mode.....	12
Automatic mode	12
Continuous mode.....	12
Speed / Shaking intensity	13
Set Speed.....	13

Cleaning and Maintenance	14
Cleaning.....	14
Maintenance	14
Troubleshooting	15
Possible errors and how to resolve	15
Montage	16
Mounting test tube tray	16
Electrical connections.....	17
Disassembly and Storage	18
Dismantling, transportation and storage	18
Dismantling	18
Transportation and storage	18
Accessories and Spare Parts	19
Scope of delivery.....	19
Accessories.....	19
Attachments	20
Technical data	20
Sevice	21
Contact / Technical Service	21
Warranty	22
Confirmation of condition of unit.....	23

➤ Version and Variants

Version

This manual describes the function, operation and maintenance of the Hei-MIX Reax top and Hei-MIX Reax control test tube shakers.

Version	Alteration date
0.3	03/2017

Variants

The units are available in different configurations. Certain features and functions apply to specific configurations only. The varieties are documented in this manual.

➤ About this manual

This operating manual is a component of the device described here.

- Please read this manual carefully and obey all safety and warning notices.
- Ensure that every operator reads this manual.
- Ensure that this manual is accessible for every operator.
- Pass on the operating manual to the subsequent owner.

For the current version of this operating manual in pdf format, please go to:

<http://www.heidolph-instruments.com/Support/Operation Manuals/Shakers & Mixers/Reax and Multi Reax/Reax top and Reax control>

In addition please observe the regional regulations.

➤ Terms and signs used

In this manual you will find standardized terms and signs to warn you of possible dangers or give important hints. Please take special notice to these signs to avoid personal injury or damage to the unit.

The following terms and type of signs (basic drawing elements) are used:

Symbol	Additional term / Description
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following terms: DANGER: Indicates a hazardous situation which, if not avoided, will result in irreversible injury or death. WARNING: Indicates a hazardous situation which, if not avoided, could result in serious injury or material damage. CAUTION: Indicates a hazardous situation which, if not avoided, may result in a minor injury or material damage.
Prohibitory signs 	Do not do: The red circle indicates a situation that should be avoided under all circumstances to avoid injury or damage.
Mandatory signs 	Must do: The blue circle indicates important information. Please obey to avoid property damage.

Other signs used

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪ ▪ ▪	List of information
a. b. c.	List of variants

➤ **General safety instructions**

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see attachment „EU Declaration of Conformity“.

The device has been constructed according to state-of-the-art technology and recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure the operating manual is available at all times.

The device may only be used under the following circumstances:

- Only operate the device, if it is in full working order.
- Ensure all operators of the device possess the necessary safety and risk awareness.
- Operate the device according to the instructions stipulated in this manual only.
- If there is something you do not understand, or certain information is missing, ask your manager or contact the manufacturer.
- Do not do anything on the device without authorization.
- Only use the device according to its intended use.

Intended use

The device is intended for use by trained and authorized personnel only.

The device is suitable for the following use:	<ul style="list-style-type: none"> • mixing • shaking • blending of solvents
The operation for the device is suitable in one of the following locations:	In any research application

Improper use

Any use which deviates from the device's intended use is considered to be improper. The manufacturer does not accept liability for any damages resulting from non-permitted uses. The risk is carried by the operator alone.

Installation / Electrical safety

- The device may only be connected when the mains voltage corresponds to the information on the type plate of the unit.
- The mains connection must be accessible at all times.
- Repairs may only be performed by a qualified electrician.
- Never operate the unit with a damaged power cord.
- Always turn the unit OFF and disconnect mains power before performing any maintenance or service.

Personnel qualification

- The device may only be operated by qualified persons.
- The device may only be operated by individuals who have been instructed in its proper use by qualified persons.
- The device may only be operated and maintained by persons who are of legal age.
- Other personnel may only work with the unit under continuous supervision of an experienced qualified operator.
- This manual must be read and understood by all persons working with the device.
- The personnel must receive special safety instructions in order to guarantee responsible and safe work procedures.

➤ Operating company's obligations

Installation site

- The device must be positioned in a suitable location.
- The device must be installed sufficiently stable.
- All screw connections must be securely tightened.
- Do not operate the device in proximity of highly inflammable and explosive substances.
- Installation and operation of the device is only permitted in facilities which are fitted with the appropriate laboratory equipment (e.g. with air extraction units).
- We recommend operation in fume hood enclosures.
- The device may only be operated in enclosed spaces and under the following environmental conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decrease linearly till max. 50 % relative humidity
Installation altitude	0 - 2,000 m (6,500 feet) height above sea level
Degree of pollution	2
Overvoltage category	II

Changes to the unit

- No unauthorized changes may be made to the unit.
- No parts may be used which have not been approved by the manufacturer.
- Unauthorized changes result in the EC Declaration of Conformity losing its validity, and the unit may no longer be operated.
- The manufacturer is not liable for any damage, danger or injuries that result from unauthorized changes or from operating the unit other than as described in this manual.

Safety for the personnel

- Ensure that only qualified personnel operates the device.
- Observe the following regulations:
 - Laboratory guidelines
 - Accident prevention regulation
 - Ordinance on Hazardous Substances
 - Other generally accepted rules of safety engineering and occupational health
 - Local regulations

➤ Safety during use

- Wear the appropriate protective clothing when working on the device (clothing, protective glasses and if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Eliminate errors immediately.
- Do not use abrasive material to clean the glass surface. Only wipe with damp cloth.
- Always switch the device off after use.
- Observe the Directive 2002-44 / 2003-10 EC regarding the exposure of workers to the risks arising from physical agents (vibration).

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device according to the appropriate national legal regulations.
- Dispose of the packaging material in accordance with the appropriate national legal regulations.

➤ Device overview

Hei-MIX Reax top

Hei-MIX Reax control

➤ Setting up device

Warning:
Risk of slipping device!

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Connect / disconnect power cord

Connect power cord

The power cord comes with a three-wire plug and a recess on the bottom side.

Unit plug

Power cord

The socket is located on the back side of the unit.

- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.


➤ Switch unit on and off

The ON/OFF switch is situated on the front panel.

Switch unit on


- Press the On/Off switch on the lower side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.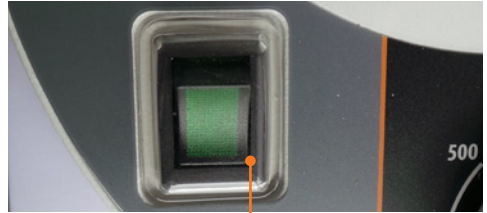

Switch On: device switched on

Switch unit off

- Press the On/Off switch on the upper side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

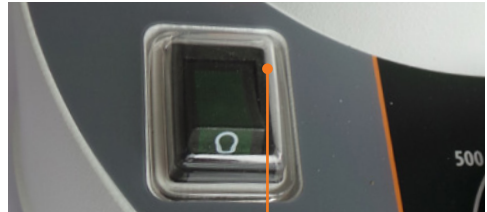Switch Off: device switched off

➤ Shaking mode

You can choose from two shaking options:

- Automatic mode: The unit starts shaking as soon as you put pressure on the test tube tray.
- Continuous mode: The unit shakes at an ongoing vibration mode until you switch off the continuous mode or the complete unit.

Warning: **Risk of injury!**

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

Automatic mode

- Turn the slide switch to the right as far as it goes.
- ✓ A delta icon ▼ appears on the left side of the slide switch.
- Place your tube in the test tube tray and push the tube with soft pressure downwards.
- ✓ The shaking operating starts as soon as you position a tube on the test tube tray.
- Remove the tube.
- ✓ The shaking operating stops.

Symbol automatic mode

Slide switch

Continuous mode

- Turn the slide switch to the left side as far as it goes.
- ✓ A circle icon ● appears on the right side of the slide switch.
- ✓ The shaking operating starts.
- Turn the slide switch to automatic mode or switch off the unit.
- ✓ The shaking operating stops.

Slide switch

Symbol continuous mode

➤ Speed / Shaking intensity

Set Speed

Warning:
Risk of splashing!

Open vessels and too high vibration frequencies may result in samples splashing.

Wear safety glasses and appropriate protective clothing.

Select low speed when starting the unit to avoid splashing from tubes. If necessary, increase the speed carefully.

Caution:
Possible risk of exertion

There is a possible risk of exertion due to physical force required to holding test tube in place.

Do not use the device any longer than recommended in the directive 2002-44 / 2003-10 EC on vibration strain.

REAX top

- ➔ Turn the speed rotation knob clockwise to the right.
- ✓ The shaking intensity increases.
- ➔ Turn the speed rotation knob counter-clockwise to the left.
- ✓ The shaking intensity decreases.

The Reax top is equipped with a guideline to set the speed.

REAX control

The Reax control is equipped with an analog dial speed control with accurate rpm numbers setting.

➤ Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Attention:
Damage to the surfaces

All surfaces can be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent including chlorine
- Ammonia
- Abrasive cleaning agents such as cleaning rags, scrubbing agents or any other agents which include metal components

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

➤ Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

Troubleshooting

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> • Check size and shape of plug and compatibility to your electrical socket • Check circuit breakers
	Light diode defective	Contact your local Heidolph Instruments distributor
No shaking movement in tilt automatic mode	Slide switch not in correct position	Push slide switch to the far right position
	Not enough pressure on the tray	Increase pressure by hand
Shaking movement stops in continuous operation mode	Pressure on tray not applied vertical	<ul style="list-style-type: none"> • Apply vertical pressure • Switch over to continuous operation mode
	<ul style="list-style-type: none"> • Motor defective • Switch defective • Control electronics defective 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

➤ Mounting test tube tray

Take off test tube tray

The test tube tray can easily be attached and pulled off the unit manually without using a tool.

- Grab the test tube tray and lift it upwards to pull it off the unit.


Attach test tube tray

A groove is located on the bottom of the test tube tray.

- Place the test tube tray with groove in line with the tongue of the unit.
- Push the test tube tray with soft pressure onto the unit.

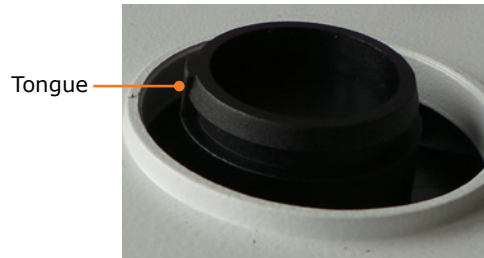Electrical connections

- Repairs of any kind are allowed by **qualified professional electricians only**. Any improper repair can result in a dangerous situation. Contact your local Heidolph Instruments distributor for any repair you may have.


Warning: Electric shock!

This unit must only be connected to a grounded electrical socket.

The unit and voltage must match. The type label on the back side of the unit provides all voltage specifics.

When shipped, the unit is grounded. In case the original plug is changed, the new plug must have a protective conductor!

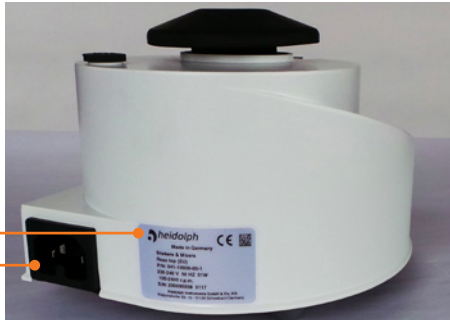Rating plate

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Color code for electrical connections:

Europe

GREEN/ YELLOW	PE: Protective conductor (Earth)
BLUE	N: Neutral conductor
BROWN	P: Phase

North America

GREEN	PE: Protective conductor (Earth)
WHITE	N: Neutral conductor
BLACK	P: Phase

The unit is connected with the power cord supplied to the electrical socket. It plugs into the housing on the rear side of the unit.

For countries utilizing a plug other than the standard one supplied:

- The plug supplied may only be changed by a professional electrician.
- If utilizing an adaptor ensure that it is grounded and approved by the local regulations.

➤ Dismantling, transportation and storage

Dismantling

**Warning:
Risk of injury!**

While the unit is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling unit ensure that the unit is switched off and the plug is pulled from the electric socket.

- ✓ The unit is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the unit.

Transportation and storage

- Store and transport the unit and its components only if they were emptied and cleaned,
- Store and transport the unit and parts of unit in the original packing material or alternatively in an adequate container to prevent damages.
- Seal the packing carefully against unauthorized or mistaken opening.
- Store the unit in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Variant	Quantity	Product number
Hei-MIX	Reax top	1	541-10000-00
	Reax control	1	541-11000-00
Tray, small		1	23-07-06-05-01
Power cord		1	14-007-003-81
Operating Manual	English / German	1	01-005-002-74
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06

➤ Accessories

Component	Quantity	Product number
Tray, large	1	549-19000-00

Additional accessories are shown in the general catalog or on our website at

- www.heidolph-instruments.com.

➤ Technical data

Hei-MIX Reax top, Hei-MIX Reax control

Standard supply voltage	230 V/50 Hz or 115 V/60 Hz
Supply power (W)	51
Protection class (DIN EN 60529)	IP22
Drive	shaded pole motor
Overheat protection	self-resetting
Speed control	electronic
Speed range (U/min)	0 - 2,500
Motion	circular vibration
Orbit (mm)	5
Operating mode	automatic or continuous
Weight (kg)	2.8
Dimensions (w x h x d) (mm)	134 x 105 x 172
Permissible ambient conditions	<ul style="list-style-type: none"> ▪ 5 - 40 °C ▪ 80 % relative humidity, no condensation at 5 - 31 °C, linear decreasing to 50 % at 40 °C

Model Hei-MIX	Reax top	Reax control
Speed setting	analog, +/- scale	analog, numbered scale
Control accuracy	unregulated	+/- 2%

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning: Danger of poisoning!

Contaminated units can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
 - take adequate protective action towards our parts receiving and service personnel
 - mark the pack in accordance with Ordinance on Hazardous Substances
-

Included at the end of this manual you will find a „Confirmation of condition of unit“.

→ Kindly copy and fill in this form and submit it prior to shipping the unit for repair.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 - 9122 - 9920-74
Fax: +49 - 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.de

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at
www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3-year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty is valid from the point of registration. The serial number is also valid without registering.

The warranty covers part and manufacturing defects.

In the case of a part or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damages incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please inform Heidolph Instruments or your local Heidolph distributor should you wish to make a warranty claim.
-

➤ Confirmation of condition of unit

→ In the case of repair, copy and complete the Confirmation of condition of unit and send it to Heidolph Instruments.

1. Details about the unit

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the unit in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the unit come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date _____ **Signature** _____

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG