

Supplement to the manual

Use of the Port Splitters RJ45-HUB

Usage

The port splitter RJ45-HUB allows you to connect your ultrasonic processor to various additional components and to use all features of your digital ultrasonic processor optimally. At the same time, the functionality of the network is retained. The RJ45-HUB is compatible with all of Hielscher's digital ultrasonic devices (featuring the touch control).

Four ports with the following functions.

- UIP port: for the connection to the Ethernet port of the ultrasonic processor
LAN port: for the connection to the user's local network
PS7 port: for the connection of the optionally available pressure sensors PS7, PS70 and PS140. The sensor can be easily activated in the setting menu of the ultrasonic device.
ON/OFF port: for the manual remote control. The port can be easily activated in the setting menu of the ultrasonic device.

The port splitter RJ45-HUB allows you to connect with various tools and gives you thereby a better process control and a higher working comfort.

<div data-bbox="662 1176 774 1272"></div> <div data-bbox="180 1234 422 1276">Please note!</div> <div data-bbox="180 1294 952 1471"><p><i>The port splitter RJ45-HUB (1) must be used when the device is operated at the network connection (TCP / IP, laptop, PC, LAN, switch, router, company network). For operation with only PS7 or only foot switch, this is not absolutely necessary.</i></p></div> <div data-bbox="180 1503 903 1574"><p><i>The cross-link adapter (2) is only required for very old network cards (10Mbit).</i></p></div>	<div data-bbox="997 1176 1332 1384"></div>
---	--

Address

Hielscher Ultrasonics GmbH
Oderstrasse 53
D-14513 Teltow
Germany

Telephone +49 (0) 33 28 / 437 3
Fax +49 (0) 33 28 / 437 444
E-Mail service@hielscher.com