





The Belly Dancer® Nesting Platform Installation

Warning: Prior to starting, ensure unit is unplugged from electrical source

To install The Belly Dancer® nesting platform you must first remove the existing standard Belly Dancer® platform. To do this first, remove the four screws attaching The Belly Dancer® platform to the flexible stabilizer legs at each corner. Use the provided hex wrench enclosed in the kit, save the detached screws for later use. Now, remove the three phillips head screws from the center of the acrylic platform. Finally, remove threaded standoffs from the bottom of acrylic platform, you may have to move them back-and-forth a few times to get them to release. Remove the standard acrylic platform.

1st

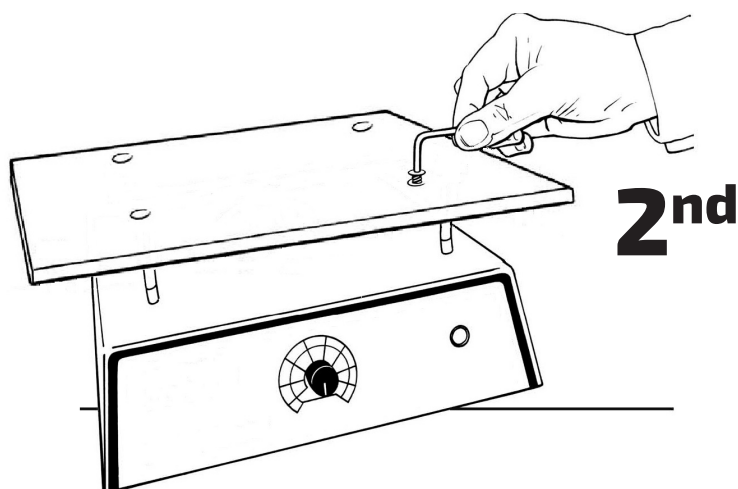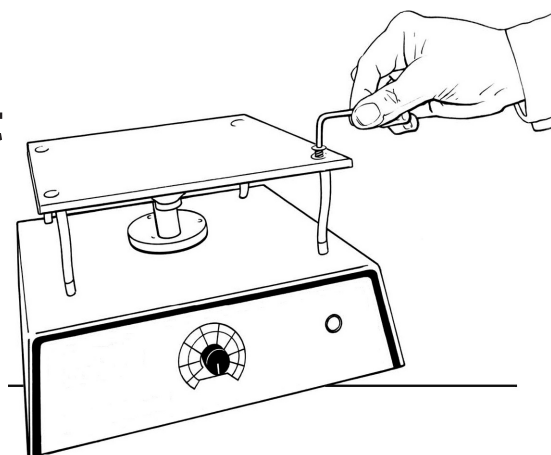

2nd

Second, install the nesting platform onto the top of the Belly Dancer® and align the (3) holes at the center of the platform with the (3) holes on the top of the tilt assembly. Replace the (3) phillips head screws previously removed from the standard platform assembly. It typically works best to start all (3) screws first, then tighten them once the platform is aligned properly. Next, align the fitting at the top of the flexible legs with each corner of the nesting platform. Insert the hex head screw and thread into the fitting until the fitting below the platform has “seated” into the acrylic and is flush with the bottom of the platform. As you tighten the screw you will feel the fitting draw up into the acrylic. Repeat this process with all (4) corners. Finally, place the new larger silicone mat onto the nesting platform. Reconnect your Belly Dancer® unit to an electrical power source and you will be ready to run.



SCIENTIFIC

7445 Chavenelle Road • Dubuque, IA 52002
(800) 253-4942 • (563) 690-0484

ibisci.com

Building
Quality
Scientific.