

## Water bath protective media AQUA STABIL

JULABO AQUA STABIL is a water disinfectant for open baths. It prevents the pollution of the bath fluid and stops the growths of algae, bacteria, and other microorganisms with its germicidal effectiveness. It increases the service life of water used for temperature control.

Application: Add 2 ml AQUA STABIL per 1 liter of water. Change the water every three months even if the blue coloration has not faded. If the coloration fades before three months, exchange the water and AQUA STABIL immediately.

Please observe the information in our safety data sheet. You can download the safety data sheet at [www.julabo.com](http://www.julabo.com) or request it from JULABO.



### Advantages:

- excellent heat conductivity
- minimum odor
- low corrosion tendency
- low toxicity

Property	Value
Working temperature °C	+20 ... +80

Order number	
8 940 006	6 bottles, 100 ml each
8 940 012	12 bottles, 100 ml each

Suitable for the following units:

- JULABO Heating Circulators (bath fluid: water)
- JULABO Water Baths
- JULABO Shaking Water Baths

-100 °C

0 °C

+100 °C