





Operating Instructions for POLYTRON[®] System PT 1200 E



Voltage



□ 100 - 120 V ± 10%, 60 Hz □ 230 V ± 10%, 50 Hz

Ensure that the supply voltage is correct and in agreement with the data on the rating plate

This is a quality product of



Luzernerstrasse 147a CH-6014 Littau- Lucerne Switzerland Tel.: +41-41-259 65 65 Fax: +41-41-259 65 75 e-mail: <u>laboratory@kinematica.ch</u>









1	INTRODUCTION	3
11	OPFRATING INSTRUCTIONS	3
11	1 RANGE OF VALIDITY	$\frac{J}{4}$
1.1	2 TARGET AUDIENCE	4
12	ORGANISATIONAL MATTERS	5
1.2	1 LOCATION OF THE OPERATING INSTRUCTIONS	5
1.2	2 MANUFACTURER AND CONTACT ADDRESS	5
1.3	WARNING NOTICES	6
2	SAFETY	7
21	SUMMARY	7
2.2	SAFETY CONCEPT	7
2.2	1 INTENDED USE OF THE EQUIPMENT	7
2.2	2 IMPROPER USE	8
2.2	3 USER ROLES	8
2.2	4 DANGER AREA	9
2.2	5 AREAS OF RESPONSIBILITY	9
2.2	.6 GENERAL SAFETY RULES	0
2.3	RESIDUAL DANGERS1	1
2.4	WARNINGS1	1
3	DESCRIPTION OF LABORATORY EQUIPEMENT	2
3.1	OVERVIEW	3
3.2	PT 1200 E DRIVE MOTOR 1	4
3.3	POLYTRON [®] dispersing equipment1	4
3.4	SPECIFICATIONS / TECHNICAL DATA	5
4	INSTALLATION	7
41	INPACKING 1	7
<u>4</u> 2	ASSEMBLING	7
4 3	COMMISSIONING 1	7
т.J		0
Э г 1		0
5.1		.8
5.2	HOUSING	.9 0
5.5	DISASSEMIBLY OF DISPERSTING UNITS	.9
6	IROUBLE SHOOTING	0
7	ACCESSORIES	0
•		1

GAT

PO

DN

PO

REA







1 INTRODUCTION

This chapter gives information on the the structure of this document. It will assist you in making use of it and show how to find the required information quickly.

1.1 OPERATING INSTRUCTIONS

Please read through these operating instructions before switching on or attempting to use the equipment. They describe the use of the POLYTRON[®] PT 1200 E, its installation and maintenance and the appropriate replacement parts and accessories. They will help you avoid erroneous use and consequent damage. Although POLYTRON[®] machines are designed for ease of service, this does not release you from the obligation to inspect your equipment carefully and to clean it thoroughly.

KINEMATICA AG is a specialist manufacturer of machines and equipment for dispersion and mixing technology.

An important objective of these operating instructions is to fully inform you, the user, about the correct and safe use of our equipment.

In order to achieve this, it is essential that you should carefully study chapter 2, "Safety", and follow the instructions in this book.

GA







1.1.1 RANGE OF VALIDITY

The information in these operating instructions relates to the $\text{POLYTRON}^{\text{®}}$ identified as follows:

Manufacturer:	KINEMATICA AG, CH-6014 Littau / Lucerne
Product name:	POLYTRON®
Type designation:	POLYTRON [®] PT 1200 E

Article number	Description
11010025	PT 1200 E, 230 V, (with EURO-plug))
11010026	PT 1200 E, 115 V, (with US-plug)
11010027	PT 1200 E, 230 V, (with CH-plug)
11010028	PT 1200 E, 230 V, (with GB-plug)
11010127	PT 1200 E, 230 V, with fixed 25'000 rpm (with EU-Plug)
11010127	PT 1200 E, 230 V, with fixed 25'000 rpm (with GB-Plug)
11030004	PT-DA 03/2EC-E050
11030012	PT-DA 05/2EC-E085 (PT-DA 1205/2EC-E)
11030031	PT-DA 07/2EC-E107 (PT-DA 1207/2EC-E)
11030062	PT-DA 12/2EC-E123 (PT-DA 1212/2EC-E)
11040006	HS 1200 E

1.1.2 TARGET AUDIENCE

These operating instructions are intended for all authorised users of our machines/equipment. We distinguish different user roles, taking account of the different demands placed on the user by the activity to be carried out.

EGATRO

You will find the definitions of user roles with the demands on the user in chapter 2, "Safety". You can fulfil one or more of these roles, provided that you meet the corresponding demands.





1.2 ORGANISATIONAL MATTERS

If you are unable to find the answer to any question in the operating instructions, please contact the equipment manufacturer directly.

1.2.1 LOCATION OF THE OPERATING INSTRUCTIONS

The operating instructions can only be of use to you if you always have them to hand. They should, therefore, always be kept at the place where the equipment is used.

1.2.2 MANUFACTURER AND CONTACT ADDRESS

KINEMATICA AG

Luzernerstrasse 147a CH-6014 Littau / Lucerne TEL: +41 41 259 65 65 FAX: +41 41 259 65 75

e-mail: <u>laboratory@kinematica.ch</u>

BA PT 1200 E English / Release 1.5 / 30.01.2009

G/1







1.3 WARNING NOTICES

Please be aware of the meaning of the following warning signs:



Safety instructions must be observed to ensure safe operation



This symbol indicates high voltage, with risk to health and environment

BA PT 1200 E English / Release 1.5 / 30.01.2009







2 SAFETY

This chapter is directed at all users of KINEMATICA laboratory equipment. It includes information on safe and optimum use.

2.1 SUMMARY

Any incorrect use of the installed equipment can be dangerous. Inadequately trained users can cause material damage and personal injury. This chapter informs you of the safety concept and the requirements for safe and optimum use of the equipment.

All those authorised to operate, service and repair the equipment are required to study chapter 2, "Safety".

2.2 SAFETY CONCEPT

The safety concept sets down the entitlement to use the equipment and the resposibilities of the individual users.

The machines and equipment are designed and constructed according to the state of the art and the recognised safety rules.

2.2.1 INTENDED USE OF THE EQUIPMENT

EGAT

The equipment is designed and constructed for the following use:

 Dispersion and homogenisation of pumpable fluid products in accordance with the technical specifications (see point 3.5) and compatibility with the materials coming into contact with the products.

If you use the equipment for any purpose other than those listed, the manufacturer cannot be held liable for any resulting damage.







2.2.2 IMPROPER USE

Any use other than the "proper use" without the written approval of the manufacturer or operation outside the technical limits of use is improper use.

2.2.3 USER ROLES

To guarantee safety, we place requirements on the users of the equipment that must be met without fail. Only persons meeting the requirements are authorised to work with the equipment.

We describe all those who work with the equipment as users. Since the requirements of these users are very much dependent on their activity, we ditinguish the following user roles.

Contract partner:

The manufacturer can impose legal obligations on the contract partner when the equipment is purchased. The contract partner is obliged to ensure that the equipment is properly used.

Operating company:

The operating company ensures that the equipment is properly used and authorises persons who are entitled to work with the equipmrnt in any one of the defined user roles. He is under the obligation to instruct the users.

Note:

Contract partner and operating company can be the same person.

EGATRON

Service technician:

The service technician is an employee of the operating company and looks after the equipment in special operating mode(s). He is a specialist with mechanical, electrical and electronic professional training. The service technician undertakes commissioning, decommissioning service and repair of the equipment. He must be appropriately trained to be able to carry out the service work required.

Operator:

The operator turns the equipment on and off. In the event of an alarm signal he informs the service technician.







2.2.4 DANGER AREA

System/equipment

The system danger area includes the whole system/equipment including the connecting lead and controls.

Proximity danger area

This refers to all areas within a defined distance of the equipment.

User danger area

This danger area includes all persons working with the equipment.

2.2.5 AREAS OF RESPONSIBILITY

In order that the system/equipment can be used safely and without risk, the users in various roles bear the responsibility for particular danger areas.

Contract partner:

The contract partner bears the responsibility for the "proximity danger area".

Operating company:

The operating company bears the responsibility for the "user danger area". Only those users may be authorised to operate the system/equipment who fulfil all requirements of the user roles concerned. In so doing, attention must be paid to the following points:

- It is to be ensured that all users of the system/equipment have fully read and understood chapter 2, "Safety" and act accordingly in a safety-conscious manner.
- It is to be ensured that no unauthorised person carries out work with the system/equipment.
- It is to be ensured that users are informed of the possible risks and dangers connected with the system/equipment.
- It is to be ensured that those being trained or engaged in general training are under the permanent supervision of a trained and authorised person.

EGATRON







Service technician:

The service technician bears the responsibility for the "system/equipment danger area". He ensures that the system/equipment is at all times free from technical faults, safe and functions correctly.

2.2.6 GENERAL SAFETY RULES

Observe the following general safety rules:

- follow these operating instructions,
- in addition, observe the legal obligations and requirements for accident prevention and environmental protection of the country in which you operate the equipment,
- do not make any modifications to the equipment without the written authorisation of the manufacturer,
- only original replacement parts may be used for repairs,
- before any service work on the equipment, it must be ensured that the electrical supply is switched off,
- after any service, maintenance or repair work has been carried out on the system/equipment, it must be given a test run by the service technician.
- depending on the place at which it is installed, circumstances may require that hearing protection is worn when remaining in the vicinity of the equipment for long periods.

EGATR







2.3 RESIDUAL DANGERS

When the system/equipment is used in accordance with rules and regulations, residual dangers are minimal.

Residual danger	Countermeasures
Tripping over feed or return lines	These should be appropriately laid
Noise	Wear hearing protection
Breakage of glass containers	Wear protective clothing (goggles etc.)
Tipping of the equipment	Stand it on a slip-proof, stable
	horizontal surface
Spitting of the product	Wear protective clothing (goggles etc.)
Rotating coupling	Never operate without attachment,
	never attempt to touch coupling or rotor

2.4 WARNINGS

•

•



- Ensure that the rated voltage of the equipment matches the supply.
- The equipment may only be opened by KINEMATICA AG authorised service specialists.
- The equipment may not be operated in areas where there is a danger of explosion.
- The PT 1200 E drive may only be operated using the PT 1200 power pack.







	•	The PT 1200 E is not qualified for continuous duty
	•	POLYTRON [®] dispersion attachments may not be operated dry – the lower sleeve bearing is cooled and lubricated by the medium being processed. Running dry will destroy the sleeve bearing.
	•	Dispersion attachments may only be changed with the power source disconnected.
	•	Ensure that the venting will never be covered by dust or any other objects. The efficiency of the motor will decrease.
	•	Ensure that dispersion attachments are thoroughly cleaned after every use.
	•	In the event that hazardous chemicals or materials that endanger health can influence the surroundings or use of the equipment, appropriate countermeasures must be taken.
	•	Ensure that before starting-up, no foreign objects are in the product .
	•	When the power supply is connected, never bring hands or tools within the vicinity of the coupling of the coupled dispersion attachment. Danger of injury by the rotating shaft and rotor.
	•	During operation, the sliding ring of the coupling may not be actuated; otherwise, the attachment may fall from the coupling.
	•	After longer use (15 min) the coupling could get very hot (over 40°C)
	•	Never turn your PT 1200 E upside down, if a dispersing unit with remains of fluid is clutched in. The fluid may damage the bearing of the motor.
	•	In the event of any power cut, the equipment must be switched off; otherwise, it may start again without proper supervision.
	•	The stator saw teeth present a danger of injury, as they are very sharp.
CE marking	•	KINEMATICA AG products comply with all the usual CE directives, carry the CE marking and are delivered with a corresponding declaration of conformity.

3 Description of laboratory equipment

ſ

BA PT 1200 E English / Release 1.5 / 30.01.2009

PO

Seite 12 von 22







POLYTRON[®] PT 1200 E Hand Unit is designed to homogenize, emulsify, blend and mix small quantities of organic and inorganic materials. The PT 1200 E conforms to all current and relevant CE regulations, carries the CE sign and is supplied with the relevant certificate of conformity.

3.1 OVERVIEW

The POLYTRON[®] PT 1200 E system produces fine dispersions, emulsions and homogenisations and is suitable for applications with working volumes up to 250 ml (depending on the product's viscosity and the dispersing equipment being used). The drive motor unit can be equipped with 4 different POLYTRON[®] dispersing units (dispersing head diameters from 3, 5, 7 and 12 mm) and they are easily exchanged via the quick release coupling.

For a workable system you require:

- the PT 1200 E drive motor with power pack
- an appropriate dispersing unit
- an electrical supply of 100 230 V, 50– 60 Hz

BA PT 1200 E English / Release 1.5 / 30.01.2009

GA







3.2 PT 1200 E DRIVE MOTOR

The PT 1200 E is fitted with a strong 24VDC-motor and a revolution setting unit.

The PT 1200 E with fixed 25'000 rpm (Ordner No. 11010127) has the setting knob too, but it's not active. The revolution is fixed according to the costumers request.

3.3 POLYTRON[®] dispersing equipment

ispersing units with diameters of 3, 5, 7 and 12 mm are available for the PT 1200 E drive motor. The illustrations below show the individual parts of a POLYTRON[®] dispersing unit. If you should have any further questions regarding the POLYTRON[®] dispersing equipment, do not hesitate to contact either the dealer of your choice or us.



PT-DA 07/2EC-E107



The unique quick release coupling makes it very easy to connect the dispersing unit to the POLYTRON[®] drive motor.

Simply slide the coupling ring upwards, push the dispersing unit into the coupling and then release the coupling ring. The dispersing unit is securely held in place.

Follow the reverse sequence to remove the unit, slide the coupling ring upwards and pull the dispersing unit out of the coupling.



Seite 14 von 22







Attention The drive motor must be disconnected from the power supply wh removing, exchanging or connecting a dispersing unit.

Ensure that the dispersing unit is well fixed in the coupling before activating the drive motor.

3.4 Specifications / Technical Data

PT 1200 E drive motor unit				
Type of motor	universal-motor with carbon brushes			
Voltage	100-120 V / 230 V ~			
Frequency	50 – 60 Hz			
Speed at no-load-operation	Approx. 25000 rpm			
Output Power	72 W			
Input Power	200 W			
Softstart	yes			
Noise Level	65 dB(A) at 25'000 min ⁻¹			
Rotatting Direction	clockwise (viewed from above)			
Ambient Temperature	-20 to 70°C			
Ambient Temperature on duty	0 to 40°C			
Relative Humidity	Max. rel. 95 %			
Protection Class	IP 20			
Max. period of continuous operation	Approx. 30 min at maximum load. Could be			
	even more, dependent upon load and			
	temperature.			
Dimensions	220 x 55 x 44 mm			
Weight	427 g			

	Dispersing units				
	PT-DA	PT-DA	PT-DA	PT-DA	
	03/2EC-E50	05/2EC-E85	07/2EC-E107	12/2EC-E123	
Shaft length,	50	85	107	123	
mm					
Stator/Rotor Ø,	2.2/3.5	5.5/3	7.8/5	12/9	
mm					
Processing	approx.	approx.	approx.	approx.	
volume	0.05 – 2 ml	0.1 – 3 ml	0.5 – 10 ml	3 – 250 ml	
Temperature	up to about 90 °C processing temperature				
Pressure	sure not pressurised				
Materials	stainless steel 1.4435 (316L) and PTFE compound				
Cleaning	can be sterilised by all the usual methods, e.g. autoclave				

BA PT 1200 E English / Release 1.5 / 30.01.2009

G

Seite 15 von 22







Product	The product to be dispersed must be pumpable and fluid and must
requirements	not contain any solid particles that might destroy the attachment.

BA PT 1200 E English / Release 1.5 / 30.01.2009

Seite 16 von 22







4 INSTALLATION

4.1 UNPACKING

Open the dispatch box and check that the contents agrees with the delivery note.



Check all parts for possible transport damage. Inform us or your dealer immediately of any disagreement or fault.

4.2 ASSEMBLING

After unpacking the PT 1200 E System is ready to use. You just have to clutch in the desired dispersing unit (see Chapter 3.4)

4.3 COMMISSIONING

Connect the drive unit to the electrical supply. **Check the supply voltage against the rating plate**.

- Immerse the dispersing unit in the product. The optimum depth of immersion of the dispersion unit is about 2/3 of the total depth of product in the container. Never immerse the upper holes of the dispersion attachment.
- Set the desired speed.

GA1







Note:

In some cases it is better to position the dispersion unit off centre in order to obtain more turbulence and better mixing. This can be achieved by tilting the drive by about 15°.

•	Never allow the POLYTRON [®] dispersion attachment to run without a liquid phase. Running dry will destroy the sleeve bearing.
•	Ensure that the dispersion attachment is thoroughly cleaned after every use.
•	Never touch the rotor of a POLYTRON [®] or BIOTRONA [®] dispersion attachment when rotating. There is great danger of injury.

5 SERVICE

Your POLYTRON[®] unit is designed for easy service. Nevertheless, it is essential to inspect your equipment carefully and to clean it thoroughly. Drawings of the separate components are to be found in appendix.



The equipment must be disconnected from the electrical supply:

- During any work on the equipment, in order to avoid any personal injury or other damage
- When changing or removing the dispersion attachment

5.1 DRIVE

Under normal working conditions, the PT 1200 E drive requires no servicing. Parts such as the motor armature and bearings are subject to natural wear.

We strongly recommend that service work and repairs should be carried out only by authorised KINEMATICA service centres or by KINEMATICA directly, where original replacement parts are available.

Any unauthorised modification or manipulation of the unit or its equipment leads to immediate annulment of the guarantee.

GA







5.2 Housing

The housing is made of styrene butadiene and meets best fire protection requirements UL-94-V0. Only suitable cleaners should be used.

5.3 Disassembly of dispersing units

PT-DA	A PT-DA PT-DA		PT-DA		
03/2EC-E050	05/2EC-E085	07/2EC-E107	12/2EC-E123		
With the univer	sal tool supplied	d, carefully	With this attachment you receive an		
knock the rotor	and attached s	haft out in the	additional tool. Use one tool to hold		
direction of the	coupling and th	en draw them	the shaft and the other to turn the		
out in the same	e direction.		rotor anti-clockwise and withdraw it.		
			You can then withdraw the shaft from		
			its tube.		
The lower sleev	ve bearing is slit	and can be	Using the tool, the sleeve bearing can		
opened and pu	lled over the sha	aft. The upper	be pushed out in the direction of the		
bearing can be	pulled off in the	direction of	coupling.		
the rotor.					
Replace defective parts. Bearings should			Replace defective parts.		
only be replace	ed in pairs.				
Reassemble in	the reverse ord	er.	Use the shaft to press the sleeve		
			bearing into the stator tube from the		
			coupling side. Screw on the rotor and		
			tighten gently.		
After every dis	assembly, and e	changing the bearings, a functional test			
	sho	ould be carried o	but in water.		

Criteria for changing the bearings:

Basically, it is the operating company that determines when and how often the bearings are to be changed. They should, however, be changed, at the latest, if:

- the rotor and stator come into contact,
- there is an increase in vibration,
- the rotor projects above the edge of the stator.

GΔ







6 TROUBLE SHOOTING

PROBLEM	CAUSE	C ORRECTIVE MEASURES
Unusual noises	Damaged drive bearings	Change ball bearings
	Damaged attachment	Change bearings
	bearings	Trace and replace defective
	Rotor/stator interference	parts (shaft, bearings)
Runs hot	Inadequate ventilation	Check ventilation slits are
	Damaged bearings	clear
	Overload	Change ball bearings
		Check manner of use
Vibrations	Bent shaft	Replace shaft
	Worn bearing(s)	Replace bearing(s)
	Defective coupling	Trace and replace defective
		parts
Drive is heating up after	Heating of electronics and	Turn the speed-controller to
longterm use.	motor	min so that drive can cool
		down.
Drive turns off during work.	Overload	Check dispersing unit.
	Electronic control causes	Check product.
	automatic power shut-off to	Turn off and on the drive.
	protect the drive.	(RESET)
Drive is heating up and	Thermal fuse has been	Send the equipment to the
turns off and cannot be	released.	nearest authorised
restarted anymore.	Overheating caused by	KINEMATICA AG service
	longterm overload use and	centre or directly to
	unsufficient cooling.	KINEMATICA AG
Drive runs stutteringly.	Carbon brushes burnt	Send the equipment to the
	Short circuit at motor anchor	nearest authorised
	Defective speed-controller	KINEMATICA AG service
		centre or directly to
		KINEMATICA AG

7 ACCESSORIES

A large selection of special homogenising containers is available, which can improve the efficiency of the processing of your product. Closed vessels to withstand vacuum or pressure can also be supplied.

Ask your dealer or contact KINEMATICA AG directly.

BA PT 1200 E English / Release 1.5 / 30.01.2009

Seite 20 von 22







8 GUARANTEE

KINEMATICA AG guarantees that the equipment that it produces will run free of any fault related to materials or manufacturing faults for **12 months**.

If thorough testing shows a fault to be due to either of the above causes, KINEMATICA AG guarantees that the equipment will be repaired or replaced free of charge.

The guarantee does not cover parts that are subject to normal wear. It is void if any person other than an employee of KINEMATICA AG or their appointed representative has made modifications to the equipment or if the damage is due to failure to comply with the operating instructions, to carelessness, accident, incorrect use or incorrect supply voltage.

KINEMATICA AG reserves the right to make technical changes to the equipment without modifying equipment delivered earlier in the same way.

In the event of technical problems, for spare parts requirements or for advice, contact our regional appointed agent, your prefered dealer or us directly at:

KINEMATICA AG

Luzernerstr. 147a CH-6014 Littau-Lucerne Switzerland Tel. +41-41-259 65 65 Fax +41-41-259 65 75 e-mail <u>laboratory@kinematica.ch</u>

BA PT 1200 E English / Release 1.5 / 30.01.2009

EGA

Seite 21 von 22





Appendix: Dimensional drawing and Art. No.



T.







Assemby Group	Main Identification	Order No.	Production-No.	Product-Identification	
	PT 1200 E, 230 V	11010025/27/28	9112017	PT1200E, 230V	
	PT 1200 E, 115 V	11010026	9112017	PT1200E, 115V	
Drive	PT 1200 E, 230 V, with	11010127	0150070	PT1200E, 230V. with	
Dive	fixed RPM	11010127	3130070	fixed RPM	
	PT 1200 E, 230 V, with	11010128	0158078	PT1200E, 230V. with	
	fixed RPM	11010120	3130070	fixed RPM	
	PT-DA 03/2EC-E050	11030004	9115386	PT-DA 03/2EC-E050	
	PT-DA 05/2EC-E085	11030012	9112005	PT-DA 05/2EC-E085	
Dispersing-Units	PT-DA 07/2EC-E107	11030031	9112007	PT-DA 07/2EC-E107	
	PT-DA 12/2EC-E123	11030062	9116012	PT-DA 12/2EC-E123	
	PT-DA 07/2SYN-E082	11030030	9753112	PT-DA 07/2SYN-E082	
Holder System	HS 1200 E	11040006	9753275	HS 1200 E	
Dispersing	GS 15 (35 ml)	11050010	9135018	GS 15 (35 ml)	
Vessels	GS 25 (150 ml)	11050011	9135019	GS 25 (150 ml)	
Furhter Accessories on request.					

BA PT 1200 E English / Release 1.5 / 30.01.2009

E

G

POL

Seite 22 von 22

PO