

Operating instructions

PLANETARY MONO MILL

PULVERISETTE 6 classic line

Valid starting with: 06.2000/03438



Read the instructions prior to performing any task!



Fritsch GmbH Milling and Sizing Industriestraße 8 D - 55743 Idar-Oberstein Telephone: +49 (0)6784/70-0 Fax: +49 (0)6784/70-11 email: info@fritsch.de Internet: www.fritsch.de



Certifications and CE conformity

Certifications and CE conformity

Certification

Fritsch GmbH has been certified by the TÜV-Zertifizierungsgemeinschaft e.V.





An audit certified that Fritsch GmbH conforms to the requirements of the DIN EN ISO 9001:2008.

CE Conformity

The enclosed Conformity Declaration lists the guidelines the FRITSCH instrument conforms to, to be able to bear the CE mark.





Table of contents

Table of contents

1	Basic structure	7
2	Safety information and use	8
	2.1 Requirements for the user	8
	2.2 Scope of application	8
	2.2.1 Operating principle	9
	2.2.2 Drive motor and speed regulation	9
	2.3 Obligations of the operator	9
	2.4 Information on hazards and symbols used in this manual 1	10
	2.5 Device safety information 1	13
	2.6 Protective equipment 1	14
	2.6.1 Opening the hood without mains connection 1	14
	2.6.2 Imbalance switch 1	14
	2.7 Hazardous points 1	16
	2.8 Electrical safety 1	17
	2.8.1 General information 1	17
	2.8.2 Protection against restart 1	17
	2.8.3 Overload protection 1	17
	2.8.4 Imbalance detection 1	17
3	Technical data 1	18
	3.1 Dimensions 1	18
	3.2 Weight 1	18
	3.3 Operating noise 1	18
	3.4 Voltage 1	18
	3.5 Current consumption 1	18
	3.6 Power consumption 1	18
	3.7 Electrical fuses 1	19
	3.8 Material 1	19
	3.9 Final fineness 1	19
4	Installation	20
	4.1 Transport 2	20
	4.2 Unpacking 2	20
	4.3 Setting up 2	21
	4.4 Ambient conditions 2	22
	4.5 Electrical connection 2	22
	4.5.1 Adjusting the mains voltage 2	23
	4.5.1.1 Adjusting the mains voltage with the rotary switch (9). 2	23
		<u>2</u> 4
		25
5	Initial start-up	26
		26
		26



Table of contents

6	Using the device	27
	6.1 Choice of grinding bowls and grinding balls	27
	6.1.1 Size of the grinding balls	29
	6.1.2 Number of balls per grinding bowl (independent of the	
	material quantity)	30
	6.1.3 Calculated weight of a ball	31
	6.2 Filling quantities of grinding bowls	31
	6.3 Filling the grinding bowl	32
	6.4 Factors with an impact on grinding	32
	6.4.1 Running time (grinding duration)	32
	6.4.2 Speed	32
	6.4.3 Reverse mode	32
	6.4.4 Number and size of the balls	32
	6.4.5 Weight of the balls (type of material)	33
	6.4.6 Dry grinding	33
	6.4.7 Wet grinding (grinding in a suspension)	33
	6.5 Clamping the grinding bowls	34
	6.5.1 Clamping with the "Safe-Lock" (5) tensioning device	34
	6.5.2 Clamping the 80 ml grinding bowls	36
	6.6 Mass balance	36
	6.7 Grinding duration	37
	6.8 Settings on the control panel	38
	6.8.1 Setting the speed	38
	6.8.2 Setting the running time	38
	6.8.2.1 Changing the time unit in setup mode	39
	6.9 Repetition of grinding / pause cycles	40
	6.10 Reverse mode	40
	6.11 Conducting a grinding operation	40
	6.11.1 Overload	41
	6.11.2 Switching off	41
	6.12 Cooling the grinding bowl	41
	6.13 Stand-by	41
7	Accessories	42
	7.1 Additional clamping system for grinding harmful sub-	
	stances or in a gas atmosphere	42
	7.1.1 Locking the additional clamping system into place	42
	7.2 Grinding in inert gas with gassing lid	45
	7.2.1 Preparation for gassing	46
	7.2.2 Gassing	47
	7.2.3 Ventilate after grinding	47
	7.2.4 Cleaning the valves	48
	7.2.5 Installing valve inserts	48
	7.3 GTM - system	49



Table of contents

8	Cleaning	50
	8.1 Grinding elements	50
	8.2 Mill	50
9	Maintenance	51
10	Repairs	53
	10.1 Checklist for troubleshooting	53
11	Examples of comminution tasks	55
12	Disposal	57
13	Guarantee terms	58
14	Exclusion of liability	60
15	Safety logbook	62
16	Index	65



Basic structure

1 **Basic structure**



- Counter weight 6
- 7 Lock

1

2

3

4

5

- 7 -

14 Support disc



2 Safety information and use

2.1 Requirements for the user

This operating manual is intended for persons assigned with operating and monitoring the Fritsch of the PULVERISETTE 6. The operating manual and especially its safety instructions are to be observed by all persons working on or with this device. In addition, the applicable rules and regulations for accident prevention at the installation site are to be observed. Always keep the operating manual at the installation site of the of the PULVERISETTE 6.

People with health problems or under the influence of medication, drugs, alcohol or exhaustion must not operate this device.

The of the PULVERISETTE 6 may only be operated by authorised persons and serviced or repaired by trained specialists. All commissioning, maintenance and repair work may only be carried out by technically qualified personnel. Qualified personnel are persons who, because of their education, experience and training as well as their knowledge of relevant standards, regulations, accident prevention guidelines and operating conditions, are authorised by those responsible for the safety of the machine to carry out the required work and are able to recognize and avoid possible hazards as defined for skilled workers in IEC 364.

In order to prevent hazards to users, follow the instructions in this manual.

Malfunctions that impair the safety of persons, the of the PULVERISETTE 6 or other material property must be rectified immediately. The following information serves both the personal safety of operating personnel as well as the safety of the products described and any devices connected to them: All maintenance and repair work may only be performed by technically qualified personnel.

This operating manual is not a complete technical description. Only the details required for operation and maintaining usability are described.

Fritsch has prepared and reviewed this operating manual with the greatest care. However, no guarantee is made for its completeness or accuracy.


Subject to technical modifications.

2.2 Scope of application

The PULVERISETTE 6 *classic line* can be applied universally for the fast, dry or wet grinding of inorganic and organic samples for analysis, quality inspection, material testing or mechanical alloying. During the synthesis, the PULVERISETTE 6 is used for mixing and homogenising dry samples, emulsions or suspensions.



2.2.1 Operating principle

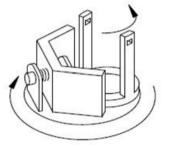

The grinding stock is crushed and ground by grinding balls in a grinding bowl. The centrifugal forces resulting from the rotation of the grinding bowls around their own axes and from the rotating support disc (14) influence the contents of the grinding bowl consisting of grinding stock and grinding balls.

The grinding bowl and the support disc (14) have opposite directions of rotation so that the influence of the centrifugal forces is alternately in the same direction and in the opposite direction. The result is that the grinding balls run down the inside of the bowl's wall as friction effect and the grinding balls hit the opposite wall of the grinding bowl as impact effect. The impact effect is amplified by the impact of the grinding balls against each other.

The loss-free comminution, even when grinding suspensions, is ensured by the hermetic seal between the grinding bowl and the lid.

- a Rotation of the grinding bowl
- b Centrifugal force
- c Movement of the support disc

2.2.2 Drive motor and speed regulation

A maintenance-free three-phase motor operated via a frequency converter is used as the drive.

2.3 Obligations of the operator

Before using the of the PULVERISETTE 6, this manual is to be carefully read and understood. The use of the of the PULVERISETTE 6 requires technical knowledge; only commercial use is permitted.

The operating personnel must be familiar with the content of the operating manual. For this reason, it is very important that these persons actually receive the present operating manual. Ensure that the operating manual is always near the device.

The of the PULVERISETTE 6 may exclusively be used within the scope of applications set down in this manual and within the framework of guidelines put forth in this manual. In case of non-compliance or improper use, the customer assumes full liability for the functional capability of the PULVERISETTE 6 and for any damage or injury arising from failure to fulfil this obligation.



By using the of the PULVERISETTE 6 the customer agrees with this and recognizes that defects, malfunctions or errors cannot be completely excluded. To prevent risk of damage to persons or property or of other direct or indirect damage, resulting from this or other causes, the customer must implement sufficient and comprehensive safety measures for working with the of the PULVERISETTE 6.

Neither compliance with this manual nor the conditions and methods used during installation, operation, use and maintenance of the of the PULVERISETTE 6 can be monitored by Fritsch GmbH. Improper execution of the installation can result in property damage and thus endanger persons. Therefore, we assume absolutely no responsibility or liability for loss, damage or costs that result from errors at installation, improper operation or improper use or improper maintenance or are in any way connected to these.

The applicable accident prevention guidelines must be complied with.

Generally applicable legal and other obligatory regulations regarding environmental protection must be observed.

2.4 Information on hazards and symbols used in this manual

Safety information

Safety information in this manual is designated by symbols. Safety information is introduced by keywords that express the extent of the hazard.



DANGER!

This symbol and keyword combination points out a directly hazardous situation that can result in death or serious injury if not avoided.



WARNING!

This symbol and keyword combination points out a possibly hazardous situation that can result in death or serious injury if not avoided.



CAUTION!

This symbol and keyword combination points out a possibly hazardous situation that can result in slight or minor injury if not avoided.



NOTICE!

This symbol and keyword combination points out a possibly hazardous situation that can result in property damage if not avoided.





ENVIRONMENT!

This symbol and keyword combination points out a possibly hazardous situation that can result in environmental damage if not avoided.

Special safety information

To call attention to specific hazards, the following symbols are used in the safety information:



DANGER!

This symbol and keyword combination points out a directly hazardous situation due to electrical current. Ignoring information with this designation will result in serious or fatal injury.



DANGER!

This symbol and keyword combination designates contents and instructions for proper use of the machine in explosive areas or with explosive substances. Ignoring information with this designation will result in serious or fatal injury.



DANGER!

This symbol and keyword combination designates contents and instructions for proper use of the machine with combustible substances. Ignoring information with this designation will result in serious or fatal injury.



WARNING!

This symbol and keyword combination points out a directly hazardous situation due to movable parts. Ignoring information with this designation can result in hand injuries.



WARNING!

This symbol and keyword combination points out a directly hazardous situation due to hot surfaces. Ignoring information with this designation can result in serious burn injuries due to skin contact with hot surfaces.



Safety information in the procedure instructions

Safety information can refer to specific, individual procedure instructions. Such safety information is embedded in the procedure instructions so that the text can be read without interruption as the procedure is being carried out. The keywords described above are used.

Example:



Close the lid carefully.

3. Tighten screw.

Tips and recommendations

Further designations

This symbol emphasises useful tips and recommendations as wells as information for efficient operation without malfunction.

To emphasise procedure instructions, results, lists, references and other elements, the following designations are used in this manual:

Designation	Explanation			
>	Step-by-step procedure instructions			
1., 2., 3				
⇔	Results of steps in the procedure			
Ŕ	References to sections in this manual and relevant documentation			
	Lists without a specific order			
[Button]	Operating elements (e.g. push button, switch), display elements (e.g. signal lamps)			
'Display'	Screen elements (e.g. buttons, function key assignment)			



2.5 Device safety information

Please observe!

- Only use original accessories and original spare parts. Failure to observe this instruction can compromise the safety of the machine.
- Accident-proof conduct is to be strictly followed during all work.
 Comply with all currently applicable pational and international accident.
- Comply with all currently applicable national and international accident prevention guidelines.



CAUTION!

Wear hearing protection!

If a noise level of 85 dB(A) is reached or exceeded, ear protection should be worn to prevent hearing damage.



WARNING!

The maximum accepted concentration (MAC) levels of the relevant safety guidelines must be observed; if necessary, ventilation must be provided or the machine must be operated under an extractor hood.



DANGER!

Explosion hazard!

- When grinding oxidizable substances, e.g. metals or coal, there is a risk of spontaneous combustion (dust explosion) if the share of fine particles exceeds a certain percentage. When grinding these kinds of substances, special safety measures must be taken and the work must be supervised from a specialist.
- The high-speed rotor mill is not explosion protected and is not designed to grind explosive materials.
- Do not remove the information signs.



NOTICE!

Immediately replace damaged or illegible information signs.

- Unauthorised alteration of the of the PULVERISETTE 6 will void Fritsch's declaration of conformity to European directives and void the guarantee.
- Only use the of the PULVERISETTE 6 when it is in proper working order, as intended and in a safety- and hazard-conscious manner adhering to the operating manual. In particular, immediately rectify any malfunctions that could pose a safety hazard.
- If, after reading the operating manual, there are still questions or problems, please do not hesitate to contact our specialised personnel.



2.6 Protective equipment



Protective equipment is to be used as intended and may not be disabled or removed.

All protective equipment is to be regularly checked for integrity and proper functioning.

For start-up, the hood (3) has to be closed.


The hood (3) is locked:

- without mains connection
- during operation



The hood (3) can only be opened, if the mill's drive is at standstill.

2.6.1 Opening the hood without mains connection

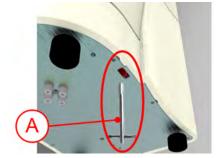

2.6.2 Imbalance switch

CAUTION!

The emergency release must not be activated while the machine is running! Disconnect the machine from the mains before the emergency release. Failure to observe this provision will void the guarantee, and releases us from liability for any resulting damage to the device or personal injury.

- **1.** Insert the included triangular key (A) into the bore hole on the bottom side of the PULVERISETTE 6 *classic line* and turn it to the right.
- 2. Unlock the latch (2) by pressing it.
- **3.** Pull the hood (3) up to open it.
- **4.** The mill can only be switched on again if the hood (3) is closed and the safety lock (7) has been enabled again by turning the triangular key to the right.

The device switches off if there is excessive imbalance. If this is the case, then the counterweight has to be readjusted.





CAUTION!

The imbalance switch can be disabled at your own risk. The Fritsch company will give no guarantee for damage resulting from disabling of the imbalance switch!



NOTICE!

Change these settings only after all work as described in & Chapter 4 'Installation' on page 20 has been carried out!




In the default setting of the imbalance switch, it is activated!

Activation / deactivation of the imbalance switch in Setup mode:

1. Press and hold the STOP button on the front control panel.



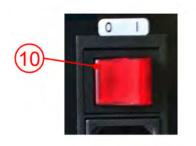


2. Switch on the device with the main switch (10) on the back side of the device, and release the STOP button.





3. If POWER SUPPLY is flashing, the device is in setup mode. If POWER SUPPLY is not flashing, repeat the procedure.

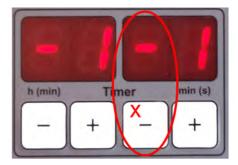



4.

NOTICE!

In the display above the right "-" button (x) in the TIMER field, there must be a minus sign. That activates the imbalance switch and prevents a drift of the device and possible damage to the device resulting from it.



5. To save and end setup mode, press the STOP button.

2.7 Hazardous points



CAUTION!

- Crushing hazard when closing the hood (3).

- Crushing hazard at the grinding bowl safe-lock tensioning device (5).
- Crushing hazard at the counterweight!

CAUTION! Risk of splashes!

During wet grinding overpressure may have been formed by the high temperature.

Wear protective goggles!





CAUTION! Risk of burning! The grinding bowls can be very hot after grinding.

Wear protective gloves!

2.8 Electrical safety

2.8.1 General information

- The main switch (10) separates the device from the mains on two poles.
- Switch off the main switch (10) if the planetary mono mill is down for a longer period of time (e.g. overnight).

2.8.2 Protection against restart

In case of power failure during operation or after switching off with the main switch (10), the hood (3) is locked. The hood lock (7) is opened when the power returns. For safety reasons, however, the mill does not restart.

2.8.3 Overload protection

- In the event of an overload, the device reduces the speed in a controlled manner. The REDUCED SPEED light is lit as a warning.
- The device switches off if the drive motor becomes too hot.
- The device switches off if the drive is blocked. (See ♦ Chapter 10 'Repairs' on page 53)

2.8.4 Imbalance detection

The device switches off if there is excessive imbalance. (See ♦ Chapter 10 'Repairs' on page 53)



Technical data

3 Technical data

3.1 Dimensions

500 x 370 x 530 mm (height x width x depth)

3.2 Weight

Net: 63 kg Gross: approx. 90 kg

3.3 Operating noise

Emissions value of workplace according to DIN EN ISO 3746:2005 is up to 85dB (A). The value fluctuates strongly, depending on the speed, the grinding stock, and the type of grinding bowl and grinding balls.

3.4 Voltage

The device can be operated in two voltage ranges:

- Single phase alternating current 100 120V ± 10%, and,
- Single phase alternating current 200 240V ± 10%.
 (See also Schapter 4.5 'Electrical connection' on page 22)
 Transient overvoltage according to overvoltage category II is permitted.

3.5 Current consumption

Depending on the mains voltage, the maximum current consumption is in the ranges:

- 100 120V→ 12 A
- 200 240V \rightarrow 7 A

3.6 Power consumption

Depending on the voltage range, the maximum power consumption is approx. 1.1 kW.



Technical data

3.7 Electrical fuses

- Circuit breaker (12): 2 x 10 A T
- Micro-fuse 10 A T in the frequency converter

3.8 Material

- Maximum feeding size approx. 10 mm
- Maximum feeding amount 225 ml

3.9 Final fineness

- **D**ry grinding up to d_{50} < 20 μ m (depending on the material)
- Wet grinding up to $d_{50} < 1 \ \mu m$ (depending on the material)



4 Installation

4.1 Transport

The device is delivered on a transport pallet with a wooden cover. We recommend using a forklift or pallet truck for transporting the packed device.



DANGER! Do not step under the transport pallet during transport.



Improper lifting can lead to personal injury or property damage. The machine is only to be lifted with suitable equipment and by qualified personnel.

The guarantee excludes all claims for damage due to improper transport.

4.2 Unpacking

- Remove the nails that fasten the wooden hood to the transport pallet.
- Lift the hood off the transport pallet.
- The pre-perforated segments can be detached so that the foam parts can be removed more easily.
- Please store the transport packaging so that it can be reused if you need to return the product. Fritsch GmbH accepts no liability for damage caused by improper packaging (packaging that is not from Fritsch).
- Compare the contents of the delivery with your order.



Grinding bowls made of tempered steel may have recesses on the surface caused during production. They do not have an impact on grinding or the grinding results and usually disappear after the first grinding operation.

These recesses on the surface, if present, are within the range of the permissible production tolerances. Complaints relating to such grinding bowls therefore cannot be accepted.



4.3 Setting up



DANGER!

Do not step under the transport pallet during transport.



CAUTION!

The weight of the planetary mono mill is approx. 67 kg!



CAUTION! Crushing hazard!


Always lift with 2 persons.

Hold the bottom edge of the housing when lifting the device.




CAUTION!

Never operate the planetary mono mill while it is standing on the transport pallet!

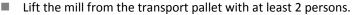

- Place the mill on a flat, stable surface. A fastening set is provided to help stabilise the mill and prevent the mill from shifting during operation. Screw the two rings diagonally into the surface using the drilling template. Place the 2 device feet of the mill into the intended rings.
- Make sure that the planetary mono mill can be easily accessed. There has to be sufficient space to reach the main switch on the back side of the device.
- Keep the air outlet on the side ventilation grate free. Risk of overheating!

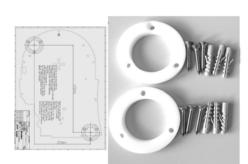



4.4 Ambient conditions



WARNING! Mains voltage!

- The device may only be operated indoors.
- The surrounding air may not carry any electrically conductive dust.
- Maximum relative humidity 80% for temperatures up to 31°C, linearly decreasing down to 50% relative humidity at 40°C.
- The room temperature has to stay between 5 40°C.
- Altitudes up to 2000 m
- Degree of pollution 2 according to IEC 664.

4.5 Electrical connection



DANGER! Provide short-circuit protection!

Risk of damage due to short-circuits.

 Make sure that the socket is connected to a mains line protected with a residual current circuit breaker.



DANGER!

Mains voltage!

Changes to the connection line may only be made by a qualified person.

CAUTION!


Ignoring the values on the type plate may result in damage to the electrical and mechanical components.

Before establishing the connection, compare the voltage and current values stated on the type plate with the values of the mains system to be used.

1. Plug the supplied power cord into the port (11) at the back of the device.

2. Then connect the device to the mains using the power cord!



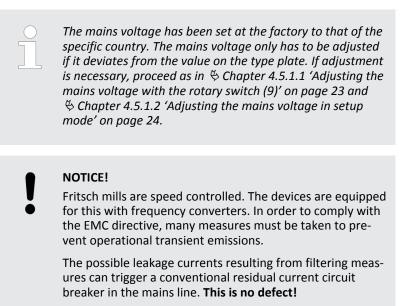

To prevent this, special residual current circuit breakers, which are adapted for operation with frequency converters, are commercially available.

Operation without a residual current switch is possible, but must be done in accordance with the relevant regulations.

4.5.1 Adjusting the mains voltage

4.5.1.1 Adjusting the mains voltage with the rotary switch (9)



CAUTION!

Only qualified personnel may change the voltage range on the device!



CAUTION!

The voltage range may only be adjusted after the mains has been disconnected. Disconnect the device from the mains!






- **1.** Disconnect the device from the mains!
- **2.** The rotary switch (9) for adjusting the mains voltage is located on the back side of the device. Rotate this switch to the required voltage.
- **3.** Connect the device to the mains.

4.5.1.2 Adjusting the mains voltage in setup mode



1. Press and hold the STOP button on the front of the control panel.

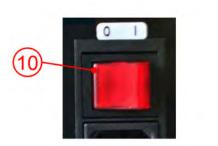

- 2. Switch on the device with the main switch (10) on the back side of the device, and release the STOP button.
- **3.** If POWER SUPPLY is flashing, the device is in setup mode. If POWER SUPPLY is not flashing, repeat the procedure.



4. Use the +/- ROTATIONAL SPEED buttons to adjust the level of the mains voltage (90-260 V) to the existing mains system.







5. To save and end setup mode, press the STOP button.

4.6 Setting device specification in setup mode



NOTICE!

"P6" must always be displayed in the REPETITIONS field. The Fritsch company will give no guarantee for damage resulting from disabling of the imbalance switch.




Initial start-up

5 Initial start-up

Perform initial start-up only after all work as described in \mathcal{G} Chapter 4 'Installation' on page 20 has been carried out.

5.1 Switching on

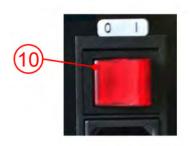

5.2 Function check

- The device must be connected to the power supply if this has not been done already.
- Switch on the device with the main switch (10) on the back side of the device.
- The POWER SUPPLY lamp lights up.



CAUTION!

Only perform the function check at a speed of 100 1/min. The counterweight must be positioned completely inwards to prevent too great an imbalance.

- Open the hood (3).
- Take out the grinding bowl tensioning device (Safe-Lock) and transport securing device (wooden block). No loose parts may remain inside the device.
- Position the counterweight completely inwards. (See 'Mass balance' on page 36).
- Close the hood (3).
- Set the speed to 100 1/min. (See <a> Chapter 6.8.1 'Setting the speed' on page 38).
- Press START on the control panel.
- The hood (3) is locked and the mill starts up at the preselected speed.





6 Using the device



CAUTION!

Before starting the machine, make sure that the grinding bowl has been tensioned correctly and that there are no loose parts inside the device. There is a risk of loose grinding bowls or parts being projected. Failure to observe this will render void the guarantee, and releases us from liability for any resulting damage to the device or personal injury.

NOTICE!

During grinding, the temperatures in the grinding bowl may get very high.

In encased grinding bowls, the inserts are glued into the casing with a two-component construction adhesive.

The adhesive is resistant to temperatures up to approx. 140 °C. Above 140 °C, the adhesive will liquefy and accumulate below the insert in the casing. When the adhesive cools down, it solidifies and pushes the casing up. That can cause irreparable damage to the insert. The grinding bowl will definitely be rendered unusable.

Above temperatures of 200 °C, the adhesive will be destroyed. The same applies for encased grinding bowl lids.



V-belt, seals, and motor require some time at the beginning before they can reach the optimal output and/or rotation. This means that initially a well-filled and heavy grinding set may be able to operate with less rotations than after an introductory phase of approx. 1 - 2 hours running time.

6.1 Choice of grinding bowls and grinding balls



CAUTION!

If the grinding elements used are not genuine accessories, we assume no guarantee and exclude all liability for damage to the device or for personal injury.





CAUTION!

The grinding element is subject to normal wear when used. Before every grinding operation, check the wall thickness of the grinding bowls. In the event of severe wear, replace the grinding bowl. If this is not done, the prevailing high centrifugal forces during grinding may cause the grinding balls to penetrate the bowl's wall and damage the mill. Failure to observe this will render void the guarantee and release us from liability for any resulting damage to the device or personal injury.

The hardness and density (specific weight) of the grinding bowl and grinding balls used must be greater than that of the material used to prevent excessive wear by abrasion.

Material (bowl and balls)	Main components of the material	Density in g/cm³ High density means high impact energy!	Abrasion resistance	Use for grinding stock
Agate	(99.9% SiO ₂)	2.65	Good	Soft to medium-hard samples
Silicon nitride	(90% Si ₃ N ₄)	3.25	Extremely good	Abrasive samples, metal-free grinding
Sintered corundum	(99.7% Al ₂ O ₃)	3.9	Fairly good	Medium-hard, fibrous samples
Zirconium oxide	(96,2% ZrO ₂)	5.7	Very good	Fibrous, abrasive samples
Stainless steel	Bowl: (17-19% Cr + 8-10% Ni) Balls: (12.5-14.5% Cr + 1% Ni)	7.8	Fairly good	Medium-hard, brittle samples
Tempered steel	Bowl: (11-12% Cr) Balls: (1.0-1.65% Cr)	7.9	Good	Hard, brittle samples
Tungsten carbide	(93% WC+6% Co)	14.9	Very good	Hard, abrasive sam- ples

The grinding bowls and grinding balls made of zirconium oxide are resistant to acids - apart from hydrofluoric acid.

Normally choose a grinding bowl and grinding balls that are made of the same material.



Exception: Tungsten carbide balls (<20 mm) may be temporarily (a few minutes) combined with grinding bowls made of tempered steel.

6.1.1 Size of the grinding balls

Type of feed material	Suitable ball diameter		
Hard samples with a maximum feed size of 10 mm	30 mm or 40 mm		
Average feed size of < 5 mm	20 mm		
Fine material < 0.5 mm	10 mm or 5 mm		
Homogenisation of dry or liquid samples	10 mm		
Homogenisation of viscous samples	20 mm		

These are reference values: The size of bowls and grinding balls may need to be determined through experimentation.



NOTICE!

It is not advisable to mix balls of different diameters. (If balls with different diameters are used, increased wear to the balls is to be expected.)



NOTICE!

Balls with a diameter of 40 mm are rarely used for grinding, because this ball size can cause damage to the grinding bowl fast if the grinding duration is too long. Only use grinding balls of this diameter for brief grinding durations.



6.1.2 Number of balls per grinding bowl (independent of the material quantity)

A higher number of balls will reduce the grinding time and the grinding result will have a smaller particle size distribution.

Ball diameter (mm)	Grinding bowl volume (ml)	80	250	500
5	Number of balls (pcs)	250 - 300	1200 - 1300	2000 - 2500
10	Number of balls (pcs)	25 -30	50 - 150	100 - 250
15	Number of balls (pcs)	10	45 - 50	70 - 100
20	Number of balls (pcs)	5	15 - 20	25 - 35
30	Number of balls (pcs)	-	5 - 6	10
40	Number of balls (pcs)	-	-	4

These are reference figures: The number of balls may need to be determined through experimentation.



CAUTION!

When grinding with a ball size of Ø 30 mm or Ø 40 mm, do not let the device run unsupervised. The vibrations may lead to shifting.



6.1.3 Calculated weight of a ball

Ball diameter in mm		5	10	15	20	30	40
Material	Density in g/cm ³	Calculated weight of a ball in g					
Agate	2.65	0,17	1,39	4,68	11,1	37,46	88,8
Silicon nitride	3.25	0,20	1,7	5,48	12,99	43,83	103,88
Sintered corundum	3.9	0,25	1,99	6,72	15,92	53,72	127,34
Zirconium oxide	5.7	0,37	2,99	10,07	23,88	80,58	191,01
Stainless steel	7.8	0,51	4,08	13,78	32,67	110,27	261,38
Tempered steel	7.9	0,52	4,14	13,96	33,09	111,68	264,73
Tungsten car- bide	14.9	0,96	7,70	25,98	61,58	207,82	492,6

To determine the weight of the required balls, the "calculated weight of a ball" is multiplied by the "number" of balls required.

Example: A 250 ml agate bowl is to be filled with 1221 agate balls with a diameter of 5 mm.

Calculation: 0.17g * 1221 St \approx 208 g

208 g of grinding balls can be weighed and inserted in the grinding bowl, thus avoiding the time required for counting the balls.

6.2 Filling quantities of grinding bowls



Never operate the mill without grinding stock! This can lead to grinding balls and grinding bowls getting damaged.


NOTICE!

If the minimum filling quantity is fallen short of, increased wear due to abrasion is to be expected. This can cause irreparable damage to the mill components.



Grinding bowl	min. sample volumes	max. sample volumes
500ml	80ml	225ml
250ml	30ml	125ml
80ml	1ml	30ml

6.3 Filling the grinding bowl

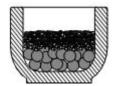

- Do not fail to comply with the following sequence:
- **1.** Place the grinding balls in the empty bowl.
- 2. Fill grinding stock onto the balls.
- **3.** Insert the seal ring.
- **4.** Place the lid on the grinding bowl.

6.4 Factors with an impact on grinding

6.4.1 Running time (grinding duration)

A longer grinding time will increase the percentage of fine material. To reduce the grinding time, you can use a grinding bowl and grinding balls with a higher density, and thus a higher impact energy.

6.4.2 Speed

Higher speeds shorten the grinding time and increase the share of fine particles.

6.4.3 Reverse mode

- Useful for mechanical alloying
- Improvement of the homogeneity of the sample

6.4.4 Number and size of the balls

Pre-grind course, hard material with large balls:

reduced percentage of fine material!

Many small balls increase the percentage of fine material during extended running time.



6.4.5 Weight of the balls (type of material)

A higher mass (specific weight) of the grinding balls accelerates grinding. (see table in \bigotimes Chapter 6.1 'Choice of grinding bowls and grinding balls' on page 27).

6.4.6 Dry grinding

Below a particle size of approx. 20 μm the surface forces prevail and the material to be ground starts to "stick".

Additional dry comminution can be achieved by adding surface-active substances to the material to be ground.

Examples (maximum amount to be added in mass%)

- Stearic acid 2-3%
- Aerosil (fine-particle silicic acid) 0.5-2%
- Quartz sand ~ 2%
- Glass powder ~ 2%

6.4.7 Wet grinding (grinding in a suspension)



DANGER!

Explosion hazard! Ignition hazard!

The device is not explosion-protected. If flammable liquids are used, make sure that the heat developing in the grinding bowl does not reach the solvent's boiling point. Program appropriate cooling phases. If the vapour pressure is too high, vapours may escape and ignite. If it can be avoided, we recommend using non-flammable liquids or liquids with a high boiling point. The boiling point should be above 80 °C and above 100 °C for a long grinding duration.

During the transition to grinding in suspension, you can add a liquid auxiliary agent with high boiling point and low vapour pressure.



6.5 Clamping the grinding bowls

6.5.1 Clamping with the "Safe-Lock" (5) tensioning device

Fritsch GmbH confirms that each Safe-Lock clamping system is produced and tested according to our internal quality standards without any exception. Fritsch GmbH measures the tension force of each individual Safe-Lock - clamping system. The tension force must be in the range 11.2 – 11.3kN +/- 0.2kN to pass our Quality Control.

Carry out the following checks for every grinding before clamping the grinding bowls:


- Has the rubber disc been inserted in the grinding bowl holder? The rough side should be facing upwards!
- Check the rubber disc in the grinding bowl holder for damage: Replace the rubber disc if it has been pressed flat!
- The Teflon flat seal (for the sealing between lid and bowl) may not be damaged or soiled.

Replace heavily deformed Teflon flat seals.

- The surfaces of the lids and bowls on which the Teflon flat seals lie must be clean!
- Check the rubber disc of the pressure piece for damage!
 Replace rubber discs that have been pressed flat and are protruding laterally from the pressure piece.

Clamping

1. Position the bracket (a) of the clamping lever vertically!

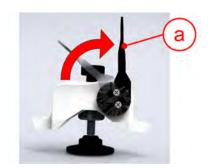

2. Hang the longer projection of the Safe-Lock tensioning device into one side of the mounting of the grinding bowl holder.





3. Hang the short projection of the Safe-Lock tensioning device into the other side.

4. Push the Safe-Lock tensioning device so that the mounting of the grinding bowl holder sits centrally in the U-shaped cut-out of the longer projection.



5. Pre-tension the setting screw (b) manually, then place the provided torque spanner on the setting screw (b) and turn until it clicks.



6. Then press the clamping lever downwards and fixate the Safe-Lock tensioning device.



DANGER!

After several minutes of grinding and in the cooling-down phases, check that the tensioning device is firmly connected.




If the Safe-Lock system is correctly tensioned, the clamping lever is automatically pulled downwards by the eccentric when horizontal.



6.5.2 Clamping the 80 ml grinding bowls

0

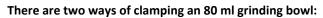


1. Clamp the 80 ml grinding bowl with mounted reducer (o) (order no. 90.1120.09).



2. Alternatively, clamp two 80 ml grinding bowls on top of one another.

6.6 Mass balance

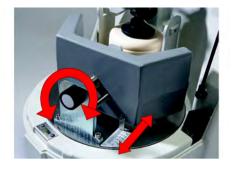

The counterweight (6) has to be positioned according to the scale in order to offset the imbalance and thus prevent the mill from shifting.

The weight indicated on the scale is the weight of the filled grinding bowl with the lid on.

NOTICE!

It may be necessary to readjust the mass balance if the bowls are heavy or the speeds are high.

NOTICE!

Extra weights, like "GTM" and "additional clamping system", have to be balanced as well.



6.7 Grinding duration



WARNING! Burn hazard!

Grinding bowls can get very hot after long grinding durations. Wear protective gloves for removal after grinding or during the grinding breaks.

Depending on the application, the grinding duration should be adapted to the development of heat in the bowls. The temperature inside the bowls is 20 - 30 °C warmer than the outer casing temperature.




CAUTION!

The maximum temperature of the grinding bowl outer casing is 100 - 110 °C (agate, max. 70 - 80 °C).

The grinding duration is therefore based on the maximum bowl temperature. The grinding duration at which the temperature is not exceeded depends on the material, ball, and speed. For this reason, the user should determine it through experimentation.

Reference value

When grinding at high speeds and with large bowls, a grinding duration of one hour (depending on the temperature) should not be exceeded. Then allow the unit to cool down for half an hour to one hour.

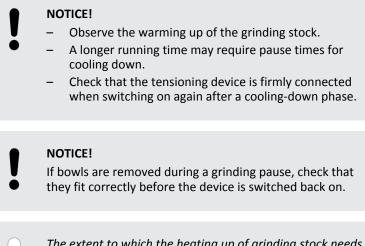



The extent to which the heating up of grinding stock needs to be observed depends of course on the sample used in each individual case. Note[®] A longer running time may also require a longer pause time for cooling down.



To reduce the grinding time, you can use a grinding bowl and grinding balls with a higher density, and thus a higher impact energy.

The mill can also run for several hours during low-speed operations for mixing and homogenisation.

Operation with an external time switch is not possible.

6.8 Settings on the control panel

6.8.1 Setting the speed



- Switch on the main switch (10) on the back side of the device (I).
- The green POWER SUPPLY ready status indicator lights up on the control panel.



ROTATIONAL SPEED control panel area

Press and hold the + or - button.

The speed can be selected in steps of 10 between 100 and 650 1/min.

The actual speed is displayed during operation. The nominal speed is briefly displayed when the + or - button is pressed.

6.8.2 Setting the running time




TIMER control panel area

- Press the MILLING button, it lights up; press the + or button and select the running time in hours (0..99) and minutes (0..59).
- Set the pause time if cooling down is necessary. Press the "Pause" button, it lights up; press the + or - button and select the pause time in hours (0..99) and minutes (0..59).

If no pause time is required, set the pause time to 0.




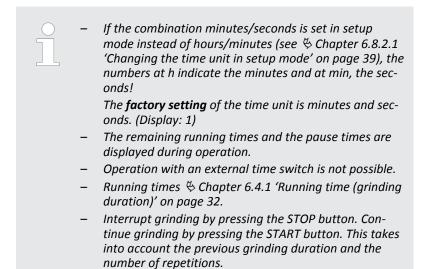

6.8.2.1 Changing the time unit in setup mode



1. When the device has been switched off, press and hold the STOP button on the front control panel.

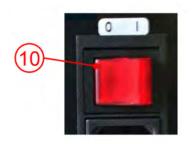

- **2.** Switch on the device with the main switch (10) on the back side of the device.
- **3.** If POWER SUPPLY is flashing, the device is in setup mode. If POWER SUPPLY is not flashing, repeat the procedure.



+ -	pulverisette
Rotational Speed	Repetitions -
	Power Supply Stand By Reduced Speed
Milling Pause Reverse	START Power On STOP

4. To perform changes, press the right "+" button (y) in the TIMER field:

Time unit, hours and minutes \rightarrow Display: -

Time unit, hours and minutes \rightarrow Display: 1 (factory setting)

Press + or - button; select the number of repetitions (0..99). The number

5. To save and end setup mode, press the STOP button.

6.9 Repetition of grinding / pause cycles



6.10 Reverse mode



 \rightarrow Press the REVERSE button. The button lights up.

of remaining cycles is displayed during operation.

REPETITIONS control panel area

After the selected running time expires, the mill will change its direction of rotation. For this, REPETITIONS has to indicate at least 1.

"Reverse" is selected when mixing dry samples, for example.

6.11 Conducting a grinding operation



- After the conditions specified in Chapter 6 'Using the device' on page 27 have all been observed, close the hood (3).
- Press the START button on the control panel.
- The hood is locked and the mill starts up.
- The mill turns at the set speed (nominal speed). If the load is too great, for example due to heavy grinding bowls, the speed is reduced (actual speed) so that the machine is not overloaded.
 → If the mill does not start, see Chapter 10 'Repairs' on page 53



While in operation, the hood (3) remains locked, even during pause times, and the fan cools the interior.



6.11.1 Overload

If the mill is overloaded, the speed is reduced and the REDUCED SPEED light flashes.

The mill switches off if the overload continues for too long; see *Chapter 10.1 'Checklist for troubleshooting' on page 53*.

6.11.2 Switching off

- Press STOP on the control panel.
- When the drive comes to a standstill, the hood is unlocked and can be opened.
- If the device is not in operation for a long time, switch off the main switch (10) on the back side.

6.12 Cooling the grinding bowl



WARNING! Burn hazard!

Grinding bowls can get very hot after long grinding durations. Wear protective gloves for removal after grinding or during the grinding breaks.

- When the hood is open (3), or
- In the programmed pause times with closed (locked) hood and the fan running.

6.13 Stand-by

If the mill is not in operation and the hood (3) is open, it switches to the energy-saving stand-by mode after one hour. The STAND-BY lamp lights.

The stand-by function is not possible when the hood is closed.



7 Accessories

7.1 Additional clamping system for grinding harmful substances or in a gas atmosphere.

The additional clamping system is used to transport a grinding bowl filled with inert gas or harmful substances from a glove box to the planetary mill and back again. This ensures that no harmful substances can be inhaled.

Grinding bowls with standard lids have to be filled in a glove box with an inert gas atmosphere and closed using the additional clamping system without being clamped in the mill.




Grinding bowls with gassing lids can also be closed and gassed outside the device using the additional clamping system.

7.1.1 Locking the additional clamping system into place




1. Place the grinding bowl in the additional clamping system.



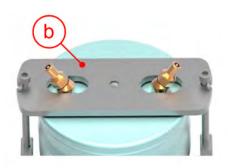

2. Position the pressure plate with rubber disc on the bowl as shown in the picture.

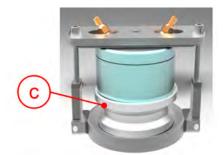

3. When using grinding bowls with a volume of 80 ml, the adapter piece (90.1120.09) has to be used as well.



NOTICE!

Under no circumstances may 2 x 80 ml bowls be clamped on top of one another. This can lead to damage to the bowl and its component seals.



4. Then secure the pressure plate on both sides with the socket-head screws shown in the picture and a hex key.




5. Place the additional clamping system with secured bowl in the grinding bowl holder.





- **6.** The pressure piece has to be removed in order to use the Safe-Lock system for tensioning.
- **7.** Clamp the additional clamping system with the Safe-Lock tensioning device into the device.

Clamping

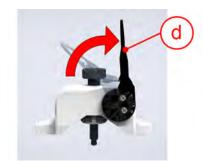

1. Position the bracket (a) of the clamping lever vertically!



2. Hang the longer projection of the Safe-Lock tensioning device into one side of the mounting of the grinding bowl holder.



3. Hang the short projection of the Safe-Lock tensioning device into the other side, and insert the setting screw of the tensioning device in the recess of the additional clamping system.





4. Push the Safe-Lock tensioning device so that the mounting of the grinding bowl holder sits centrally in the U-shaped cut-out of the longer projection.

- e
- 5. Pre-tension the setting screw (b) manually, and screw it tight using the provided torque spanner. Place the torque spanner on the setting screw (b) and turn until it clicks.



6. Then press the clamping lever downwards. Re-tighten the additional clamping system with a hex key. The additional clamping system may be loosened due to the clamping of the Safe-Lock using the clamping lever.



DANGER!

After several minutes of grinding and in the cooling-down phases, check that the tensioning devices are still firmly connected.



If the Safe-Lock system is correctly tensioned, the clamping lever is automatically pulled downwards by the eccentric when horizontal.

7.2 Grinding in inert gas with gassing lid



NOTICE!

Observe imbalance offsetting! (See \Leftrightarrow Chapter 6.6 'Mass balance' on page 36).



Wa Th ap bu spa

We carry out a worker water bath test on all gassing lids. The part to be tested is sealed, a pressure of 5.5 bar is applied and it is immersed in a water bath. If there is a leak, bubbles will develop. The air bubbles that develop within a specific interval are evaluated by the worker/tester.

Only gassing lids with a leak rate of $<10^{-4}$ [mbar l/s] are approved.

When grinding in inert gas, the same conditions apply regarding clamping and composition of grinding set and balls, as apply for standard grinding.

Two valves are screwed onto the gassing lid through which you can feed in inert gas (e.g. nitrogen) before switching on the mill. A Viton flat seal is used instead of a Teflon one.

7.2.1 Preparation for gassing



Grinding bowls outside the mill can be filled with gas using the additional clamping system.

- Gassing hose
- b Hose clamp
- c Coupling
- d Valves

а

- e Ventilation attachment
- Fill the grinding bowl with grinding balls and grinding stock.
- Put on lid with seal (use Viton seal).
- Insert the grinding bowl (closed with an additional clamping system) into the grinding bowl holder (8).
- Lock the grinding bowl with gassing lid or grinding bowl with additional clamping system in the device.

Using the additional clamping system, the following steps can also be completed in the glove box with subsequent locking into the planetary mono mill:

- Connect the gassing hose (a) to the inert gas supply using the provided hose clamp (b).
- Screw the ventilation attachment (e) onto one of the two valves (d).
- Place the coupling (c) of the gassing hose on the free valve. When doing so, press the lever of the coupling and push the coupling along the neck of the valve until it goes no further. Release the lever.



7.2.2 Gassing

- Slowly open the inert gas supply.
- Press a thin object (e.g. hex key) onto the top of the ventilation attachment (e) so that the air can escape from the grinding bowl.
- The inert gas now purges the air from the grinding bowl.
- The duration of purging has to be determined through experimentation.

It depends on grinding bowl size, filling, and gas supply, among other factors.

- To end purging, close the inert gas supply and release the ventilation attachment.
- Screw off the ventilation attachment.
- Remove the gassing hose. To do so, press the lever.



CAUTION!

Only switch on the device when both coupling and ventilation attachment have been removed.

Overpressure may occur during grinding!

7.2.3 Ventilate after grinding




CAUTION!

Always let the bowl cool down before ventilation. Hot gases and sample material may escape from the bowl during pressure equalisation, leading to serious burns.

- \rightarrow Use protective gloves when ventilating!
- When the bowl has cooled down, remove the additional clamping system with bowl from the device.
- Screw the ventilation attachment onto the valve.
 Each valve can be used for aerating or ventilating.
- For pressure equalisation (of the overpressure occurring because of grinding), carefully press on the ventilation attachment with a thin object (e.g. hex key).
- Only now should you release the additional clamping system (5)!



7.2.4 **Cleaning the valves**

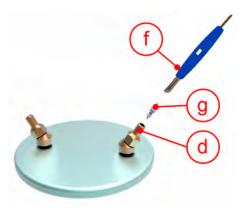

- Valve screwdriver f
- g d Valve insert
- Valves

Both valves (d) should be cleaned after every grinding!

- Insert the thin end of the valve screwdriver (f) from above into the valve (d) and turn anti-clockwise.
- Screw out the valve insert (g).
- Depending on the soiling, clean the valve insert (g) with compressed air, or place it in a small glass container filled with alcohol and clean in an ultrasonic cleaner (LABORETTE 17) and then dry carefully.
- After the two valve inserts have been removed, the two valve holders can be cleaned with compressed air from above the lid.

7.2.5 Installing valve inserts

- Insert the valve insert (g) into the valve (d) with the spring pointing upwards.
- With the valve screwdriver (f), screw the valve insert clockwise.

The following gassing lids for the grinding sets are available, each with two valves and a soft sealing ring:

Material	Order number
Hardmetal tungsten carbide 250 ml	50.8600.00
Tempered chrome steel 80 ml	50.8700.00
Tempered chrome steel 250 ml	50.8500.00
Tempered chrome steel 500 ml	50.8400.00
Stainless steel 80 ml	50.8800.00
Stainless steel 250 ml	50.8300.00
Stainless steel 500 ml	50.8200.00
Agate 250 ml	50.8100.00
Agate 500 ml	50.8000.00



NOTICE!

The soft black sealing rings made of "Viton" can endure temperatures of approx. 200 °C.

The valves (d) can endure temperatures of approx. 180 °C for one hour at most.



NOTICE!

The grinding parts made of agate are only designed for temperatures of up to 100 °C. After this point, they need to be slowly and carefully cooled down.

7.3 GTM - system

The GTM system is available as an accessory for recording pressure and temperatures in the grinding bowl during grinding.

Instructions are included with the GTM system.



Cleaning

8 Cleaning



DANGER!

Mains voltage!

- Before beginning with cleaning work, disconnect the mains plug and protect the device against being unintentionally switched back on!
- Do not allow any liquids to flow into the device.
- Indicate cleaning work with warning signs.
- Put safety equipment back into operation after cleaning work.



When cleaning the entire device, adhere to the guidelines of the Accident Prevention Regulation (BGV A3) - especially when the device has been set up in a dusty environment or when the grinding stock processed produces dust.

8.1 Grinding elements

- Clean the grinding bowl and grinding balls each time after using them: Clean them, e.g., under running water using a brush and a commercially available cleaning agent.
- Half fill the grinding bowl with grinding balls, sand and water, and run for 2 to 3 minutes (correctly tensioned) in the planetary mono mill.
- Cleaning with an ultrasonic cleaner is permitted.
- For sterilisation in the heat cabinet, only heat the grinding elements up to 100 °C.


8.2 Mill

The planetary mono mill can be wiped down with a damp cloth when it is switched off.



Maintenance

9 Maintenance

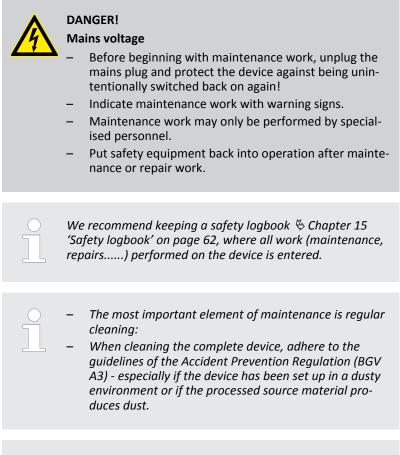



NOTICE!

Safe-lock - clamping systems might lose their tension force over time of usage. To watch the right tension force of each Safe-lock - clamping system Fritsch GmbH stipulates a scheduled maintenance interval of one year. Please send back all Safe-lock - clamping systems for regular checks to Fritsch GmbH. This service is liable to pay the costs.

Functional part	Task or description	Test	Maintenance interval
Safety lock	Hood lock (7)	Is the closed hood (3) locked in place when the main switch is off? If this test is failed, do not continue to work until the fault has been rectified.	Before each use
Rotating bearings	Permanent lubrication	Bearing clearance	Every 2,000 h or annually



Maintenance

Functional part	Task or description	Test	Maintenance interval	
Drive motor	Permanent lubrication	Bearing clearance	Every 4,000 h or annually	
V-belt	Motor planetary disc	Check belt tension Disconnect the device from the mains. Screw off the (top) rear cover plate. The belt must not slacken by more than 10 mm when pushed with your thumb.	Once a year	
Fan, ventilation slots	Grinding chamber cooling and electronics	Clean functional parts when soiled	Twice a year	
Grinding bowl tensioning device, Safe-Lock (5)	Rubber of the pressure piece and rubber disc in the grinding bowl holder	Signs of use; replace if pressed flat and no longer elastic	After every 1,000 h	
	Indicated surfaces of Safe- Lock are round Original height, 12.5 mm Minimum height, 12 mm (y)	Safe-Lock has frequently been tensioned too loosely	After every 200 h If the height is under the minimum height, the corre- sponding part has to be replaced.	
Grinding bowl lid seal	nding bowl lid seal Seal the grinding bowl Replace seal if dirt has penetrated		After every 100 h	
Grinding bowl holder (8)	Keep tension Original height, 14mm	Minimum measurement at grinding bowl holder, 11 mm (x) If the height is under minimum height, the sponding part has to replaced.		
(y)	min. 12mm	x min. 7	I1mm	

Abb. 1

Abb. 2



Repairs

10 Repairs



DANGER!

Mains voltage!

- Before beginning with repair work, unplug the mains plug and protect the device against being unintentionally switched back on.
- Indicate repair work with warning signs.
- Repair work may only be performed by specialised personnel.
- Put safety equipment back into operation after maintenance work.

10.1 Checklist for troubleshooting

Fault description	Cause	Remedy
Mill does not start up	No mains connection	Plug in mains plug.
	Main switch (10) at 0 (OFF)	Switch on main switch.
	Device fuse	Check device fuse; fuse insert on the back side of device $2 \times 10 \text{ A T}$ (12).
"START" button is pressed but	Check if POWER SUPPLY is lit	See 🖔 Chapter 4 'Installation' on page 20.
mill does not start up	Pause time active	Wait for end of pause or press "STOP".
	Safety lock (7) was opened man- ually	See $\stackrel{<}{\Rightarrow}$ Chapter 2.6.1 'Opening the hood without mains connection' on page 14.
Mill reduces speed automati- cally	If REDUCED SPEED is lit: Over- load	Reduce load or accept automatically set speed.
Mill stops running	Switched off due to thermal overload of the drive	Allow device to cool down and select a lower speed.
	Imbalance of machine too high	Improve mass balance. (See & <i>Chapter 6.6</i> ' <i>Mass balance' on page 36</i>).
	Drive was blocked	Rectify malfunction in grinding chamber.
	Motor's V-belt loose or ripped	Check V-belt, replace if necessary.
	Speed sensor is defective	Contact customer service.
The hood cannot be opened	The button (1) on the front of the hood was not actuated when opening.	Actuate button when opening.



Repairs

Fault description	Cause	Remedy	
	No mains connection	Plug in mains plug.	
	Main switch (10) at 0 (OFF)	Switch on main switch (I).	
	Device fuse	Check device fuse; fuse insert on the back side of device $2 \times 10 \text{ A T}$ (12).	
Grinding stock escapes	Tensioning device (5) loose	Check and re-tension if necessary.	
	Seal ring is soiled or defective	Clean or replace seal ring.	
Runs unevenly with strong vibrations	Masses balanced inadequately	Set mass balance more adequately: See	



Examples of comminution tasks

11 Examples of comminution tasks

Material			
Feeding amount	Material of grinding bowl and balls	Grinding balls, St. x diam- eter	Result
Feed size	Grinding bowl size	Grinding duration	Final fineness
Ruby (stone)			
140 g	Cr-Ni steel	6 x 30 mm	100%
12 mm	250 ml	3 min	< 250 μm
Titanium dioxide TiO ₂ (dry a	nd wet grinding in water)		
40 g	Cr-Ni steel	6 x 30 mm	100%
2 mm	250 ml	30 min	< 40 µm
Titanium dioxide TiO ₂ (wet g	rinding in water)		
40 g / 50 ml water	Cr-Ni steel	6 x 30 mm	100%
2 mm	250 ml	60 min	< 10 µm
Coal (dry and wet grinding in	n water)		
5 g	Zirconium oxide	5 x 20 mm	100%
0.5 mm	80 ml	120 min	< 15 µm
Aluminium oxide / silicon ox	ide		
100 g	WC + Co	15 x 20 mm	90%
0.1 mm	250 ml	90 min	< 20 μm
Ferrovanadium			
70 g	WC + Co	5 x 30 mm	70%
3 mm	250 ml	20 min	< 100 µm
Glass			
50 g	Agate	15 x 20 mm	100%
4 mm	250 ml	15 min	< 90 µm
Silicon carbide (dry and wet	grinding in water)		
15 g	WC + Co	5 x 20 mm	100%
3 mm	80 ml	30 min	< 150 μm
Silicon carbide (dry and wet	grinding in water)		
15 g / 5 ml water	WC + Co	5 x 20 mm	100%
3 mm	80 ml	45 min	< 71 µm



Examples of comminution tasks

Raw phosphate					
40 g	Cr steel	15 x 20 mm	100%		
3 mm	250 ml	2 min	< 250 μm		
Manganese dioxide MnO ₂ (v	Manganese dioxide MnO ₂ (wet grinding in water)				
50 g / 40 ml water	WC + Co	15 x 20 mm	100%		
0.1 mm	250 ml	60 min	< 20 µm		
Sludge (dry)					
180 g	Al ₂ O ₃	10 x 30 mm	100%		
8 mm	500 ml	30 min	< 250 μm		
Active carbon (wet grinding	in water)				
150 ml	Cr-Ni steel	15 x 20 mm	100%		
0.025 mm	250 ml	30 min	< 5 μm		
Plaster					
300 g	Cr steel	10 x 30 mm	100%		
10 mm	500 ml	20 min	< 200 µm		
Protein					
50 g	Sintered corundum 1	6 x 30 mm	90%		
20 mm	250 ml	90 min	< 50 µm		
Grains (barley)					
100 g	Sintered corundum 1	3 x 40 mm	100%		
3 mm	500 ml	20 min	< 150 µm		
Dough products					
100 g	Sintered corundum 1	10 x 30 mm	100%		
5 mm	500 ml	3 min	< 250 μm		
Sugar (wet grinding in alcoh	ol)				
200 g	Agate	10 x 30 mm	100%		
1 mm	500 ml	45 min	< 10 µm		



Disposal

12 Disposal

It is hereby confirmed that FRITSCH has implemented the directive 2002/95/EC of the European Parliament and Council from 27th January 2003 for the limitation of the use of certain dangerous substances in electrical and electronic devices.

FRITSCH has registered the following categories according to the German electrical and electronic equipment act, section 6, paragraph 1, clause 1 and section 17, paragraphs 1 and 2:

Mills and devices for the preparation of samples have been registered under category 6 for electrical and electronic tools (except for large stationary industrial tools).

Analytical devices have been registered under category 9, monitoring and control instruments.

It has been accepted that FRITSCH is operating only in the business-tobusiness area. The German registration number for FRITSCH is WEEE reg. no. DE 60198769

FRITSCH WEEE coverage

Since the registration of FRITSCH is classified for bilateral transactions, no legal recycling or disposal process is described. FRITSCH is not obliged to take back used FRITSCH devices.

FRITSCH declares it is prepared to take back used FRITSCH devices for recycling or disposal free of charge whenever a new device is purchased. The used FRITSCH device must be delivered free of charge to a FRITSCH establishment.

In all other cases FRITSCH takes back used FRITSCH devices for recycling or disposal only against payment.



Guarantee terms

13 Guarantee terms

Guarantee period	As manufacturer, FRITSCH GmbH provides – above and beyond any guar- antee claims against the seller – a guaranty valid for the duration of two years from the date of issue of the guarantee certificate supplied with the device.
	Within this guarantee period, we shall remedy all deficiencies due to material or manufacturing defects free of charge. Rectification may take the form of either repair or replacement of the device, at our sole discre- tion. The guarantee may be redeemed in all countries in which this FRITSCH device is sold with our authorisation.
Conditions for claims against the guar- antee	This guarantee is subject to the condition that the device is operated according to the instructions for use / operating manual and its intended use.
	Claims against the guarantee must include presentation of the original receipt, stating the date of purchase and name of the dealer, together with the complete device type and serial number.
	For this guarantee to take effect, the answer card entitled "Securing of Guarantee" (enclosed with the device) must be properly filled out and
	despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information.
Reasons for loss of the guarantee	despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <i>online registration</i> must be
Reasons for loss of the guarantee	despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information.
Reasons for loss of the guarantee	 despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information. The guarantee will not be granted in cases where: Damage has arisen due to normal wear and tear, especially for wear parts, such as: Crushing jaws, support walls, grinding bowls, grinding balls, sieve plates, brush strips, grinding sets, grinding disks, rotors, sieve rings, pin inserts, conversion kits, sieve inserts, bottom sieves, grinding inserts, cutting tools, sieve cassettes, sieve and measuring
Reasons for loss of the guarantee	 despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information. The guarantee will not be granted in cases where: Damage has arisen due to normal wear and tear, especially for wear parts, such as: Crushing jaws, support walls, grinding bowls, grinding balls, sieve plates, brush strips, grinding sets, grinding disks, rotors, sieve rings, pin inserts, conversion kits, sieve inserts, bottom sieves, grinding inserts, cutting tools, sieve cassettes, sieve and measuring cell glasses. Repairs, adaptations or modifications were made to the device by
Reasons for loss of the guarantee	 despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information. The guarantee will not be granted in cases where: Damage has arisen due to normal wear and tear, especially for wear parts, such as: Crushing jaws, support walls, grinding bowls, grinding balls, sieve plates, brush strips, grinding sets, grinding disks, rotors, sieve rings, pin inserts, conversion kits, sieve inserts, bottom sieves, grinding inserts, cutting tools, sieve cassettes, sieve and measuring cell glasses. Repairs, adaptations or modifications were made to the device by unauthorized persons or companies. The device was not used in a laboratory environment and/or has
Reasons for loss of the guarantee	 despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information. The guarantee will not be granted in cases where: Damage has arisen due to normal wear and tear, especially for wear parts, such as: Crushing jaws, support walls, grinding bowls, grinding balls, sieve plates, brush strips, grinding sets, grinding disks, rotors, sieve rings, pin inserts, conversion kits, sieve inserts, bottom sieves, grinding inserts, cutting tools, sieve cassettes, sieve and measuring cell glasses. Repairs, adaptations or modifications were made to the device by unauthorized persons or companies. The device was not used in a laboratory environment and/or has been used in continuous operation. Damage is present due to external factors (lightning, water, fire or
Reasons for loss of the guarantee	 despatched without delay after receipt of the device and be received by us within three weeks or alternatively, <u>online registration</u> must be carried out with the above-mentioned information. The guarantee will not be granted in cases where: Damage has arisen due to normal wear and tear, especially for wear parts, such as: Crushing jaws, support walls, grinding bowls, grinding balls, sieve plates, brush strips, grinding sets, grinding disks, rotors, sieve rings, pin inserts, conversion kits, sieve inserts, bottom sieves, grinding inserts, cutting tools, sieve cassettes, sieve and measuring cell glasses. Repairs, adaptations or modifications were made to the device by unauthorized persons or companies. The device was not used in a laboratory environment and/or has been used in continuous operation. Damage is present due to external factors (lightning, water, fire or similar) or improper handling. Damage is present that only insubstantially affects the value or



Costs not covered by the guarantee	This guarantee excludes any costs for transport, packaging or travel that accrue in the event the product must be sent to us or in the event that one of our specialist technicians is required to come to your site. Any servicing done by persons not authorised by us and any use of parts tha are not original FRITSCH accessories and spare parts will void the guar- antee.	
Further information about the guar- antee	The guarantee period will neither extend nor will a new period of guar- antee begin in the event that a claim is placed against the guarantee.	
	Please provide a detailed description of the type of error or the com- plaint. If no error description is enclosed, we shall interpret the shipment as an assignment to remedy all recognisable errors or faults, including those not covered by the guarantee. Errors or faults not covered by the guarantee shall in this case be rectified at cost.	
	We recommend reading the operating manual before contacting us or your dealer, in order to avoid unnecessary inconvenience.	
	Ownership of defective parts is transferred to us with the delivery of the replacement part; the defective part shall be returned to us at buyer's expense.	
	NOTICE! Please note that in the event that the device must be returned, the device must be shipped in the original Fritsch packaging. Fritsch GmbH denies all liability for any damage due to improper packaging (packaging not from Fritsch).	

Any enquiries must include a reference to the serial number imprinted on the type plate.



Exclusion of liability

14 Exclusion of liability

Before using the product, be sure to have read and understood this operating manual.

The use of the product requires technical knowledge; only commercial use is permitted.

The product may be used exclusively within the scope of applications set down in this operating manual and within the framework of guidelines put forth in this operating manual and must be subject to regular maintenance. In case of non-compliance, improper use or improper maintenance, the customer assumes full liability for the functional capability of the product and for damage or injury arising from violating these obligations.

The contents of this operating manual are subject in entirety to copyright law. This operating manual and its contents may not be copied, further distributed or stored in any form, in part or in whole, without the prior written consent of Fritsch.

This operating manual has been prepared to the best of our knowledge and checked for accuracy at the time of printing. FRITSCH GMBH assumes no guarantee or liability whatsoever for the accuracy or completeness of the contents of this operating manual, including but not limited to the implied warranties of merchantability and fitness for a particular purpose, unless liability is expressly prescribed by applicable laws or jurisprudence.

FRITSCH GMBH expressly reserves the right to modify and/or update this operating manual without prior notice. The same applies to modifications and improvements to the products described in this operating manual. It is the responsibility of the user to ensure that they have the current version of this operating manual. For more information, please contact your local FRITSCH GMBH distributor or Fritsch GmbH, Industriestr. 8, D-55473 Idar-Oberstein.

Not all parts shown here are necessarily installed in the product. The buyer is not entitled to delivery of these parts. If interested, please contact your local FRITSCH GMBH distributor or Fritsch GmbH, Industriestr. 8, D-55743 Idar-Oberstein.

FRITSCH GMBH takes the greatest care to ensure that the quality, reliability and safety of your products are continuously improved and adapted to the state of the art. The supplied products as well as this operating manual conform to the current state of the art when they leave the sphere of influence of FRITSCH GMBH.

By using the product the customer agrees with this and recognizes that defects, malfunctions or errors cannot be completely excluded. To prevent risk of damage to persons or property or of other direct or indirect damage, resulting from this or other causes, the customer must implement sufficient and comprehensive safety measures for working with the product.



Exclusion of liability

Fritsch GmbH excludes any liability, warranty, or other obligation to compensate for damages, regardless of whether this liability, warranty, or other obligation is explicit or implicit, contractual or arising from unlawful acts or prescribed contractually, by law, or otherwise. In no event shall the buyer be entitled to any compensation from Fritsch GmbH for any special, direct, indirect, coincidental or consequential damage, including but not limited to lost profits, lost savings, lost sales or financial loss of any kind or for compensation of third parties, for downtimes, for lost goodwill, for damage to or replacement of equipment and property, for costs or restoration of materials or goods related to the product or the use of our products, for other damage or injury to persons (including fatal injuries) or similar. The above exclusion of liability is limited by mandatory liability as prescribed by laws or jurisprudence. Liability for negligence is excluded in all cases.

No permission is given expressly, implicitly or otherwise for the use of patents, brands or other copyrights. We also assume no liability for copyright infringements or infringements of the rights of third parties arising from the use of this product.

Neither compliance with this operating manual nor the conditions and methods used during installation, operation, use and maintenance of the product can be monitored by Fritsch GmbH. Improper execution of the installation can result in property damage and thus endanger persons. Therefore, we assume absolutely no responsibility or liability for loss, damage or costs that result from errors at installation, improper operation or improper use or improper maintenance or are in any way connected to these.



Safety logbook

15 Safety logbook

Date	Maintenance / Repair	Name	Signature



Safety logbook

Date	Maintenance / Repair	Name	Signature



Safety logbook

Date	Maintenance / Repair	Name	Signature



Index

16 Index

Α

Accident prevention
Additional clamping system
Adjusting to the mains voltage
Authorised persons

С

Clamping the 80 ml grinding bowls36Clamping with the Safe-Lock (5) tensioning device34Cleaning the valves48Conducting a grinding operation40Cooling the grinding bowl41Current consumption18

D

Dimensions	
Disposal	
Dry grinding	

Ε

Electrical safety
Examples of comminution tasks
Exclusion of liability
Explanation of signs
Explanation of symbols

F

Filling grinding bowls
Filling quantities of the grinding bowls
Final fineness

G

Grinding duration
Grinding in inert gas
GTM - system
Guarantee terms

Н

Hazardous points
Imbalance detection17Initial start-up26Installing valve inserts48

Μ

Maintenance	
Material	

Ν

Number of balls per bowl	
--------------------------	--

0

Opening the hood without mains connection 14
Operating noise
Operating principle
Overload
Overload protection

Ρ

Power consumption	3
Preparation for gassing	5
Protection against restart	/
Protective equipment	ł

R

Repairs	,
Repetition of grinding / pause cycles 40	l
Requirements for the user	
Reverse mode	l
Running time	

S

Safety information 10 Safety logbook 62, 63, 64
Scope of application
Selecting the grinding balls
Selecting the grinding bowl
Setting the running time
Setting the speed
Setting up
Size of the grinding balls
Skilled workers
Speed
Stand By - Mode

Т

Troubleshooting checklist .			•	•	•	•			. !	53

U

Unpacking	 •				•	 		•		•	•	•	 	2	0
v															

Ventilate after grinding 47 Voltage 18

W

Warning information	
WEEE	



Index

Wet grinding	33
--------------	----





© 2014 Fritsch GmbH Milling and Sizing Industriestraße 8 D - 55743 Idar-Oberstein Telephone: +49 (0)6784/ 70-0 Fax: +49 (0)6784/ 70-11 email: info@fritsch.de Internet: www.fritsch.de