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Extent of delivery

- schuett petriturn-E
- Power supply
- Operating instructions
- Warranty Card with reply card

Warranty

The manufacturer grants a one-year guarantee on the schuett petriturn-E.

To establish your right for warranty coverage, send in the reply card accompanying the Warranty Card to the manufacturer within 14 days.

Device faults and malfunctions resulting from improper use are not covered by the warranty.

schuett petriturn-E

The schuett petriturn-E is an electrically driven petri-dish turntable for uniform distribution of liquid suspensions of micro organisms onto agar plates. The turning speed is continuously adjustable from 20 to 120 rpm to fit the needs of the user's particular application. The turntable is started manually by either a touch-free sensor or by a foot-switch (optionally). The turning time may be adjusted from 2 s to 3 min or continuously.

Instructions for set-up

Carefully take the schuett petriturn-E out of its shipping package and place the device in its intended location. Compare the data on the power supply type plate with the available line/mains voltage. Connect the power supply with the line/mains.

Controls

All controls are located on the front of the schuett petriturn-E.

The turning knob on the right-hand side marked "on/off – speed" is suitable for adjusting the speed continuously.

The turning knob on the left-hand-side marked "running – time" is suitable for adjusting the turning time and activating the sensor.

The jacks for the power supply and the foot-switch are located at the back of the device.

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Putting into operation

Please follow the order of installation.

Connect the cable of the power supply to the corresponding jack at the schuett petriturn-E (backside). Connect the power supply with the line/mains. Insert the requested petri dish for inoculation into the turntable of the schuett petriturn-E.

Sensor operation (without foot-switch):

For sensor operation, the foot-switch should not be connected.

1. Make sure both turning knobs are in the "0" position.
2. Adjust the needed turn time by moving the turning knob "running – time" clockwise (6 = approx. 3 minutes).
3. Turn the turning knob "on/off – speed" to the right in order to start the device.
4. Activate the turntable by moving your hand in front of the sensor.
5. Adjust the requested speed with the turning knob "on/off – speed".

Continuous operation (without foot-switch):

For continuous operation:

1. Make sure both turning knobs are in the "0" position.
2. Turn the turning knob "on/off – speed" to the right in order to start the device.
3. Adjust the requested speed with the same turning knob.
4. Leave the turning knob "running – time" in the "0" position.

Operation with foot switch:

Tip: In order to limit the quantity of cables on your working place, connect the cable of the power supply with the corresponding jack at the foot-switch and connect the cable of the foot-switch to the corresponding jack at the schuett petriturn-E (backside).

1. Make sure both turning knobs are in the "0" position.
2. Adjust the needed turn time by moving the turning knob "running – time" clockwise (6 = approx. 3 minutes).
3. Turn the turning knob "on/off – speed" to the right.
4. Press the foot-switch. The turntable of the schuett petriturn-E starts turning at the set speed as long as the foot-switch is pressed.
5. The speed may be adjusted continuously by using the turning knob "on/off – speed".

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Operation with petri dishes of up to 150 mm Ø

For operating the schuett petriturn-E with petri dishes of up to 150 mm diameter insert the corresponding turntable attachment (optional) to the turntable.

Cleaning

For cleaning the schuett petriturn-E, you may use standard, soft cloths and detergents. The silicone disk in the turntable may be removed and autoclaved. Also the adapter may be autoclaved. The turntable itself may be removed from the base frame for easier cleaning by pulling upright.

Accessories	Cat.-No.
schuett petriturn-E foot-switch made of stainless steel	3.361 302
schuett petriturn-E turntable attachment for petri-dishes of up to 150 mm Ø (autoclavable)	3.361 402
inoculating hook, stainless steel	3.362 152
inoculating hook, glass	3.362 162

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Trouble shooting

Description of fault	Measures to eliminate fault
----------------------	-----------------------------

A simple fault may often be the reason for a malfunction, and as such can easily be corrected by the user himself.

After turning the knob „on/off – speed“ to the right hand side and activating the sensor, the turntable does not start turning.

- Check line/mains connections and voltage.

When operated with a foot-switch the turntable does not turn although the foot-switch is pressed.

- Check the cable connection of the power supply and the foot-switch.
- If the schuett petriturn-E works properly in the sensor mode, replace the foot-switch.

If the measures listed above do not restore the device to its operational state or if the fault is not listed above, please contact the manufacturer.

Technical Data

Dimensions (Ø x H)	160 x 70 mm
Weight	approx. 1 kg
Revolutions direct adjustable	20-120 rpm
Power	12 VDC, 5 VA